



Mounting the transfer connection to the K-377

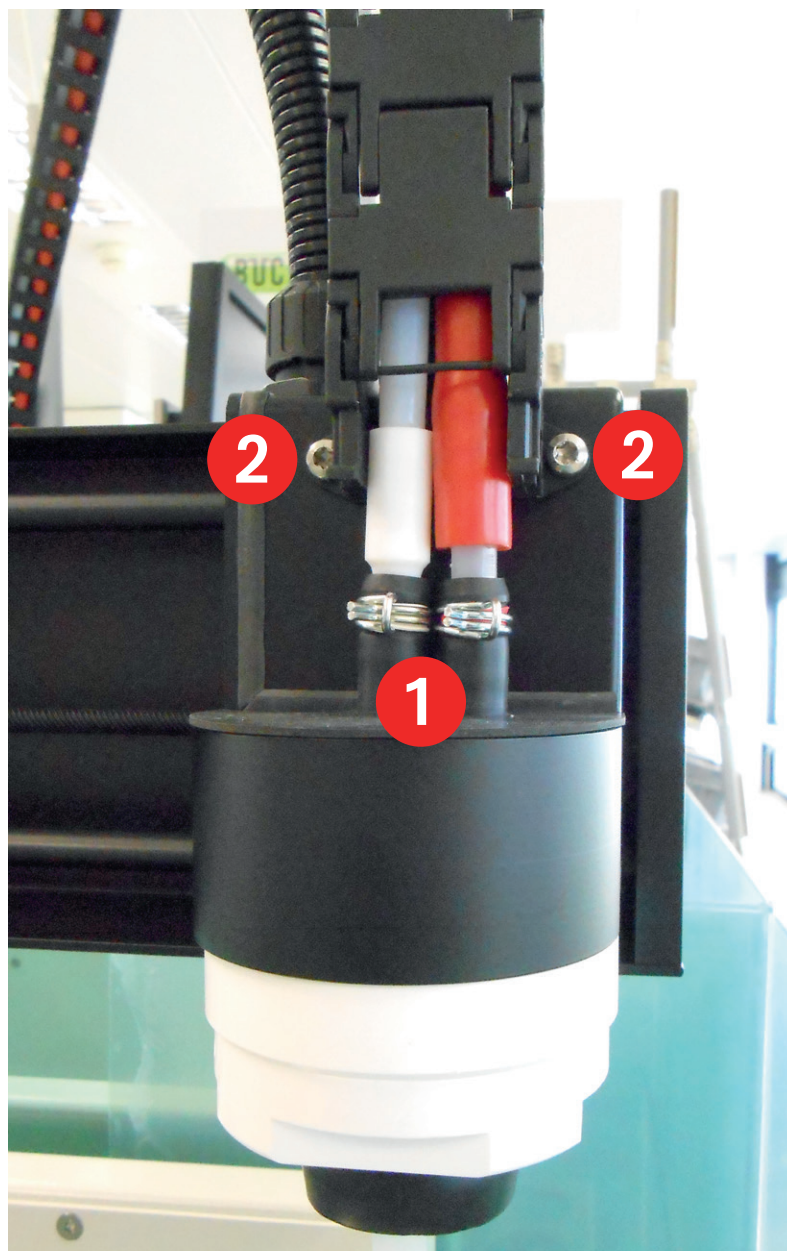

Picture 1: Attaching the transfer connection to the K-377 KjelSampler

- Connect the transfer hose and the steam hose to the two fittings on top of the sealing cap on the sampler arm. Secure both connections with hose clamps ①.
The red steam hose has to be fixed to the first position (marked with a red ring) - pointing to the front of the instrument!
- Fix the plastic cable channel with the two provided screws ② to the sampler arm.

Mounting the transfer connection to the K-375

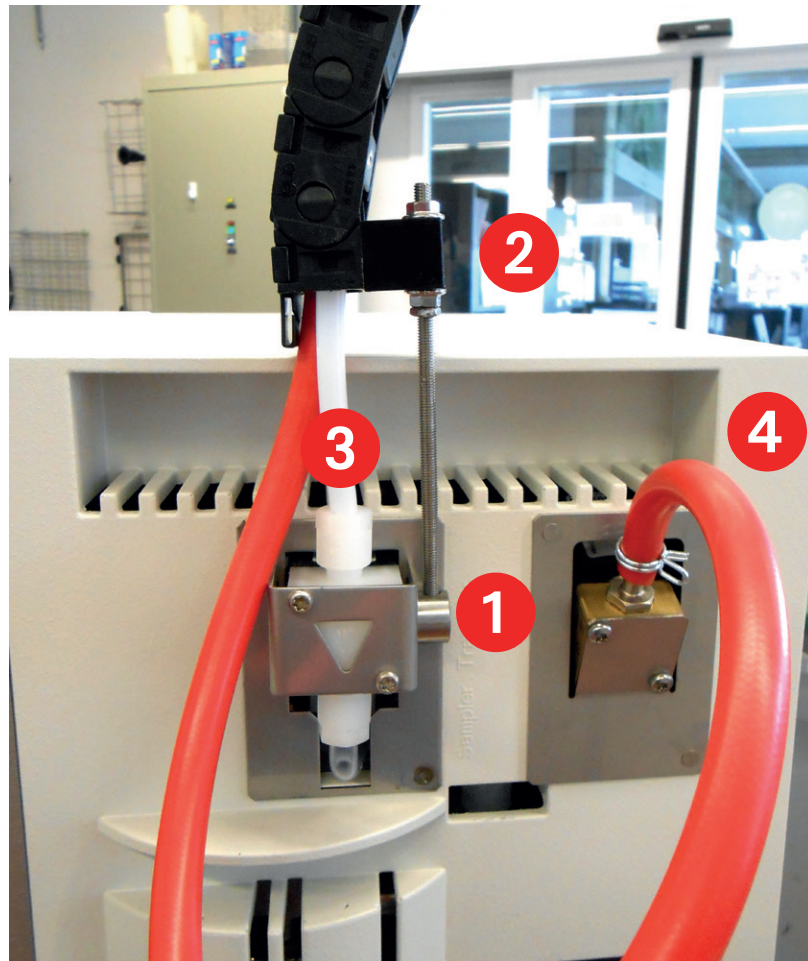

Picture 2: Attaching the transfer connection to the K-375 KjelMaster

- Remove the tightening screw from the holder on the valve **1**.
(The screw is not required for connection to K-377.)
- Slide the ring of the chain fastener onto the holder on the valve **1** and fix it by tightening the threaded bar.
- Slide the the plastic holder of the transfer connection onto the threaded bar **2**. Hold it in place by screwing the second nut hand-tight on top.
- Mount the white transfer hose on the valve using the provided screw connection **3**.
- Mount the red steam hose on the steam valve and secure it with a hose clamp **4**.

BÜCHI Labortechnik AG
9230 Flawil 1/Switzerland
T +41 71 394 63 63
F +41 71 394 65 65

www.buchi.com

Quality in your hands

