


Mounting the level sensor to the secondary condenser

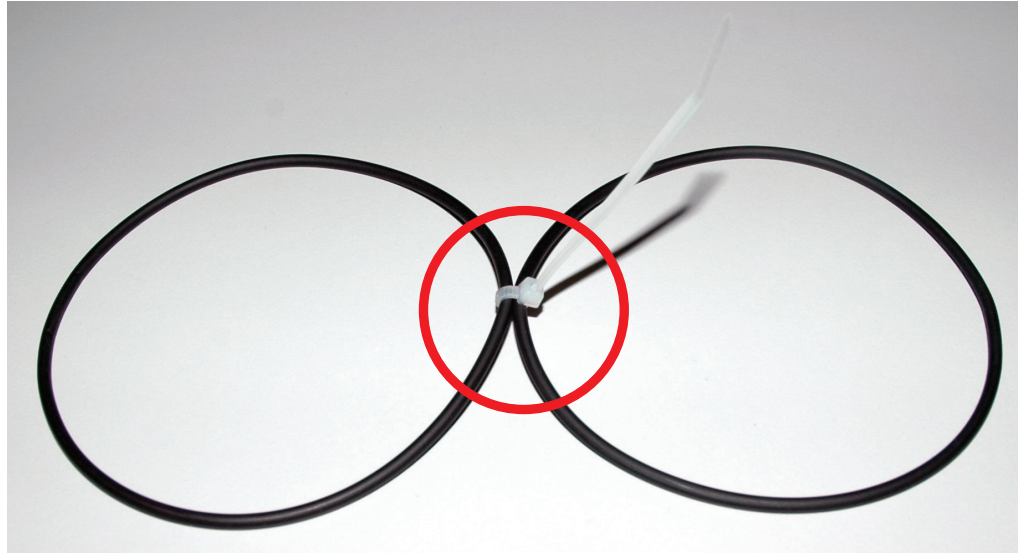

Picture 1: Tying up the two O-rings



Picture 2: Attaching the level sensor to the flask

- Tie the two O-rings up, using the provided cable strap. (see picture 1)
 - Attach the level sensor to the secondary condenser flask using the prepared O-rings. (see picture 2)
- (The cross on the level sensor has to point towards the inside of the flask!)**

BÜCHI Labortechnik AG
9230 Flawil 1/Switzerland
T +41 71 394 63 63
F +41 71 394 65 65

www.buchi.com

Quality in your hands

