



Bath shield (11055364)

Installation instruction

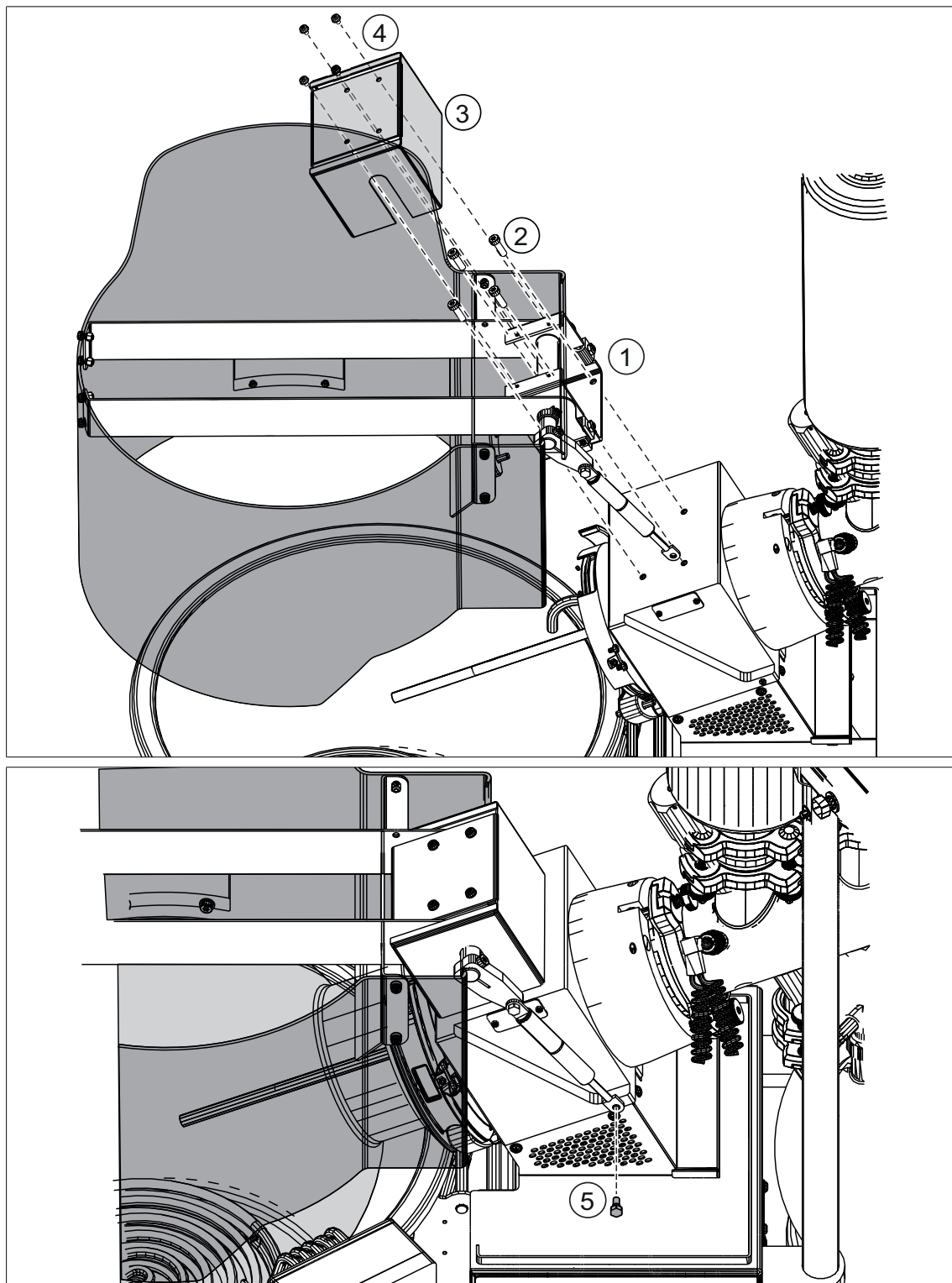

Activate bath shield R-220SE

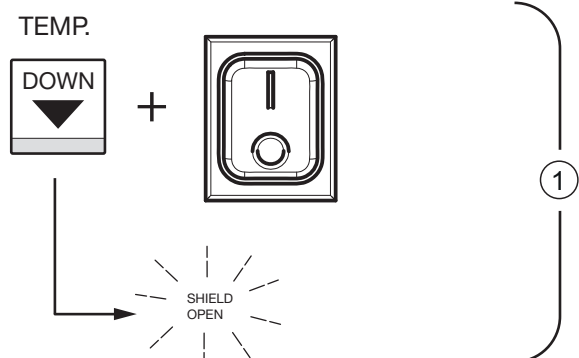


- Keep the **TEMP DOWN** button pressed while switching on the Rotavapor to enter the setup menu.
⇒ **SHIELD OPEN** flashes.




- Press the **TEMP DOWN** button to select **on** (activate bath shield)
- Press **AERATE** button to store the selected setting.

Activate bath shield R-220Pro

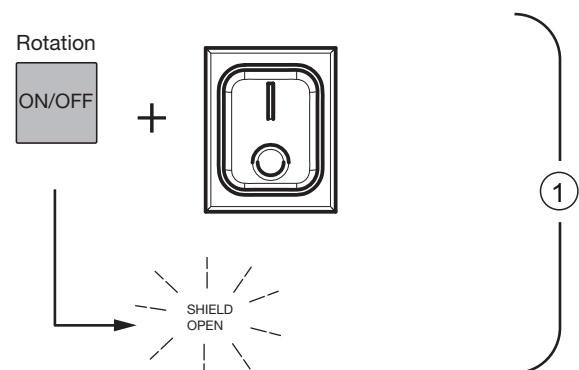

- Keep the button **Rotation ON/OFF** pressed to pressed while switching on the Rotavapor to enter the setup menu.
⇒ **SHIELD OPEN** flashes.

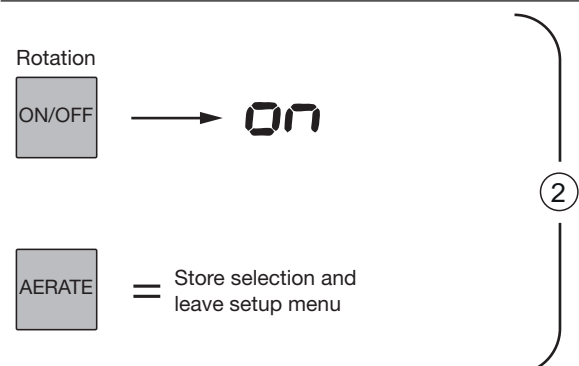

- Press the button **Rotation ON/OFF** to select **on** (activate bath shield)
- Press **AERATE** button to store the selected setting.