

Operation Manual

Interface I-80



Imprint

Product Identification:
Operation Manual (Original) Interface I-80
11594493

Publication date: 03.2024

Version A

BÜCHI Labortechnik AG
Meierseggstrasse 40
Postfach
CH-9230 Flawil 1
E-Mail: quality@buchi.com

BUCHI reserves the right to make changes to the manual as deemed necessary in the light of experience, especially with respect to structure, illustrations and technical details.
This manual is copyrighted. Information from it may neither be reproduced, distributed, or used for competitive purposes, nor made available to third parties. The manufacture of any component with the aid of this manual without prior written agreement is also prohibited.

Table of contents

| | | |
|----------|---|-----------|
| 1 | About this document | 5 |
| 1.1 | Mark-ups and symbols..... | 5 |
| 1.2 | Trademarks..... | 5 |
| 1.3 | Connected instruments..... | 5 |
| 2 | Safety | 6 |
| 2.1 | Intended use..... | 6 |
| 2.2 | Use other than that intended..... | 6 |
| 2.3 | Staff qualification..... | 6 |
| 2.4 | Personal protective equipment..... | 7 |
| 2.5 | Warning notices in this document..... | 7 |
| 2.6 | Warning symbols..... | 7 |
| 2.7 | Residual risks..... | 7 |
| 2.7.1 | Faults during operation..... | 7 |
| 2.8 | Modifications..... | 7 |
| 3 | Product description | 9 |
| 3.1 | Description of function..... | 9 |
| 3.2 | Configuration..... | 9 |
| 3.2.1 | Front view..... | 9 |
| 3.2.2 | Rear view..... | 10 |
| 3.2.3 | Connections..... | 10 |
| 3.2.4 | Display layout..... | 11 |
| 3.2.5 | Display symbols..... | 11 |
| 3.3 | Scope of delivery..... | 11 |
| 3.4 | Type plate..... | 11 |
| 3.5 | Technical data..... | 12 |
| 3.5.1 | Interface I-80..... | 12 |
| 3.5.2 | Ambient conditions..... | 12 |
| 3.5.3 | Materials..... | 13 |
| 3.5.4 | Installation site..... | 13 |
| 4 | Transport and storage | 14 |
| 4.1 | Transport..... | 14 |
| 4.2 | Storage..... | 14 |
| 4.3 | Lifting the instrument..... | 14 |
| 5 | Installation | 15 |
| 5.1 | Installing the interface..... | 15 |
| 5.1.1 | Installing on the Rotavapor®..... | 15 |
| 5.1.2 | Installing on the vacuum pump..... | 16 |
| 5.1.3 | Installing on a laboratory stand..... | 17 |
| 5.2 | Connecting the BUCHI communication cable..... | 17 |
| 5.3 | Connecting the vacuum pump..... | 18 |

| | | |
|-----------|---|-----------|
| 6 | Operation | 19 |
| 6.1 | Main functions..... | 19 |
| 6.1.1 | Start/Stop the vacuum pump | 19 |
| 6.1.2 | Control the vacuum pump speed | 19 |
| 6.1.3 | Stop the vacuum pump | 19 |
| 6.2 | Adjusting the needle valve..... | 20 |
| 6.2.1 | Vacuum regulation | 20 |
| 6.3 | Aerating the system | 21 |
| 6.4 | Settings..... | 21 |
| 6.4.1 | Setting the pump speed | 21 |
| 6.4.2 | Setting to default settings..... | 22 |
| 6.5 | Advanced settings | 22 |
| 7 | Cleaning and servicing | 23 |
| 7.1 | Maintenance work..... | 23 |
| 7.2 | Cleaning the housing | 23 |
| 8 | Help with faults | 24 |
| 8.1 | Troubleshooting | 24 |
| 8.1.1 | Error codes | 24 |
| 8.1.2 | Customer service | 25 |
| 9 | Taking out of service and disposal | 26 |
| 9.1 | Taking out of service..... | 26 |
| 9.2 | Disposal | 26 |
| 9.3 | Returning the instrument | 26 |
| 10 | Appendix | 27 |
| 10.1 | Spare parts and accessories | 27 |
| 10.1.1 | Spare parts | 27 |
| 10.1.2 | Wear parts | 27 |
| 10.1.3 | Accessories..... | 28 |

1 About this document

This operation manual is applicable for all variants of the instrument. Read this operation manual before operating the instrument and follow the instructions to ensure safe and trouble-free operation.

Keep this operation manual for later use and pass it on to any subsequent user or owner.

BÜCHI Labortechnik AG accepts no liability for damage, faults and malfunctions resulting from not following this operation manual.

If you have any questions after reading this operation manual:

- ▶ Contact BÜCHI Labortechnik AG Customer Service.

<https://www.buchi.com/contact>

1.1 Mark-ups and symbols



NOTE

This symbol draws attention to useful and important information.

- ☑ This character draws attention to a requirement that must be met before the instructions below are carried out.
- ▶ This character indicates an instruction that must be carried out by the user.
- ⇒ This character indicates the result of a correctly carried out instruction.

| Mark-up | Explanation |
|---------------------------|--|
| <i>Window</i> | Software Windows are marked-up like this. |
| <i>Tab</i> | Tabs are marked-up like this. |
| <i>Dialog</i> | Dialogs are marked-up like this. |
| <i>[Button]</i> | Buttons are marked-up like this. |
| <i>[Field names]</i> | Field names are marked-up like this. |
| <i>[Menu / Menu item]</i> | Menus or menu items are marked-up like this. |
| Status | Status is marked-up like this. |
| Signal | Signals are marked-up like this. |

1.2 Trademarks

Product names and registered or unregistered trademarks that are used in this document are used only for identification and remain the property of the owner in each case.

1.3 Connected instruments

In addition to this operation manual, follow the instructions and specifications in the documentation for the connected instruments.

2 Safety

2.1 Intended use

The instrument is intended for regulating and indicating vacuum within an operating range of 0 mbar to atmospheric pressure. The instrument has been designed and built as an item of laboratory equipment and can be used in conjunction with the following devices:

- Distillation apparatus, especially rotary evaporators (Rotavapor®)
- Vacuum-drying cabinets
- Vacuum pumps

2.2 Use other than that intended

Use of any kind other than that described in the section Chapter 2.1 "Intended use", page 6 and any application that does not comply with the technical specifications (see Chapter 3.5 "Technical data", page 12) constitutes use other than that intended.

In particular, the following applications are not permissible:

- Use of the instrument in an environment with a potential risk of explosion or areas which require explosion-safe apparatus.
- Use of the instrument with gases with unknown chemical composition.
- Use of the instrument at pressure above atmospheric.
- Use of the instrument for calibration of other equipment.

Damage or hazards attributable to use of the product other than as intended are entirely at the risk of the user alone.

2.3 Staff qualification

Unqualified persons are unable to identify risks and are therefore exposed to greater dangers.

The instrument must be operated by suitably qualified laboratory staff.

The instrument must be commissioned and serviced by suitably qualified technicians.

These operating instructions are aimed at the following target groups:

Users

The users are persons that meet the following criteria:

- They have been instructed in the use of the instrument.
- They are familiar with the contents of these operating instructions and the applicable safety regulations and apply them.
- They are able, on the basis of their training or professional experience, to assess the risks associated with the use of the instrument.

Operator

The operator (generally the laboratory manager and qualified technicians) are responsible for the following aspects:

- The instrument must be correctly installed, commissioned, operated and serviced.
- Only suitably qualified staff must be assigned the task of performing the operations described in these operating instructions.
- The staff must comply with the local applicable requirements and regulations for safe and hazard-conscious working practices.
- Safety-related incidents that occur while using the instrument should be reported to the manufacturer (quality@buchi.com).

BUCHI service technicians

Service technicians authorized by BUCHI have attended special training courses and are authorized by BÜCHI Labortechnik AG to carry out special servicing and repair measures.

2.4 Personal protective equipment

Depending on the application, hazards due to heat and/or corrosive chemicals may arise.

- ▶ Always wear appropriate personal protective equipment such as safety goggles, protective clothing and gloves.
- ▶ Make sure that the personal protective equipment meets the requirements of the safety data sheets for all chemicals used.

2.5 Warning notices in this document

Warning notices warn you of dangers that can occur when handling the instrument. There are four danger levels, each identifiable by the signal word used.

| Signal word | Meaning |
|-------------|---|
| DANGER | Indicates a danger with a high level of risk which could result in death or serious injury if not prevented. |
| WARNING | Indicates a danger with a medium level of risk which could result in death or serious injury if not prevented. |
| CAUTION | Indicates a danger with a low level of risk which could result in minor or medium-severity injury if not prevented. |
| NOTICE | Indicates a danger that could result in damage to property. |

2.6 Warning symbols

The following warning symbols are displayed in this operation manual or on the instrument.

| Symbol | Meaning |
|---|-----------------|
|  | General warning |

2.7 Residual risks

The instrument has been developed and manufactured using the latest technological advances. Nevertheless, risks to persons, property or the environment can arise if the instrument is used incorrectly.

Appropriate warnings in this manual serve to alert the user to these residual dangers.

2.7.1 Faults during operation

If an instrument is damaged, sharp edges, glass splinters, moving parts or exposed electrical wires can cause injuries.

- ▶ Regularly check instruments for visible damage.
- ▶ If faults occur, switch off the instrument immediately, unplug the power cord and inform the operator.
- ▶ Do not continue to use instruments that are damaged.

2.8 Modifications

Unauthorized modifications can affect safety and lead to accidents.

- ▶ Use only genuine BUCHI accessories, spare parts and consumables.
- ▶ Carry out technical changes only with prior written approval from BUCHI.

- ▶ Only allow changes to be made by BUCHI service technicians.

BUCHI accepts no liability for damage, faults and malfunctions resulting from unauthorized modifications.

3 Product description


3.1 Description of function

The instrument is designed for indicating, adjusting and controlling the vacuum.
The instrument can be used to:

- Measure and display the pressure
- Regulate the vacuum pump speed
- Start/Stop the vacuum pump
- Manually regulate the system leakage using a needle valve

3.2 Configuration


3.2.1 Front view

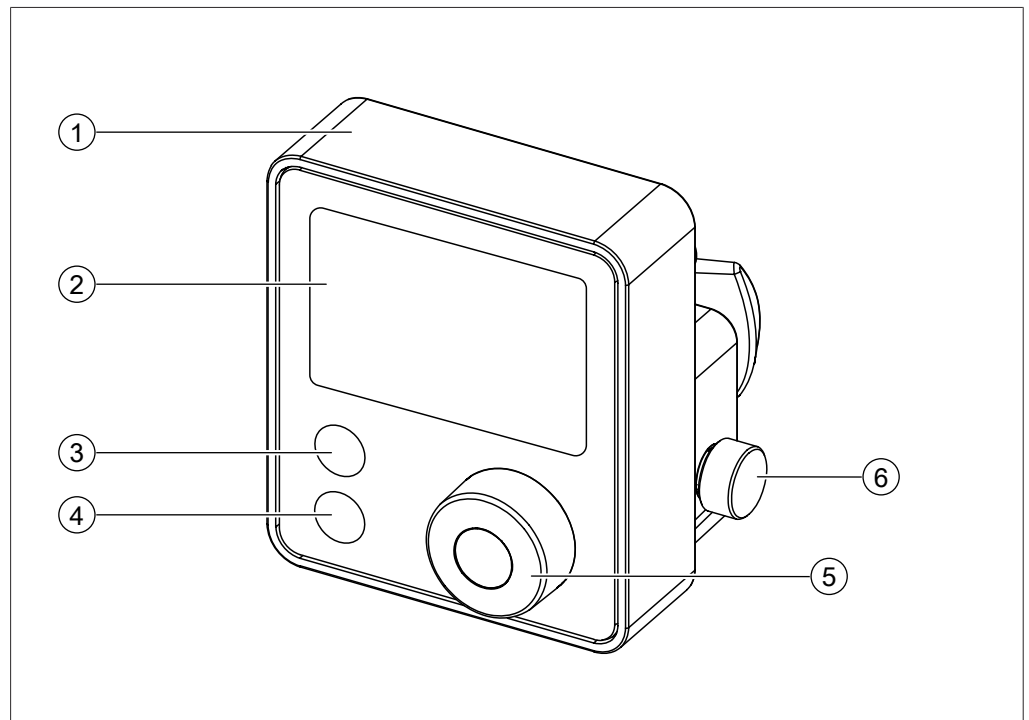

- 1 Interface cover
3 Button **SET**
5 **Navigation control**


- 2 Display
4 Button **STOP**
6 Needle valve control

3.2.2 Rear view

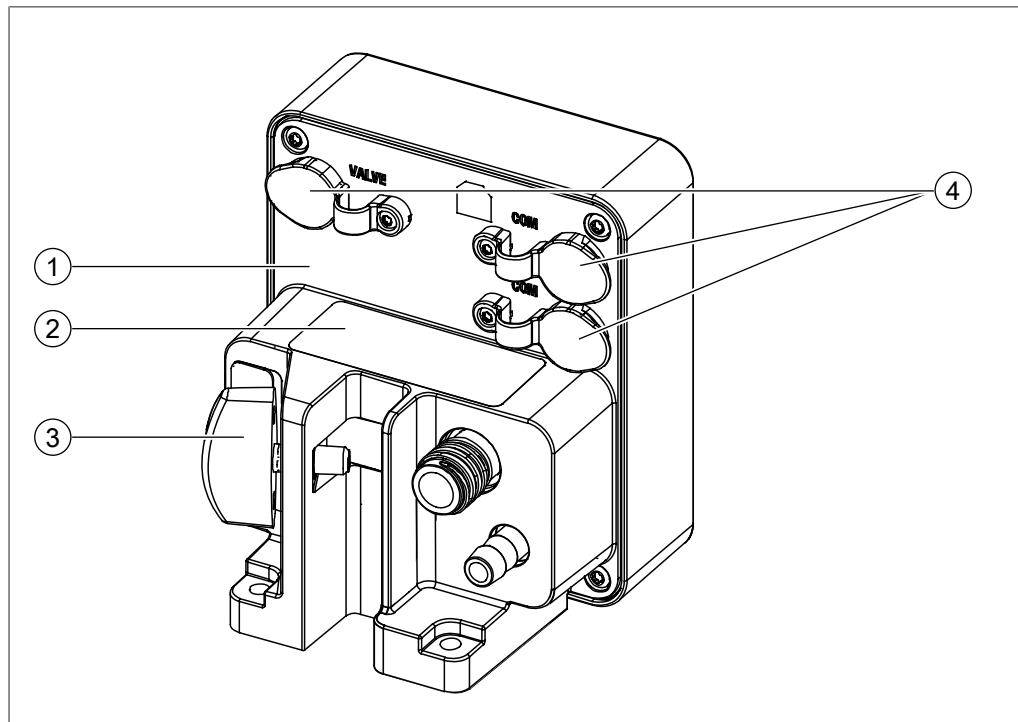


- | | | | |
|---|----------------------|---|------------|
| 1 | Interface back cover | 2 | Type plate |
| 3 | Mounting fixture | 4 | Plug |

3.2.3 Connections

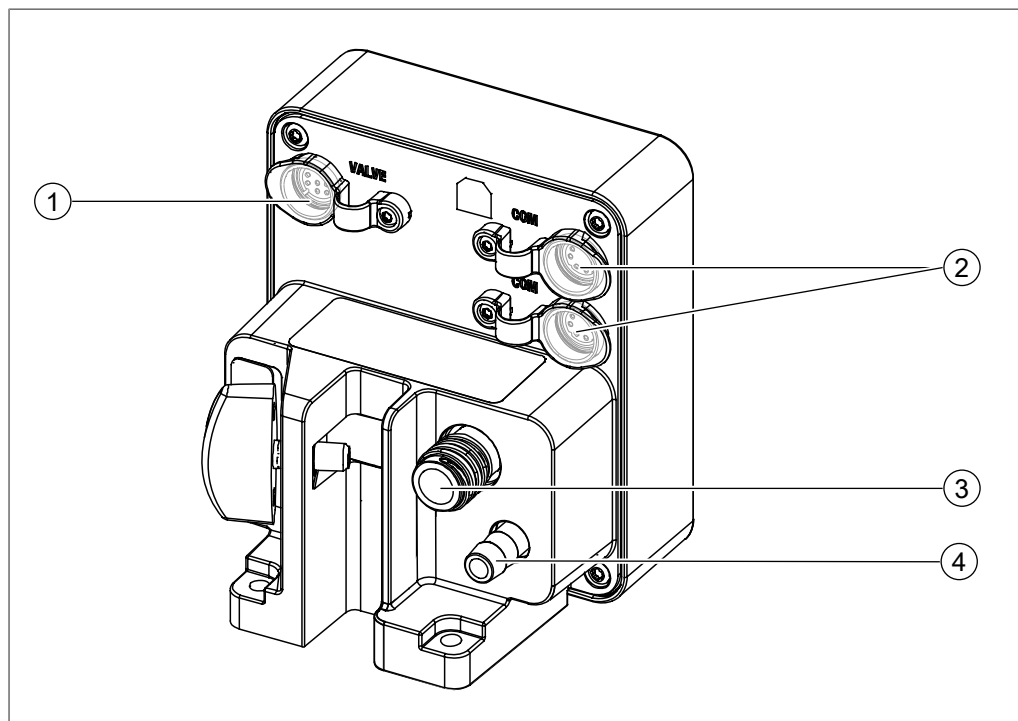

- | | | | |
|---|-------------------------|---|--------------------------|
| 1 | Valve unit VALVE | 2 | Communication COM |
| 3 | Vacuum | 4 | Inert gas |

3.2.4 Display layout

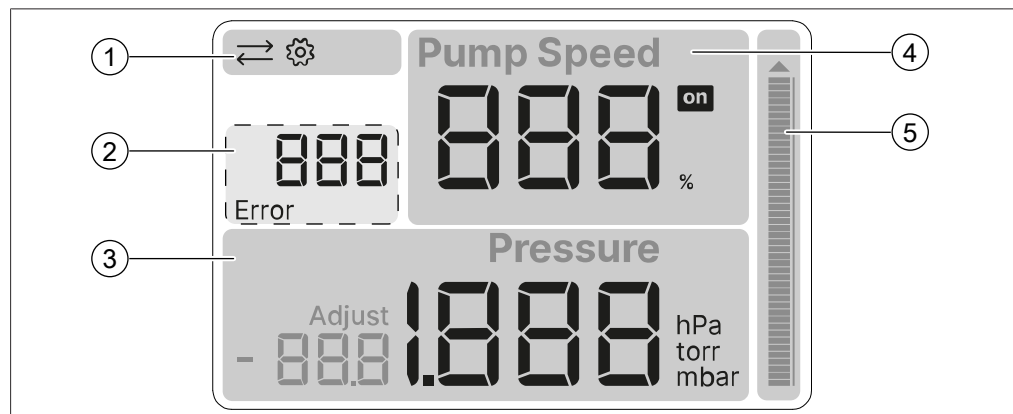

- 1 Status bar
- 2 Error code
- 3 Pressure display
- 4 Pump control
- 5 Pressure indicator

3.2.5 Display symbols

| Symbol | Description |
|--------|-----------------------------|
| | BUCHI COM connected |
| | Settings |
| Error | Error occurrence |
| | Pump ON |
| Adjust | One point calibration value |

3.3 Scope of delivery



NOTE


The scope of delivery depends on the configuration of the purchase order.

Accessories are delivered as per the purchase order, order confirmation, and delivery note.

3.4 Type plate

The type plate identifies the instrument. The following type plate is an example. For more details refer to the type plate on the instrument.

The type plate is located at the rear of the instrument.

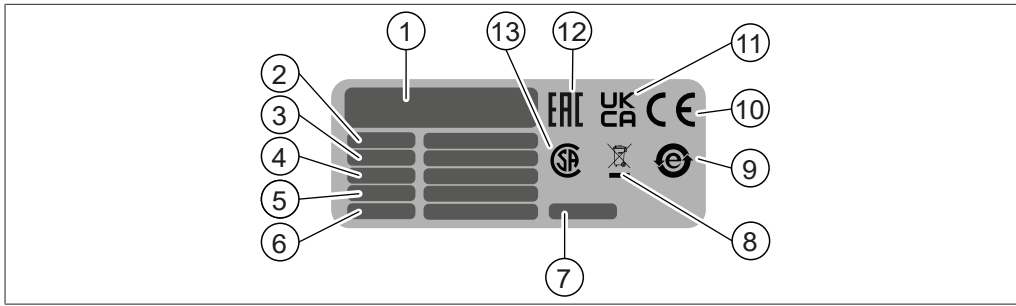

- | | |
|--|---|
| 1 Company name and address | 2 Instrument name |
| 3 Serial number | 4 Input voltage range |
| 5 Frequency | 6 Power consumption maximum |
| 7 Year of manufacture | 8 Symbol for "Do not dispose of as household waste" |
| 9 Symbol for "electronics recycling" | 10 Symbol for "CE conformity" |
| 11 Symbol for "UK Conformity Assessed" | 12 Symbol for "Eurasian Conformity" (optional) |
| 13 Symbol for "CSA certified" (optional) | |

3.5 Technical data

3.5.1 Interface I-80

| Specification | Value |
|--------------------------------|---|
| Dimensions (W × D × H) | 95 mm × 103 mm × 103 mm |
| Weight | 370 g |
| Connection voltage | 30 VDC +5% / -15% |
| Power consumption | 5 W |
| Measurement range | 1,400 – 0 mbar |
| Measurement accuracy | ± 2 mbar (after calibration at constant temperature) |
| Vacuum connection | GL14 |
| IP Code | IP44 |
| Minimum clearance on all sides | none |
| Display type | 3" Dark segment display |
| Certificate | CB, CE, UL/CSA |

3.5.2 Ambient conditions

For indoor use only.

| Specification | Value |
|---------------------------------|---|
| Max. altitude above sea level | 2,000 m |
| Ambient and storage temperature | 5 – 40 °C |
| Max. relative humidity | 80% for temperatures up to 31 °C decreasing linearly to 50% relative humidity at 40 °C |

3.5.3 Materials

| Component | Material |
|----------------------|------------------------------------|
| Casing | PBT |
| Vent tube connection | PP |
| Pressure sensor | Al ₂ O ₃ 96% |

3.5.4 Installation site

- The installation site meets the safety requirements. See Chapter 2 "Safety", page 6
- The installation site meets the specifications according to the technical data (e.g. weight, dimension, etc.). See Chapter 3.5 "Technical data", page 12
- The installation site has no obstacles (e.g. water taps, drains, etc.).
- The installation site is not exposed to external thermal loads, such as direct solar radiation.
- The installation site has enough space that cables / tubes can be routed safely.
- The installation site meets the requirements for the connected devices. See related documentation.
- The installation site is in a well ventilated area.
- The instrument is required to be installed on a Rotavapor®, vacuum pump or a laboratory stand.

4 Transport and storage

4.1 Transport



NOTICE

Risk of breakage due to incorrect transportation

- ▶ Make sure that the instrument is fully dismantled.
- ▶ Pack all instrument components properly to prevent breakage. Use the original packaging whenever possible.
- ▶ Avoid sharp movements during transit.


-
- ▶ After transporting, check the instrument and all glass components for damage.
 - ▶ Damage that has occurred in transit should be reported to the carrier.
 - ▶ Keep packaging for future transportation.

4.2 Storage

- ▶ Make sure that the ambient conditions are complied with (see Chapter 3.5 "Technical data", page 12).
- ▶ Wherever possible, store the instrument in its original packaging.
- ▶ After storage, check the instrument, all seals and tubing for damage and replace if necessary.

4.3 Lifting the instrument

- ▶ Lift the instrument at the points indicated.

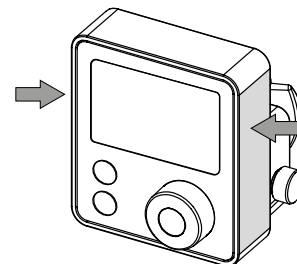

5 Installation

5.1 Installing the interface


The interface is designed to be mounted on BUCHI instruments or a laboratory stand.

5.1.1 Installing on the Rotavapor®


**NOTE**

The interface is designed to be mounted on a Rotavapor® R-80.


- ▶ Remove and discard the two screws on the Rotavapor® R-80 tower.

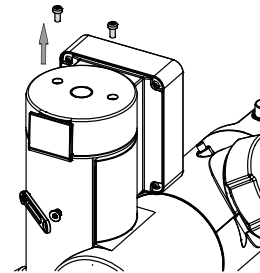


- ▶ Remove the mounting fixture.

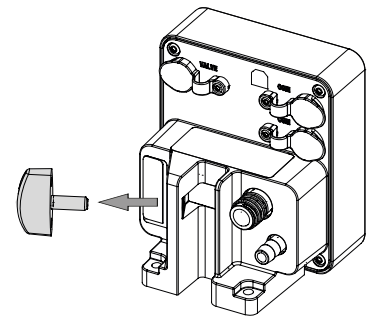


- ▶ Position the interface on the Rotavapor® R-80 tower.

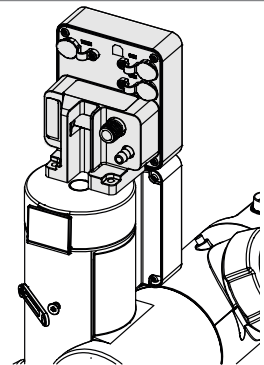

- ▶ Mount the interface with the long screws from the screw set shipped with the interface.
- ▶ Tighten the screws hand-tight. (Max. **0.8 Nm**)

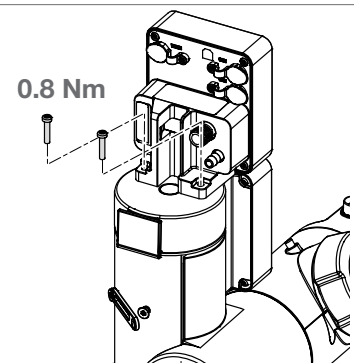

- ▶ Re-install the mounting fixture.

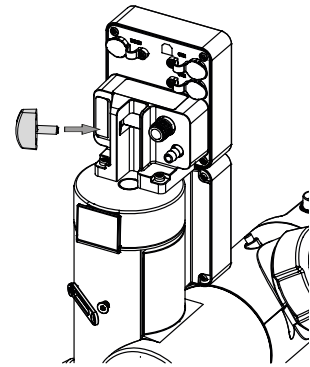

- ▶ Connect the instrument. See additional chapter according to the delivered components.


5.1.2 Installing on the vacuum pump




NOTE

The interface is designed to be mounted on a Vacuum Pump V-80 or Vacuum Pump V-180.

- ▶ Position the interface holder.

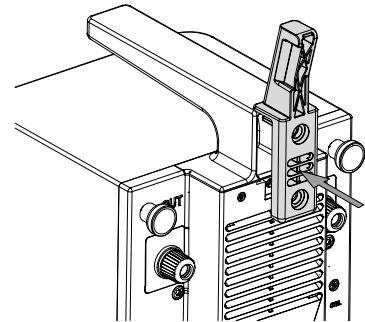


Vacuum Pump V-180

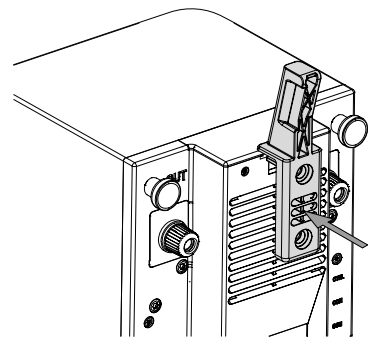


Vacuum Pump V-80


- ▶ Tighten the knurled-head screw.

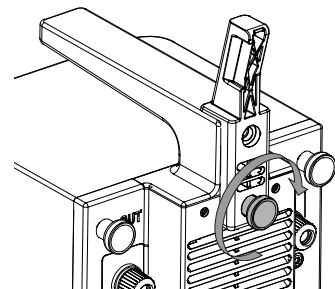

- ▶ Slide the interface onto the interface holder.
- ▶ Position the interface.

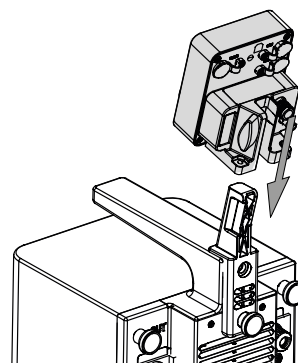


- ▶ Tighten the mounting fixture.

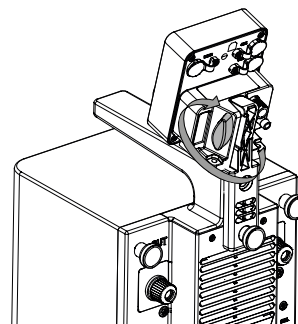

- ▶ Connect the instrument. See additional chapter according to the delivered components.

5.1.3 Installing on a laboratory stand

- ▶ Slide the interface onto the laboratory stand.
- ▶ Position it in place.
- ▶ Tighten the mounting fixture.

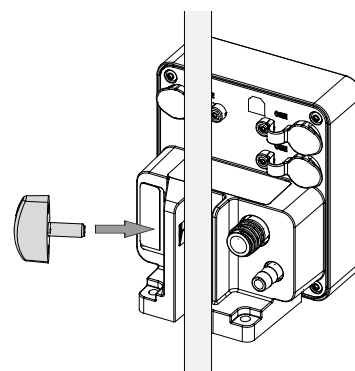


- ▶ Connect the instrument. See additional chapter according to the delivered components.


5.2 Connecting the BUCHI communication cable

Precondition:

- The Interface I-80 / I-180 is installed.
- ▶ Plug the communication cable into the interface.
- ▶ Connect the communication cable to the Rotavapor® or the vacuum pump.

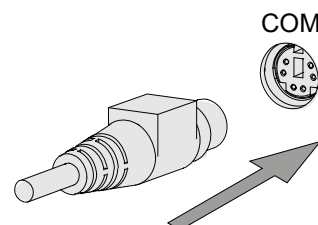

- ▶ Clamp the cable through the fixture.

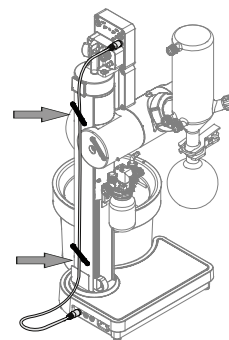

5.3 Connecting the vacuum pump

- ▶ Connect the vacuum pump. See additional manual according to purchase order.




6 Operation

6.1 Main functions


6.1.1 Start/Stop the vacuum pump

- ▶ Push the *navigation control*.
- ⇒ Activates the function.

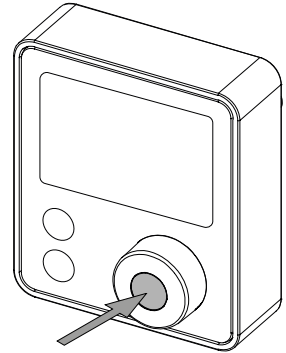


6.1.2 Control the vacuum pump speed

- ▶ Turn the *navigation control*.
- ⇒ Changes the value.

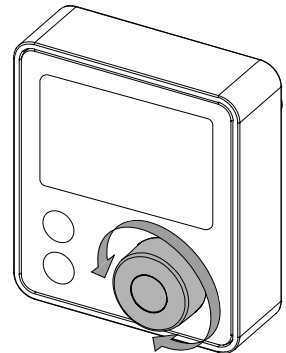


6.1.3 Stop the vacuum pump

- ▶ Touch the **STOP** button.
- ⇒ Stops the instrument.

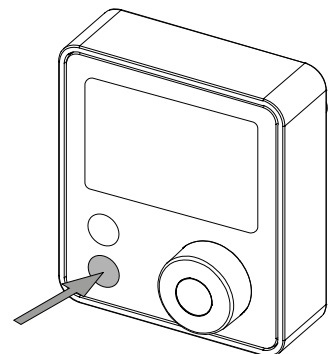

6.2 Adjusting the needle valve

- ▶ Turn the **Needle valve control**.
- ⇒ The needle valve leakage gets regulated.

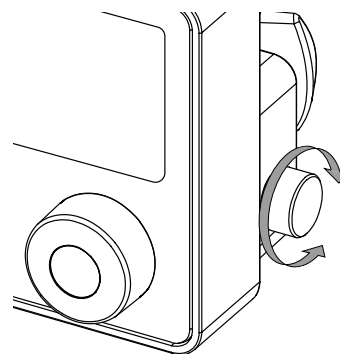

6.2.1 Vacuum regulation

The vacuum regulation (mbar) is achieved with the combination of:

- The pump speed (%)
- The needle valve leakage (rotation)



NOTE

Grey values in the table are not recommended.


Interface I-80 vacuum regulation matrix (mbar) with approximate values with a Vacuum Pump V-80 are indicated. Please consider deviations due to regional differences in atmospheric pressure.

| | | Pump speed | | | | | | | | | |
|----------------------------------|-----|------------|-----|-----|-----|-----|-----|-----|-----|-----|------|
| | | 10% | 20% | 30% | 40% | 50% | 60% | 70% | 80% | 90% | 100% |
| Rotation of needle valve control | 0 | 15 | 15 | 15 | 15 | 15 | 15 | 15 | 15 | 15 | 15 |
| | ¼ | 20 | 20 | 15 | 15 | 15 | 15 | 15 | 15 | 15 | 15 |
| | ½ | 340 | 210 | 160 | 120 | 110 | 100 | 90 | 90 | 80 | 80 |
| | ¾ | 660 | 460 | 350 | 270 | 240 | 210 | 180 | 170 | 170 | 160 |
| | 1 | 810 | 660 | 540 | 440 | 380 | 330 | 300 | 280 | 260 | 250 |
| | 1 ¼ | 860 | 750 | 690 | 580 | 520 | 460 | 390 | 380 | 360 | 340 |
| | 1 ½ | 880 | 800 | 780 | 700 | 640 | 550 | 500 | 490 | 460 | 430 |
| | 1 ¾ | 890 | 820 | 830 | 770 | 720 | 640 | 610 | 600 | 560 | 530 |
| | 2 | 900 | 840 | 890 | 850 | 820 | 790 | 740 | 720 | 700 | 680 |
| | 2 ¼ | 910 | 890 | 900 | 880 | 860 | 840 | 800 | 790 | 770 | 760 |
| | 2 ½ | 930 | 930 | 920 | 890 | 870 | 850 | 830 | 810 | 790 | 780 |
| | 2 ¾ | 930 | 930 | 920 | 900 | 880 | 850 | 840 | 820 | 800 | 790 |
| | 3 | 930 | 930 | 920 | 900 | 880 | 850 | 840 | 820 | 810 | 790 |

6.3 Aerating the system

Precondition:


- ☑ The vacuum pump is not in operation.
- ▶ Turn the aeration cap on the Rotavapor® condenser.
- ⇒ System is aerating.

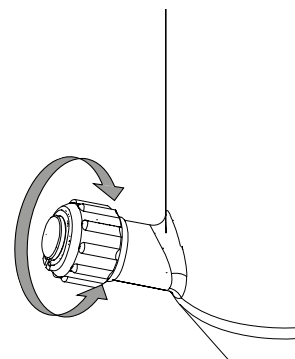


6.4 Settings

6.4.1 Setting the pump speed


Navigation path

→  → Set pump speed


- ▶ Touch the **SET** button.
- ⇒ **Settings** symbol appears.
- ⇒ Blinking value is active.

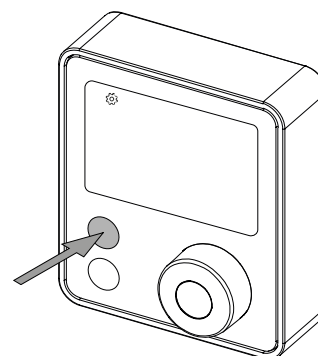

- ▶ Turn the **navigation control**.
- ⇒ Changes the symbol.

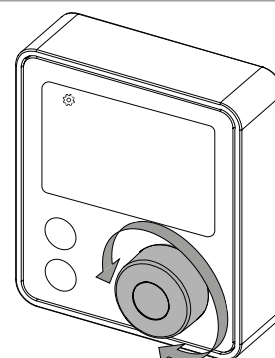


- ▶ Push the **navigation control**.
- ⇒ Exits the settings.

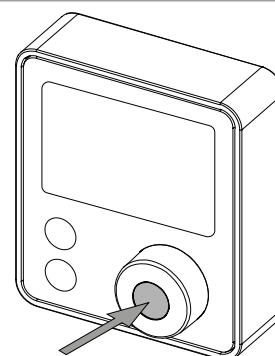

6.4.2 Setting to default settings

Precondition:

- ☑ All connected instruments are switched off.
- ▶ Press and hold the **navigation control**.
- ▶ Switch on a connected instrument.
- ▶ Wait until the indication bar is fully loaded.
- ⇒ Interface is reset to default settings.

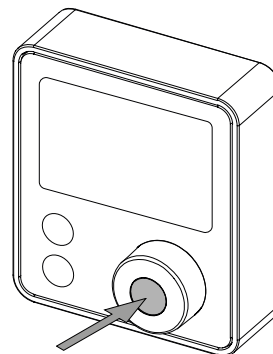


6.5 Advanced settings

| Navigation path | Symbol | Description |
|----------------------|------------------------|---|
| Pressure unit | mbar, torr, hPa | Changing the pressure unit. |
| Pressure calibration | Adjust | Setting an offset for the pressure calibration. |


- ▶ Touch the **SET** button and the **navigation control**.

- ⇒ **Settings** symbol appears.
- ⇒ Blinking symbol or value is active.

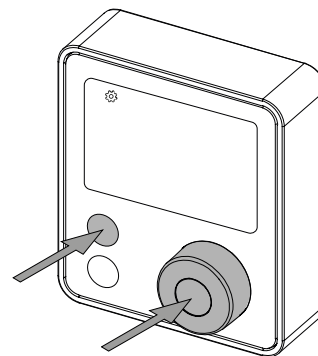


- ▶ Turn the **navigation control**.

- ⇒ Changes the symbol or value.

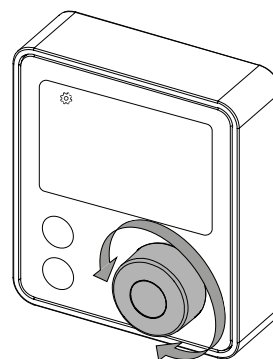

- ▶ Touch the **SET** button to navigate through the settings.

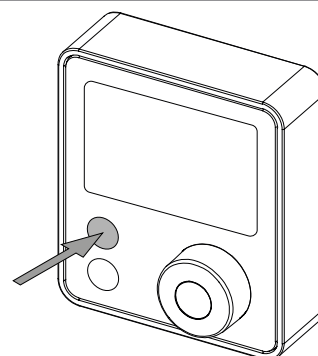

7 Cleaning and servicing



NOTE

- ▶ Carry out only the service and cleaning operations described in this section.
- ▶ Do not carry out any servicing and cleaning operations that involve opening the housing.
- ▶ Use only genuine BUCHI spare parts in order to ensure correct operation and preserve the warranty.
- ▶ Carry out the service and cleaning operations described in this section to extend the lifetime of the instrument.

7.1 Maintenance work

| Action | Weekly | Additional information |
|-----------------------------|---------------|------------------------|
| 7.2 Cleaning the housing | 1 | |
| 1 - Operator | | |

7.2 Cleaning the housing

- ▶ Wipe down the housing with a damp cloth.
- ▶ If heavily soiled, use ethanol or a mild detergent.
- ▶ Wipe down the display with a damp cloth.

8 Help with faults

8.1 Troubleshooting

| Problem | Possible cause | Action |
|---|--|---|
| Instrument does not work | BUCHI communication cable is not connected | ▶ Plug-in the communication cable. See Chapter 5.2 "Connecting the BUCHI communication cable", page 17. |
| | Main switch of the connected instrument with power supply is off | ▶ Switch on the main switch of the connected instrument with power supply. |
| Displayed pressure is wrong | Wrong calibration | ▶ Do the one-point calibration in the advanced setting. |
| Desired vacuum level can not be reached | Needle valve sealing is damaged | ▶ Unscrew the needle valve. |
| | | ▶ Replace the needle valve sealing. |
| | | ▶ Re-install the needle valve. |

8.1.1 Error codes

| Error code | Description | Action |
|------------|--------------------------------|---|
| 140 | Supply voltage too low | ▶ Check the power supply. ⇒ If the error code is still indicated. ▶ Contact BUCHI Customer Service. |
| 150 | Maximum pressure exceeded | ▶ Check possible causes of the system overpressure. |
| 180 | Pressure sensor not calibrated | ▶ Contact BUCHI Customer Service. |
| 181 | Pressure sensor defective | ▶ Contact BUCHI Customer Service. |
| 199 | Initialization error | ▶ Contact BUCHI Customer Service. |
| 391 | Valve unit driver error | ▶ Contact BUCHI Customer Service. |

Error codes from connected BUCHI Vacuum Pump

| Error code | Description | Action |
|------------|--|---|
| 116 | Fill level sensor removed during operation | ▶ Check the sensor connection. |
| 117 | Fill level sensor reached | ▶ Empty the secondary condenser receiving flask. |
| 650 | No fan response | ▶ Make sure the fan can rotate. ⇒ If the error code is still indicated. |
| | | ▶ Contact BUCHI Customer Service. |
| 651 | Housing open | ▶ Make sure the housing is closed. ⇒ If the error code is still indicated. |
| | | ▶ Contact BUCHI Customer Service. |

| Error code | Description | Action |
|-------------------|----------------------------------|---|
| 681 | Electronic circuitry over-heated | <ul style="list-style-type: none"> ▶ Switch off the instrument. ▶ Let the instrument cool down. ▶ Clean the air intake. ▶ Switch on the instrument. ⇒ If the error code is still indicated. ▶ Contact BUCHI Customer Service. |
| 682 | No reliable rotation measurement | ▶ Contact BUCHI Customer Service. |
| 683 | Wrong turning direction | ▶ Contact BUCHI Customer Service. |
| 685 | Motor blocked | ▶ Contact BUCHI Customer Service. |
| 686 | Supply voltage too low | <ul style="list-style-type: none"> ▶ Check the power supply. ⇒ If the error code is still indicated. ▶ Contact BUCHI Customer Service. |

8.1.2 Customer service

Only authorized service personnel are allowed to perform repair work on the instrument which is not described in this manual. Authorization requires a comprehensive technical training and knowledge of possible dangers which might arise when working at the instrument. Such training and knowledge can only be provided by BUCHI.

The customer service and support offers the following support:

- Spare part delivery
- Repairs
- Technical advice

Addresses of official BUCHI customer service offices can be found on the BUCHI website.

www.buchi.com

9 Taking out of service and disposal

9.1 Taking out of service

- ▶ Switch off the instrument.
- ▶ Disconnect the instrument from the power supply.
- ▶ Remove all connections.
- ▶ Clean the instrument.

9.2 Disposal

The operator is responsible for proper disposal of the instrument.

- ▶ When disposing the equipment observe the local regulations and statutory requirements regarding waste disposal.
- ▶ When disposing, observe the disposal regulations of the materials used. For the used materials see Chapter 3.5 "Technical data", page 12 or the material labeling on the parts.

9.3 Returning the instrument

Before returning the instrument, contact the BÜCHI Labortechnik AG Service Department.

<https://www.buchi.com/contact>

10 Appendix

10.1 Spare parts and accessories




Use only genuine BUCHI consumables and spare parts in order to ensure correct, safe and reliable operation of the system.







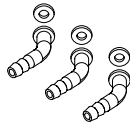
NOTE

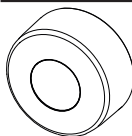Any modifications of spare parts or assemblies are only allowed with the prior written permission of BUCHI.

10.1.1 Spare parts





| | Order no. | Image |
|---|-----------|--|
| Navigation control knob | 11074581 |  |
| Receiving vessel, GL 40, 125 mL, P+G For Woulff bottle | 047233 |  |
| Woulff bottle holder | 11075161 |  |

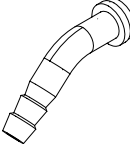
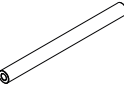
10.1.2 Wear parts

| | Order no. | Image |
|---|-----------|---|
| Seal Woulff bottle | 047165 |  |
| O-ring Needle valve, NBR | 003560 |  |
| Hose barb, bent, GL 14, incl. silicone seal | 018916 |  |
| Hose barbs, set. 2 pcs, bent (1), straight (1), GL 14, silicone seal Content: Hose barbs, cap nuts, seals | 041939 |  |
| Hose barbs, set. 3 pcs, bent, GL 14, silicone seal Content: Hose barbs, seals. | 041987 |  |






| | Order no. | Image |
|---|-----------|---|
| Hose barbs, set. 4 pcs, bent GL 14, silicone seal Content: Hose barbs, cap nuts, seals | 037287 |  |
| Hose barbs, set. 4 pcs, bent, GL 14, EPDM seal Content: Hose barbs, cap nuts, seals. | 043129 |  |
| Hose barbs, set. 4 pcs, bent, GL 14, FPM seal Content: Hose barbs, cap nuts, seals. | 040295 |  |
| Hose barbs, set. 4 pcs, straight, GL 14, EPDM seal Content: Hose barbs, cap nuts, seals. | 043128 |  |
| Hose barbs, set. 4 pcs, straight, GL 14, FPM seal Content: Hose barbs, cap nuts, seals. | 040296 |  |
| Hose barbs, set. 4 pcs, straight, GL 14, silicone seal Content: Hose barbs, cap nuts, seals | 037642 |  |
| Hose barbs, set. 6 pcs, bent (4), straight (2), GL 14, silicone seal Content: Hose barbs, cap nuts, seals. | 038000 |  |

10.1.3 Accessories
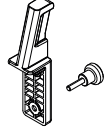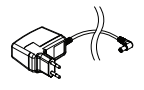
| | Order no. | Image |
|--|-----------|---|
| Stand. V stand with rod, 600 mm Used as a holder for interface when it cannot be mounted on an instrument. | 048891 |  |
| Woulff bottle, 125 mL, P+G, incl. holder For trapping particles and droplets and for pressure equalization. | 11075622 |  |
| Interface holder To install the Interface I-80 / I-180 on the Vacuum pump V-80 / V-180 | 11075588 |  |
| Power adapter. 30 V, 30 W, frequency 50/60 Hz In combination with Mini-DIN to DC-Jack adapter to power the interface. | 11060669 |  |

| | Order no. | Image |
|---|------------------|---|
| Mini-DIN to DC-Jack, 300 mm | 11062266 |  |
| Tubing, synthetic rubber, Ø6/13 mm, black, per m Use: Vacuum | 11063244 |  |

Communication cable
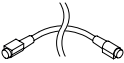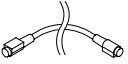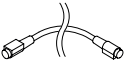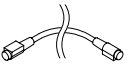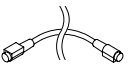
| | Order no. | Image |
|---|------------------|---|
| Communication cable. BUCHI COM, 0.3 m, 6p | 11058705 |  |
| Communication cable. BUCHI COM, 0.9 m, 6p | 11070540 |  |
| Communication cable. BUCHI COM, 1.8 m, 6p | 11058707 |  |
| Communication cable. BUCHI COM, 5.0 m, 6p | 11058708 |  |
| Communication cable. BUCHI COM, 15 m, 6p | 11064090 |  |



11594493 | A en

We are represented by more than 100 distribution partners worldwide.
Find your local representative at:

www.buchi.com

Quality in your hands
