

# Interface I-80 / I-180 on the Rotavapor® R-80

Unlock more insights and guidance through our media resources:



[www.buchi.com/r-80/installation](http://www.buchi.com/r-80/installation)



## NOTE

Reading and following the instructions in this document does not replace reading and following the instructions in the operation manual.

## Tools

	Order no.	Image
Torx size 20	none	


## Installing on the Rotavapor®




## NOTE

The interface is designed to be mounted on a Rotavapor® R-80.


- ▶ Remove and discard the two screws on the Rotavapor® R-80 tower.

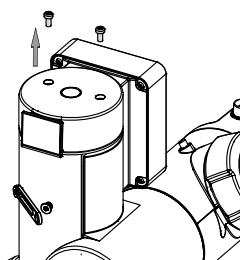


- ▶ Remove the mounting fixture.

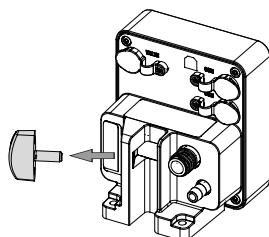


- ▶ Position the interface on the Rotavapor® R-80 tower.

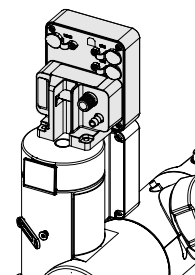

- ▶ Mount the interface with the long screws from the screw set shipped with the interface.
- ▶ Tighten the screws hand-tight. (Max. 0.8 Nm)

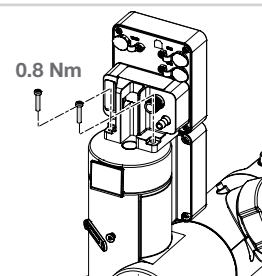

- ▶ Re-install the mounting fixture.

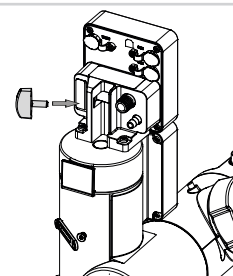

- ▶ Connect the instrument. See additional chapter according to the delivered components.

