

# Installation Guide Scrubber OneScrub

## 1. Scope of Delivery

Product	Article No.
Scrubber OneScrub	114150010

## 2. Installation

1. Place the 3L condensate vessel on the black ground plate on the left side and place the vessel cover on top of it.

2. Mount the absorption vessel on the cover. Make sure it is filled with activated charcoal granules for absorption of fumes.

3. Place the pump Vacuum pump on the black ground plate on the right side and connect it to an electrical plug.

4. Connect the black hose coming from the SpeedDigester to the inlet of the neutralization vessel and the black hose from the outlet of the adsorption vessel to the vacuum pump.

5. Before starting operation during digestion, please make sure that the neutralization vessel is filled with sufficient level of buffer solution.

