

Operation Manual

K-365 Dist Line



Imprint

Product Identification:
Operation Manual (Original) K-365 Dist Line
11594193

Publication date: 04.2026

Version F

BÜCHI Labortechnik AG
Meierseggrasse 40
Postfach
CH-9230 Flawil 1
E-Mail: quality@buchi.com

BUCHI reserves the right to make changes to the manual as deemed necessary in the light of experience, especially with respect to structure, illustrations and technical details.
This manual is copyrighted. Information from it may neither be reproduced, distributed, or used for competitive purposes, nor made available to third parties. The manufacture of any component with the aid of this manual without prior written agreement is also prohibited.

Table of contents

1	About this document	8
1.1	Mark-ups and symbols	8
1.2	Trademarks	8
1.3	Connected devices	8
2	Safety	9
2.1	Proper use	9
2.2	Use other than that intended	9
2.3	Staff qualification	9
2.4	Personal protective equipment	9
2.5	Warning notices in this document	10
2.6	Warning and directive symbols	10
2.7	Protection devices	11
2.8	Residual risks	11
	2.8.1 Hot surfaces	11
	2.8.2 Faults during operation	11
	2.8.3 Glass breakage	12
	2.8.4 Not suitable titrators	12
	2.8.5 Dosing pump defect	12
	2.8.6 Acidic liquids	12
2.9	Modifications	12
2.10	Quality of the repair	13
3	Product description	14
3.1	Description of function	14
3.2	Configuration	14
	3.2.1 Front view	14
	3.2.2 Rear view	16
	3.2.3 Receiving area	17
	3.2.4 Communication connections	18
	3.2.5 Connections on the rear side	20
3.3	Scope of delivery	23
3.4	Type plate	24
3.5	Technical data	24
	3.5.1 K-365 Dist Line	24
	3.5.2 Ambient conditions	27
	3.5.3 Materials	27
	3.5.4 Installation site	27
4	Transport and storage	28
4.1	Transport	28
4.2	Storage	28
4.3	Lifting the instrument	28

5	Installation	29
5.1	Before installation.....	29
5.2	Establishing electrical connections	29
5.3	Securing against earthquakes.....	29
5.4	Installing the cooling water supply	30
5.4.1	Installing a tap water supply (option)	30
5.4.2	Installing the recirculating chiller on the condenser (option)	30
5.5	Installing the drain tubing (option).....	31
5.6	Installing the waste pump connection (MultiDist option only).....	31
5.7	Installing the sample waste pump connection (BasicDist option, MultiDist only)	32
5.8	Installing the H ₂ O supply for steam generation.....	32
5.9	Installing NaOH supply (MultiDist only).....	33
5.10	Installing H ₂ BO ₃ supply (MultiDist only).....	33
5.11	Installing the H ₂ O supply for the sample tube (MultiDist only)	34
5.12	Installing the level sensor on the canister (option).....	35
5.13	Installing a titrator with LAN (MultiDist option with Eco Titrator only)	35
5.14	Installing a titrator with RS232 (MultiDist option only).....	36
5.15	Installing a dispenser (MultiDist only).....	37
5.16	Installing the reaction detection sensor (MultiDist option only)	37
5.17	Installing a SD card (Interface pro only).....	37
5.18	Installing the OnLevel sensor (option).....	37
5.19	Installing the stirrer (MultiDist option only)	38
6	Description of the interface (EasyDist and BasicDist)	39
6.1	Layout of the interface.....	39
6.2	Function bar interface	40
6.3	Menu bar interface	40
6.3.1	Home menu	41
6.3.2	Manual control menu	41
6.3.3	Method menu	41
6.3.4	Configuration menu.....	41
6.3.5	Determination data menu.....	43
6.4	Status bar interface	43
7	Description of the interface pro (BasicDist and MultiDist)	44
7.1	Layout of the interface pro	44
7.2	Function bar interface pro	45
7.3	Menu bar interface pro	45
7.3.1	Home menu	46
7.3.2	Manual control menu	46
7.3.3	Method menu	47
7.3.4	Configuration menu.....	47
7.3.5	Determination data menu.....	49
7.4	Status bar interface pro	49

8	Preparations for a determination.....	50
8.1	Editing the preheating function.....	50
8.2	Editing the priming function.....	50
8.3	Editing the single determination.....	50
8.3.1	Changing the name of the single determination.....	50
8.3.2	Selecting the determination type for a single determination.....	51
8.3.3	Selecting the method for a single determination.....	51
8.3.4	Changing the sample quantity for single determination.....	52
8.3.5	Changing the unit for a single determination.....	52
8.4	Editing a series.....	53
8.4.1	Creating a new series.....	53
8.4.2	Changing the name of a series.....	53
8.4.3	Adding a determination to a series.....	54
8.4.4	Deleting a determination from a series.....	54
8.4.5	Changing the determination name for a series.....	54
8.4.6	Selecting the determination type within a series.....	55
8.4.7	Selecting the method for a determination within a series.....	55
8.4.8	Changing the sample quantity for determination within a series.....	56
8.4.9	Changing the amount unit for a determination within a series.....	56
8.5	Editing a template.....	57
8.5.1	Creating a new template.....	57
8.5.2	Changing the name of a template.....	58
8.6	Editing a method.....	58
8.6.1	Creating a new method.....	58
8.6.2	Deleting a method.....	58
8.6.3	Changing the name of a method.....	59
8.6.4	Changing the reaction detection for a method (option for MultiDist only).....	59
8.6.5	Changing the H ₂ O volume for a method (MultiDist only).....	59
8.6.6	Changing the H ₂ SO ₄ volume for a method (MultiDist only).....	60
8.6.7	Changing the NaOH volume for a method (MultiDist only).....	60
8.6.8	Changing the acid volume for a method (BasicDist, MultiDist only).....	61
8.6.9	Changing the reaction time for a method (BasicDist, MultiDist only).....	61
8.6.10	Changing the number of steam steps for a method.....	62
8.6.11	Changing the time a steam step is carried out for a method.....	63
8.6.12	Changing the steam power for a method.....	63
8.6.13	Changing the level detection sensor settings (option) for a method.....	64
8.6.14	Changing the distillation time for a method.....	64
8.6.15	Changing the distillation stirring speed for a method (MultiDist only).....	65
8.6.16	Changing the titration type for a method (MultiDist option only).....	65
8.6.17	Changing the H ₃ BO ₃ volume for a method (MultiDist only).....	65
8.6.18	Changing the dosing unit volume for a method (MultiDist option only).....	66
8.6.19	Changing the titration stirring speed for a method (MultiDist option only).....	66
8.6.20	Changing the titration start time for a method (MultiDist option only).....	67
8.6.21	Changing the aspiration time for the sample tube for a method (BasicDist, MultiDist only).....	67
8.6.22	Changing the aspiration time for the receiving vessel for a method (MultiDist only).....	68
8.7	Editing user setting.....	68
8.7.1	Creating a new user.....	68
8.7.2	Deleting a user.....	69
8.7.3	Changing the name of a user.....	69
8.7.4	Selecting a user.....	69
8.8	Installing the sample tube.....	70
8.9	Installing a receiving vessel.....	70
8.10	Preparing the end stop for the sample tube size.....	71
8.11	Preparing the OnLevel sensor (option).....	71
8.12	Selecting the operation mode.....	71

8.13	Dosing manually.....	72
8.13.1	Dosing H ₂ O manually	72
8.13.2	Dosing NaOH manually	72
8.13.3	Dosing H ₃ BO ₃ manually	72
8.14	Aspirating manually.....	72
8.14.1	Aspirating the sample tube manually	72
8.14.2	Aspirating the receiving vessel manually	73
8.15	Changing the stirring speed manually	73
8.16	Switching valve adjustment (MultiDist only).....	73
9	Carrying out a determination	75
9.1	Preparing the instrument.....	75
9.2	Starting a determination	75
9.3	Ending a determination	75
9.4	Shutting down the instrument.....	76
9.5	Filtering determination data.....	76
9.6	Showing determination data.....	76
9.7	Deleting determination data	76
9.7.1	Deleting one determination	76
9.7.2	Deleting all determination data	76
9.8	Exporting determination data (Interface pro only).....	77
9.8.1	Exporting one determination (Interface pro only).....	77
9.8.2	Exporting all determination data (Interface pro only).....	77
10	Cleaning and servicing	78
10.1	Regular maintenance work	78
10.2	Cleaning and servicing the sample tube	79
10.3	Cleaning the glass components	79
10.4	Cleaning and servicing the hoses and hose connections	80
10.5	Cleaning and servicing the sample tube seal.....	80
10.6	Cleaning and servicing the splash protector	80
10.7	Cleaning and servicing the steam inlet and condenser outlet tube	80
10.8	Cleaning the housing	80
10.9	Cleaning and servicing the warning and directive symbols.....	80
10.10	Cleaning and servicing the bridge splash protector to condenser	80
10.11	Cleaning and servicing the dosing pumps	81
10.12	Cleaning and servicing the condenser	81
10.13	Cleaning and servicing the steam generator.....	81
10.14	Calibrating the pumps	81
10.14.1	Calibrating the H ₂ O pump.....	81
10.14.2	Calibrating the acid pump	82
10.14.3	Calibrating the NaOH pump.....	82
10.14.4	Calibrating the H ₃ BO ₃ Pump	83
10.15	Cleaning and servicing the waste pumps.....	83
10.16	Decalcifying the instrument	83
10.17	Replacing the splash protector.....	84
10.18	Replacing the condenser	85
10.19	Attaching and removing the protection shield	86
10.19.1	Removing the protection shield.....	86
10.19.2	Attaching the protection shield.....	86
10.20	Cleaning the drip tray	86
10.21	Rinsing a pump	86
10.22	Replacing the sample tube seal	87

11	Help with faults	88
11.1	Troubleshooting digestion	88
11.2	Troubleshooting distillation.....	89
11.3	Troubleshooting instrument.....	91
11.4	Tighten the sample tube sealing	92
12	Taking out of service and disposal	93
12.1	Taking out of service	93
12.2	Disposal	93
12.3	Returning the instrument.....	93
12.4	Removing water from the steam generator	93
13	Appendix	94
13.1	Spare parts and accessories.....	94
13.1.1	Accessories.....	94
13.1.2	Splash protector spare parts.....	95
13.1.3	Tank spare parts	97
13.1.4	Condenser spare parts	98
13.1.5	Titration related parts	98
13.1.6	Sample tubes	100
13.1.7	Cable and tubing.....	101
13.1.8	Other spare parts	102
13.1.9	Consumables	103
13.1.10	Maintenance kit.....	104
13.1.11	Upgrade kits	104
13.1.12	Documentation.....	104

1 About this document

This operation manual is applicable for all variants of the instrument.

Read this operation manual before operating the instrument and follow the instructions to ensure safe and trouble-free operation.

Keep this operation manual for later use and pass it on to any subsequent user or owner.

BÜCHI Labortechnik AG accepts no liability for damage, faults and malfunctions resulting from not following this operation manual.

If you have any questions after reading this operation manual:

- ▶ Contact BÜCHI Labortechnik AG Customer Service.

<https://www.buchi.com/contact>

1.1 Mark-ups and symbols



NOTE

This symbol draws attention to useful and important information.

- ☑ This character draws attention to a requirement that must be met before the instructions below are carried out.
- ▶ This character indicates an instruction that must be carried out by the user.
- ⇒ This character indicates the result of a correctly carried out instruction.

Mark-up	Explanation
<i>Window</i>	Software Windows are marked-up like this.
<i>Tab</i>	Tabs are marked-up like this.
<i>Dialog</i>	Dialogs are marked-up like this.
<i>[Button]</i>	Buttons are marked-up like this.
<i>[Field names]</i>	Field names are marked-up like this.
<i>[Menu / Menu item]</i>	Menus or menu items are marked-up like this.
Status	Status is marked-up like this.
Signal	Signals are marked-up like this.

1.2 Trademarks

Product names and registered or unregistered trademarks that are used in this document are used only for identification and remain the property of the owner in each case.

1.3 Connected devices

In addition to these operating instructions, follow the instructions and specifications in the documentation for the connected devices.

2 Safety

2.1 Proper use

The instrument is designed for steam distillation.

The instrument can be used in laboratories for the following tasks:

- Distillation of steam-volatile substances.

2.2 Use other than that intended

The use of the instrument other than described in proper use and specified in technical data is use other than that intended.

The operator is responsible for damages or hazards that are caused by use other than that intended.

Specially the following uses are not permitted:

- Use of the instrument in areas which require explosion-safe instruments.
- Use of samples, which can explode or inflame (example: explosives, etc.) due to shock, friction, heat or spark formation.
- Use of the instrument with other than original BUCHI glassware.

2.3 Staff qualification

Unqualified persons are unable to identify risks and are therefore exposed to greater dangers.

The device may only be operated by suitably qualified laboratory staff.

These operating instructions are aimed at the following target groups:

Users

Users are persons that meet the following criteria:

- They have been instructed in the use of the device.
- They are familiar with the contents of these operating instructions and the applicable safety regulations and apply them.
- They are able on the basis of their training or professional experience to assess the risks associated with the use of the device.

Operator

The operator (generally the laboratory manager) is responsible for the following aspects:

- The device must be correctly installed, commissioned, operated and serviced.
- Only suitably qualified staff may be assigned the task of performing the operations described in these operating instructions.
- The staff must comply with the local applicable requirements and regulations for safe and hazard-conscious working practices.
- Safety-related incidents that occur while using the device should be reported to the manufacturer (quality@buchi.com).

BUCHI service technicians

Service technicians authorized by BUCHI have attended special training courses and are authorized by BÜCHI Labortechnik AG to carry out special servicing and repair measures.

2.4 Personal protective equipment

Depending on the application, hazards due to heat and/or corrosive chemicals may arise.

- ▶ Always wear appropriate personal protective equipment such as safety goggles, protective clothing and gloves.
- ▶ Make sure that the personal protective equipment meets the requirements of the safety data sheets for all chemicals used.








2.5 Warning notices in this document

Warning notices warn you of dangers that can occur when handling the device. There are four danger levels, each identifiable by the signal word used.

Signal word	Meaning
DANGER	Indicates a danger with a high level of risk which could result in death or serious injury if not prevented.
WARNING	Indicates a danger with a medium level of risk which could result in death or serious injury if not prevented.
CAUTION	Indicates a danger with a low level of risk which could result in minor or medium-severity injury if not prevented.
NOTICE	Indicates a danger that could result in damage to property.

2.6 Warning and directive symbols

The following warning and directive symbols are displayed in this operation manual or on the instrument.

Symbol	Meaning
	Hot surface
	Corrosive
	General warning
	Device damage
	Dangerous electrical voltage
	Wear protective gloves
	Wear safety goggles

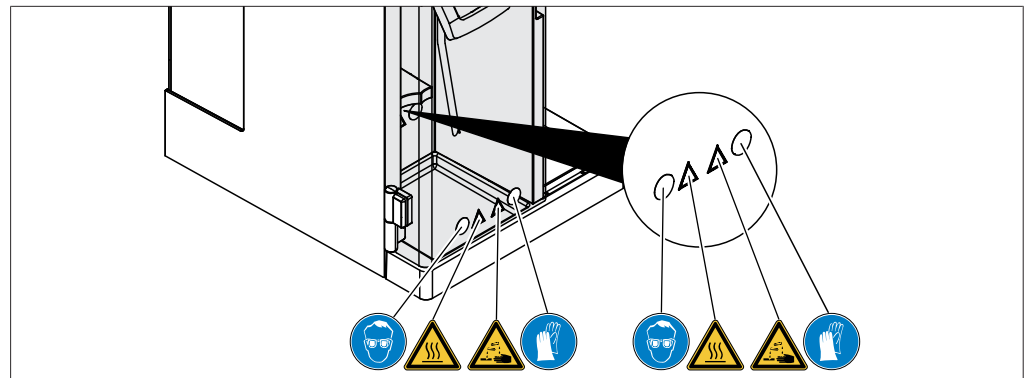


Fig. 1: Location of the warning and directive symbols

2.7 Protection devices

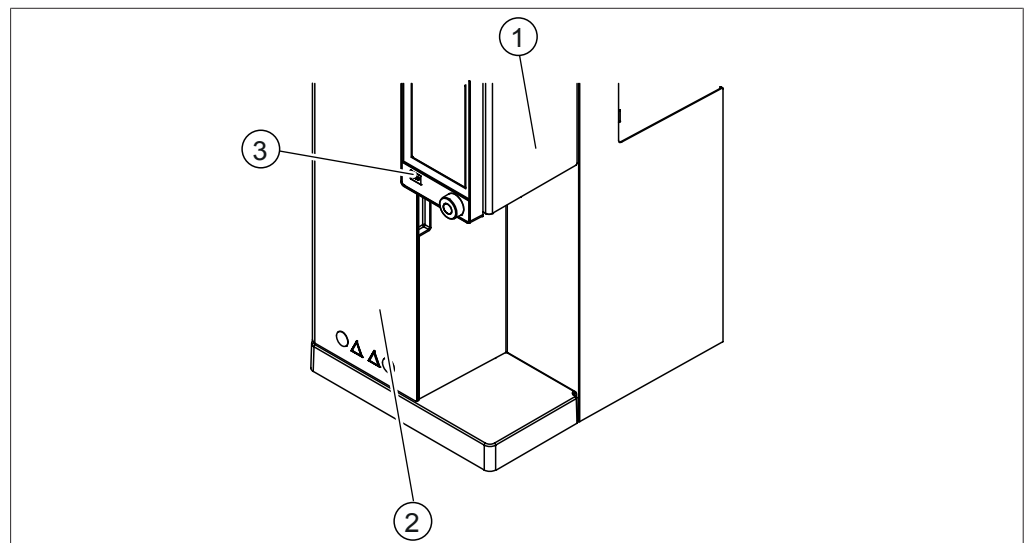


Fig. 2: Protection devices

- | | | | |
|---|-------------------|---|-------------------|
| 1 | Protection shield | 2 | Protection shield |
| 3 | Stop button | | |

2.8 Residual risks

The device has been developed and manufactured using the latest technological advances. Nevertheless, risks to persons, property or the environment can arise if the device is used incorrectly.

Appropriate warnings in this manual serve to alert the user to these residual dangers.

2.8.1 Hot surfaces

The surfaces of the device can become very hot. If touched they can cause skin burns.

- ▶ Do not touch hot surfaces or else wear suitable protective gloves.

2.8.2 Faults during operation

If a device is damaged, sharp edges, moving parts or exposed electrical wires can cause injuries.

- ▶ Regularly check device for visible damage.
- ▶ If faults occur, switch off the device immediately, unplug the power cord and inform the operator.
- ▶ Do not continue to use devices that are damaged.

2.8.3 Glass breakage

Broken glass can cause severe cuts.

Minor damage to the ground joints impairs the sealing effect and may therefore diminish suction capacity.

- Handle the glass components carefully and do not drop them.
- Always place the glassware in a suitable holder when they are not in use.
- Always visually inspect glass components for damage every time they are to be used.
- Do not continue to use glass components that are damaged.
- Always wear protective gloves when disposing of broken glass.

2.8.4 Not suitable titrators

Not suitable titrators can lead to malfunction and wrong results.

- ▶ Use recommended titrators only.

2.8.5 Dosing pump defect

A defect dosing pump can spill hazardous liquids on the installation site.

- ▶ Set the On / Off master switch to Off.
- ▶ Disconnect the power supply.
- ▶ Use personal protective equipment to remove the liquid.
- ▶ Dispose of the remains of the liquid according to the local legal regulations.

2.8.6 Acidic liquids

Accumulation of liquid on surfaces of unknown origin might contain acid.

- ▶ Use personal protective equipment to remove the liquid.
- ▶ Dispose of the remains of the liquid according to the local legal regulations.

2.9 Modifications

Unauthorized modifications can effect safety and lead to accidents.

- ▶ Use only genuine BUCHI accessories, spare parts and consumables.
- ▶ Carry out technical changes only with prior written approval from BUCHI.
- ▶ Only allow changes to be made by BUCHI service technicians.

BUCHI accepts no liability for damage, faults and malfunctions resulting from unauthorized modifications.

2.10 Quality of the repair

Quality of products and service is the basis of a good relationship between the customer and manufacturer for BUCHI. As a help to maintain a good quality service level, some general rules must be followed:


- Before starting any work, make sure the device and parts are decontaminated and clean.
- Document the functional status / problem and system condition upon arrival.
- Discharge yourself from electrostatic charges using the Electro Static Discharge (ESD) service kit (included in the tool kits available from BUCHI upon request) before touching any electronic components.
- If a replacement Printed Circuit Board (PCB) gets damaged immediately after a successful replacement, check for short circuits and overvoltage before exchanging the part again.
- Handle all parts with care. In particular don't bend or twist items to avoid cracks and mechanical stress to components on the Printed Circuit Boards (PCB's).
- Carry out all necessary calibrations and functional tests after replacement of components, Printed Circuit Boards (PCB's) or subassemblies.
- Check for electrical safety after service.
- Inform the customer about the functional status and system condition after service.

3 Product description

3.1 Description of function

The instrument is suitable for determining steam-volatile substances (e.g. of alcohol, sulfur dioxide, volatile acids) and nitrogen by using Kjeldahl and Devarda methods (MultiDist).

- Steam is introduced into the sample solution to drive out volatile components.
- After condensation in condenser the condensate is collected in a receiver solution.

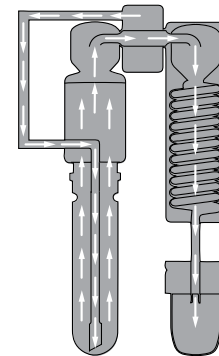

3.2 Configuration

3.2.1 Front view



NOTE

The interface differs according to the configuration:

- ▶ Interface see Chapter 6 “Description of the interface (EasyDist and BasicDist)”, page 39
⇒ EasyDist, BasicDist
- ▶ Interface Pro see Chapter 7 “Description of the interface pro (BasicDist and MultiDist)”, page 44
⇒ BasicDist (option), MultiDist

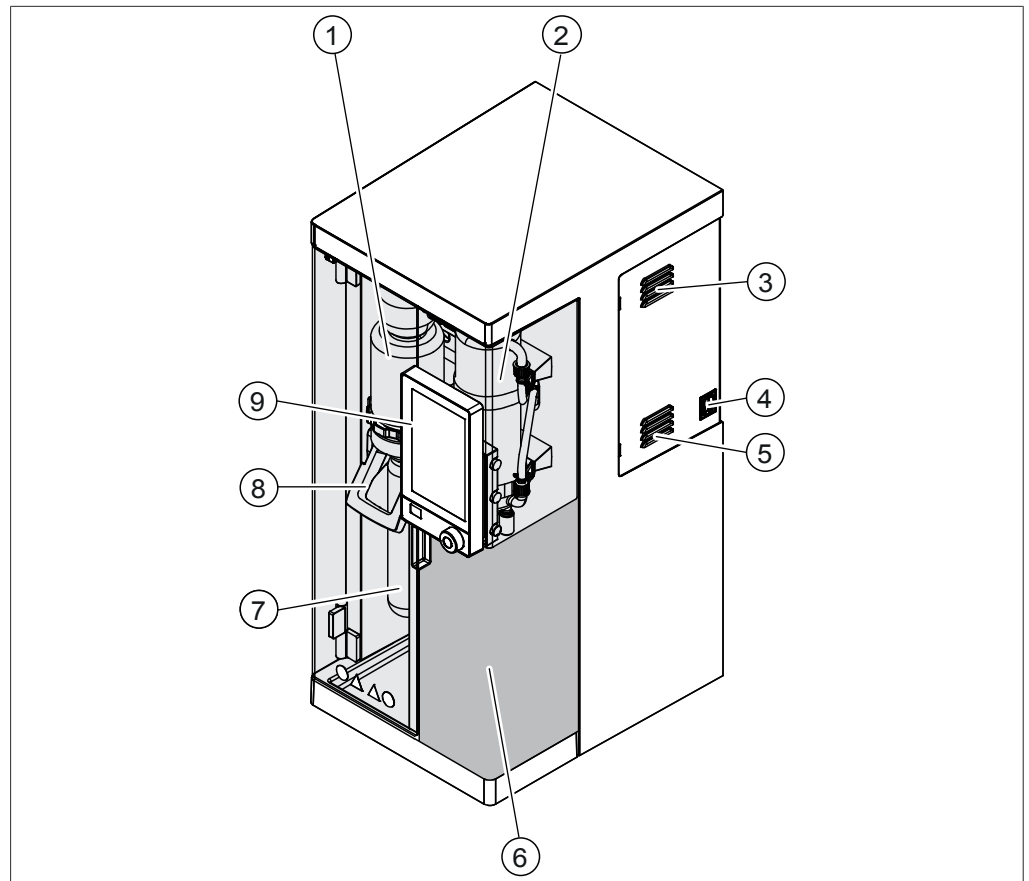


Fig. 3: Front view

- | | | | |
|---|---|---|--|
| 1 | Splash protector | 2 | Condenser |
| 3 | Ventilation slots | 4 | On/Off master switch |
| 5 | Ventilation slots | 6 | Receiving area
See Chapter 3.2.3 "Receiving area",
page 17 |
| 7 | Sample tube | 8 | Handle |
| 9 | Interface
(according to the instrument
configuration) | | |

3.2.2 Rear view

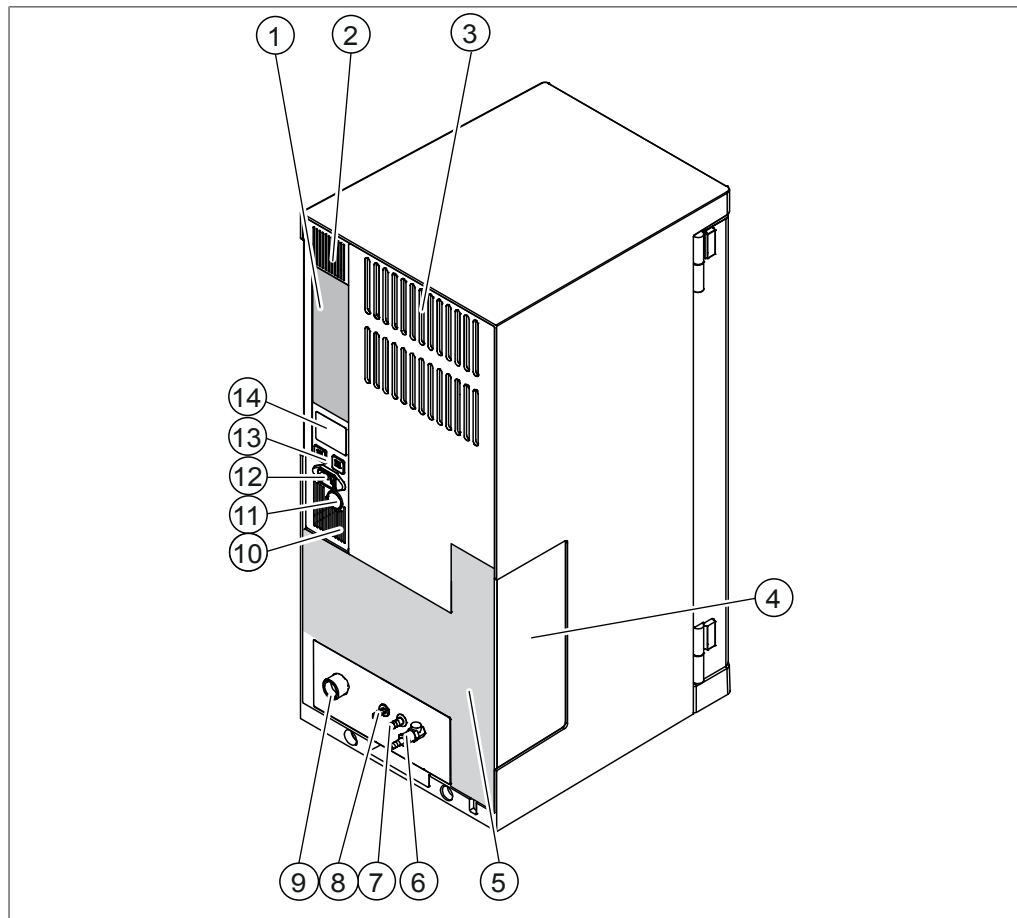


Fig. 4: Rear view

- | | | | |
|----|---|----|---|
| 1 | Communication connections
See Chapter 3.2.4 “Communication connections”, page 18 | 2 | Ventilation slots |
| 3 | Ventilation slots | 4 | Dosing pumps window |
| 5 | Connections on the rear side
See Chapter 3.2.5 “Connections on the rear side”, page 20 | 6 | Steam valve |
| 7 | Cooling water out | 8 | Drain connection
(according to the instrument configuration) |
| 9 | Cooling water in | 10 | Ventilation slots |
| 11 | Rear cable duct | 12 | Power supply connection |
| 13 | Fuses
(resettable) | 14 | Type plate
See Chapter 3.4 “Type plate”, page 24 |

3.2.3 Receiving area

The receiving area is different for each instrument configuration.

Receiving area

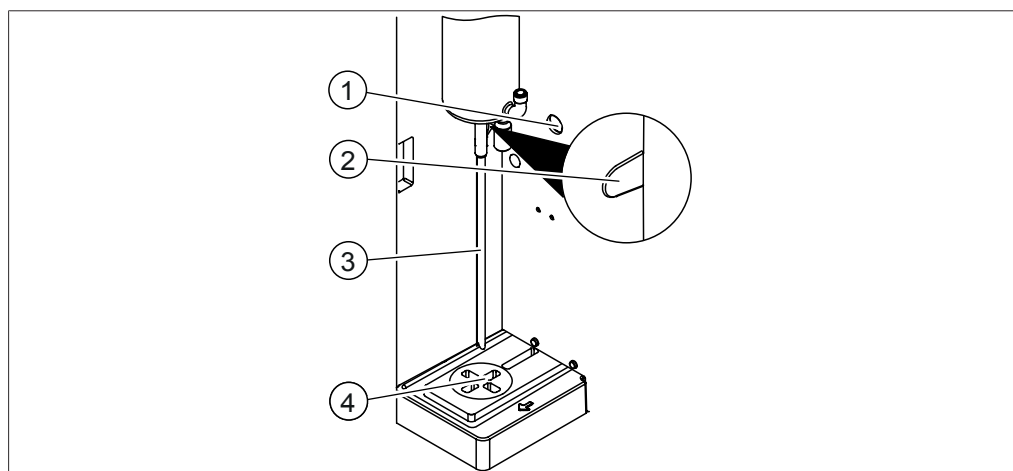


Fig. 5: Receiving area

- | | | | |
|---|------------------|---|--------------------------------------|
| 1 | Rear cable duct | 2 | Cable duct reaction detection sensor |
| 3 | Condenser outlet | 4 | Receiving vessel area |

Receiving area (MultiDist option only)

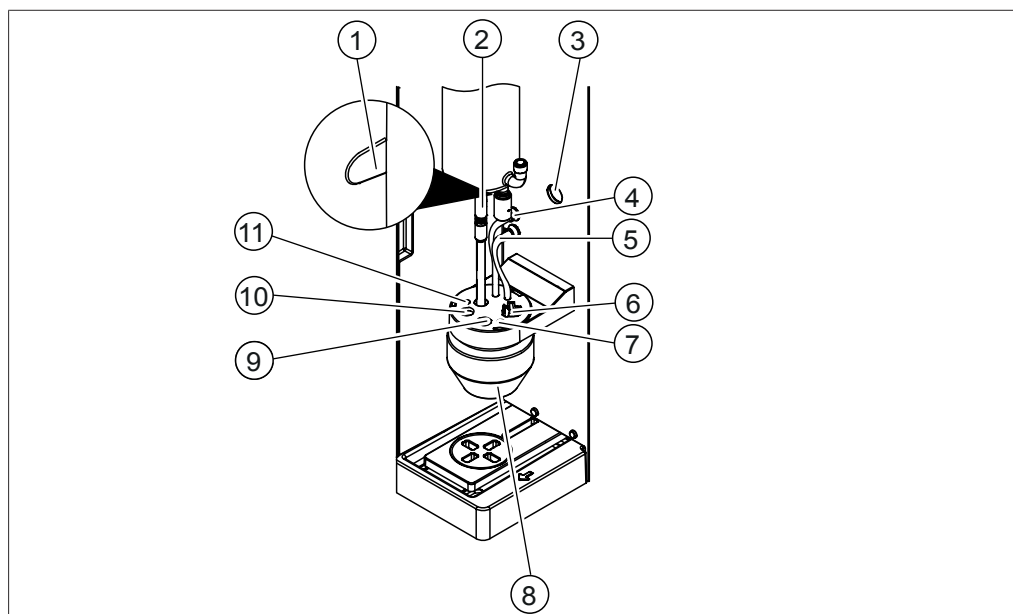


Fig. 6: Receiving area

- | | | | |
|----|--------------------------------------|----|----------------------------------|
| 1 | Cable duct reaction detection sensor | 2 | Condenser outlet tubing |
| 3 | Rear cable duct | 4 | Tubing for boric acid aspiration |
| 5 | Boric acid dosing | 6 | Retainer stirrer |
| 7 | Retainer dosing tip titration | 8 | Receiving vessel |
| 9 | Retainer titration sensor | 10 | Retainer temperature sensor |
| 11 | Retainer dosing tip back titration | | |

3.2.4 Communication connections

Communication connection EasyDist

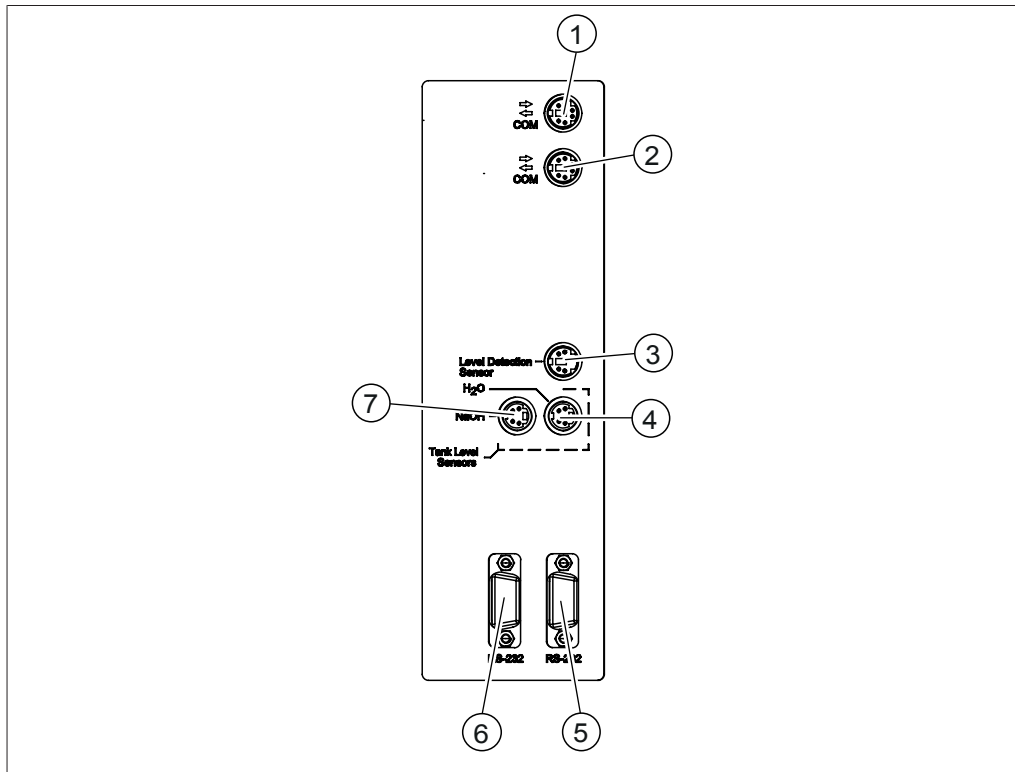


Fig. 7: Communication connections

- | | | | |
|---|---|---|--|
| 1 | BUCHI standard communication port
(marked <i>COM</i>) | 2 | BUCHI standard communication port
(marked <i>COM</i>) |
| 3 | Level detection sensor port
(marked <i>Level Detection</i>) | 4 | H ₂ O tank level sensor port
(marked <i>H₂O</i>) |
| 5 | Communication port
(marked <i>RS-232</i>) | 6 | Communication port
(marked <i>RS-232</i>) |
| 7 | NaOH tank level sensor port
(marked <i>NaOH</i>) | | |

Communication connection BasicDist

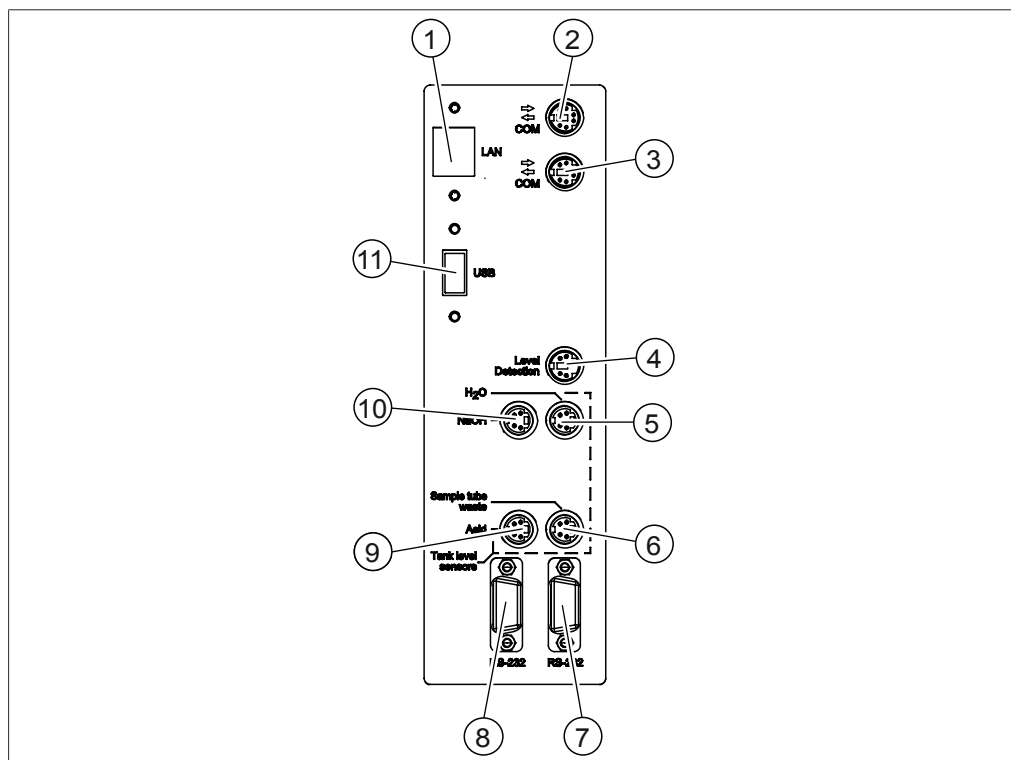


Fig. 8: Communication connection

- | | |
|--|--|
| 1 LAN port
(marked <i>LAN</i>) | 2 BUCHI standard communication port
(COM)
(marked <i>COM</i>) |
| 3 BUCHI standard communication port
(COM)
(marked <i>COM</i>) | 4 OnLevel sensor port
(marked <i>Level Detection</i>) |
| 5 H ₂ O tank level sensor port
(marked <i>H₂O</i>) | 6 Waste tank level sensor port
(marked <i>Sample Tube Waste</i>) |
| 7 Communication port
(marked <i>RS-232</i>) | 8 Communication port
(marked <i>RS-232</i>) |
| 9 Acid tank level sensor port
(marked <i>Acid</i>) | 10 NaOH tank level sensor port
(marked <i>NaOH</i>) |
| 11 USB port
(marked <i>USB</i>) | |

Communication connection MultiDist

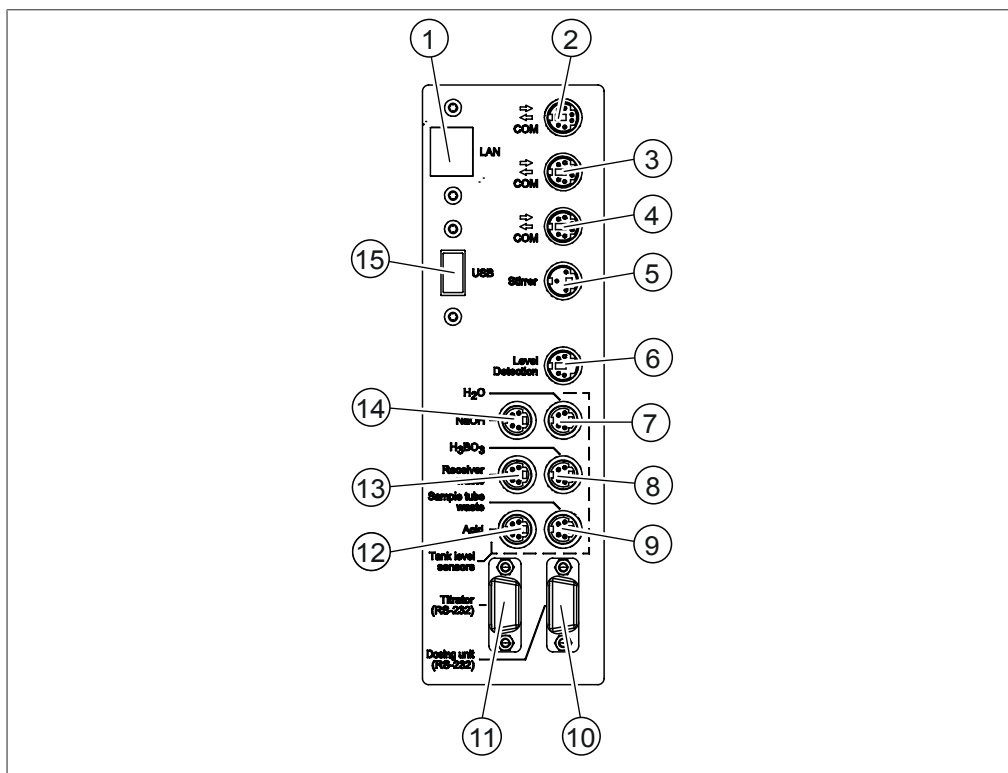


Fig. 9: Communication connections

- | | |
|--|---|
| 1 LAN port
(marked <i>LAN</i>) | 2 BUCHI standard communication port
(marked <i>COM</i>) |
| 3 BUCHI standard communication port
(marked <i>COM</i>) | 4 BUCHI standard communication port
(marked <i>COM</i>) |
| 5 Stirrer port
(marked <i>Stirrer</i>) | 6 OnLevel sensor port
(marked <i>Level Detection</i>) |
| 7 H ₂ O tank level sensor port
(marked <i>H₂O</i>) | 8 H ₃ BO ₃ tank level sensor port
(marked <i>H₃BO₃</i>) |
| 9 Waste tank level sensor port
(marked <i>Sample Tube Waste</i>) | 10 Dispenser port
(marked <i>Dosing Unit</i>) |
| 11 Titrator port
(marked <i>Titrator</i>) | 12 Acid tank level sensor port
(marked <i>Acid</i>) |
| 13 Waste tank level sensor port
(marked <i>Receiver Waste</i>) | 14 NaOH tank level sensor port
(marked <i>NaOH</i>) |
| 15 USB port
(marked <i>USB</i>) | |

3.2.5 Connections on the rear side

The connections on the rear side are depending on the version of the instrument.

Connections on the rear side EasyDist

The connections are located at the rear side of the instrument. See Chapter 3.2.2 “Rear view”, page 16

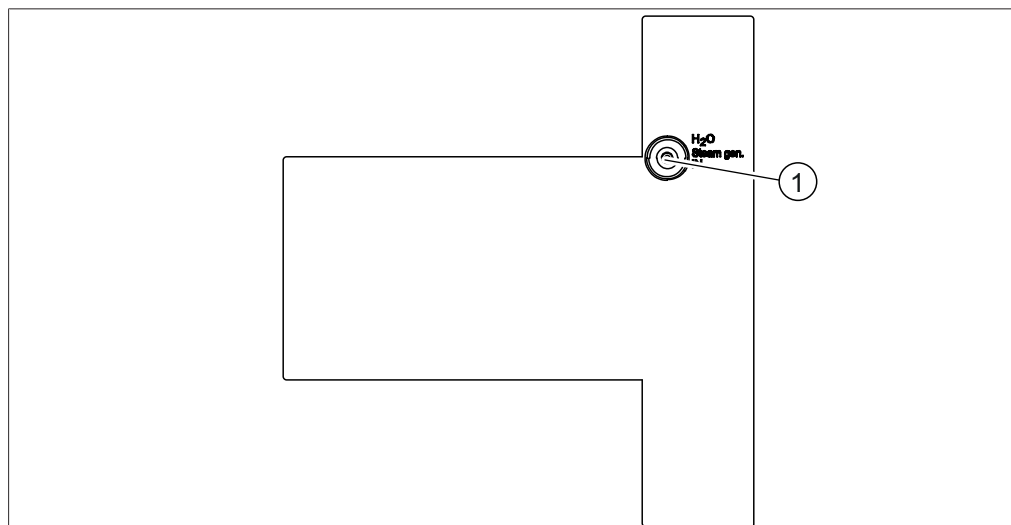


Fig. 10: Connections on the rear side

- 1 H₂O supply for steam generation

Connections on the rear side BasicDist (base unit)

The connections are located at the rear side of the instrument. See Chapter 3.2.2 “Rear view”, page 16

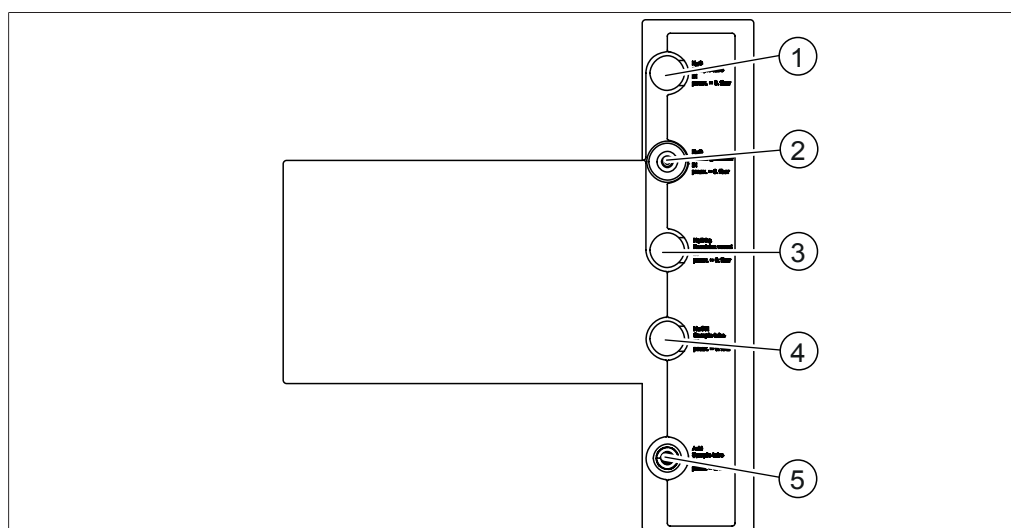


Fig. 11: Connections on the rear side

- | | |
|-------------------------------|--|
| 1 Reserve | 2 H ₂ O supply for steam generation |
| 3 Reserve | 4 Reserve |
| 5 Acid supply for sample tube | |

Connections on the rear side BasicDist (option)

The connections are located at the rear side of the instrument. See Chapter 3.2.2 “Rear view”, page 16

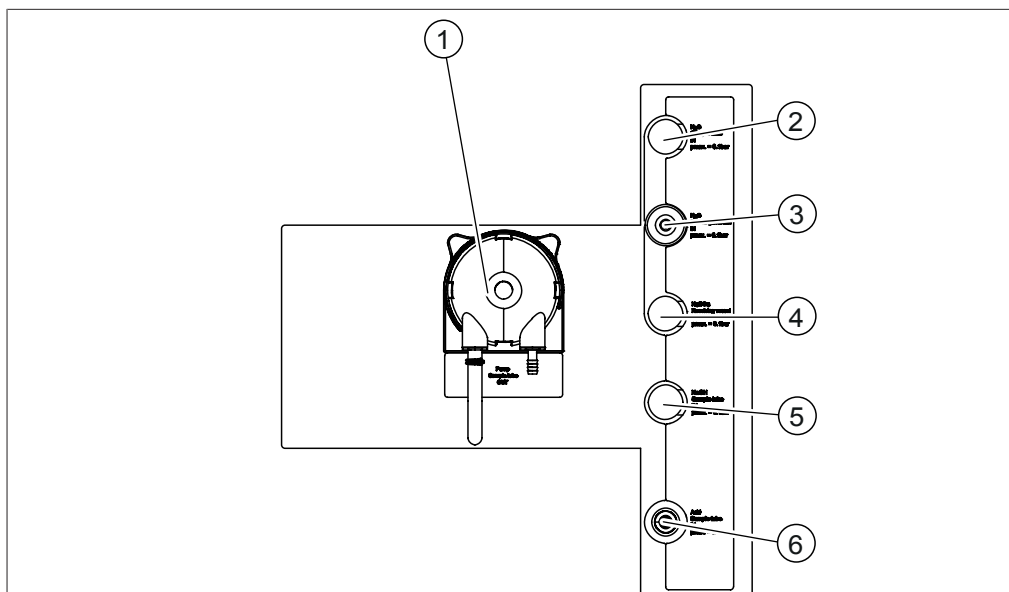


Fig. 12: Connections on the rear side

- | | |
|--|-------------------------------|
| 1 Waste pump for sample tube | 2 Reserve |
| 3 H ₂ O supply for steam generation | 4 Reserve |
| 5 Reserve | 6 Acid supply for sample tube |

Connections on the rear side MultiDist (base unit)

The connections are located at the rear side of the instrument. See Chapter 3.2.2 “Rear view”, page 16

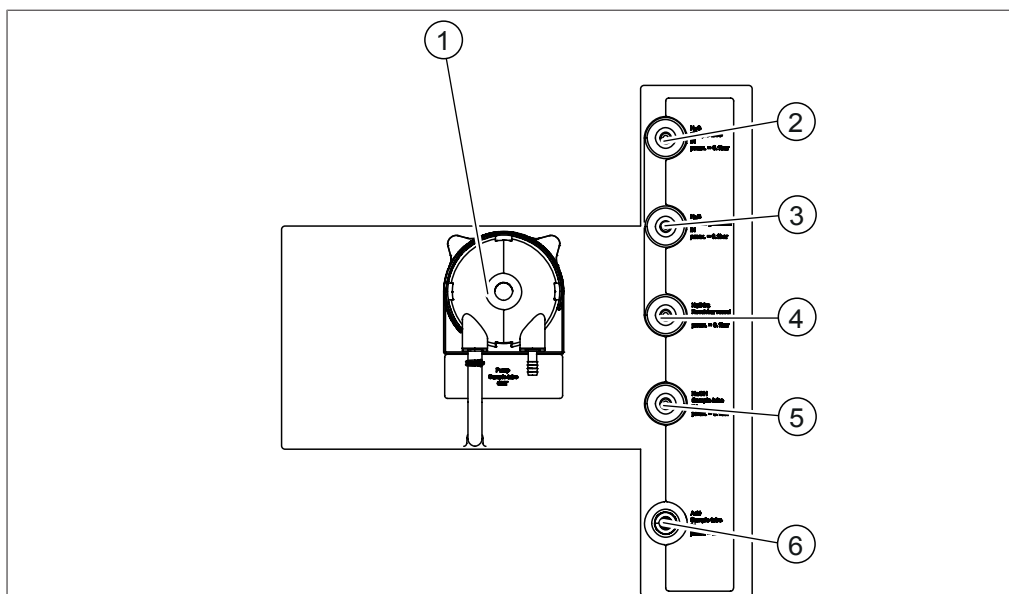


Fig. 13: Connections on the rear side

- | | |
|--|---|
| 1 Waste pump for sample tube | 2 H ₂ O supply for sample tube |
| 3 H ₂ O supply for steam generation | 4 Boric acid supply |
| 5 NaOH supply for sample tube | 6 Acid supply for sample tube |

Connections on the rear side MultiDist (option)

The connections are located at the rear side of the instrument. See Chapter 3.2.2 “Rear view”, page 16

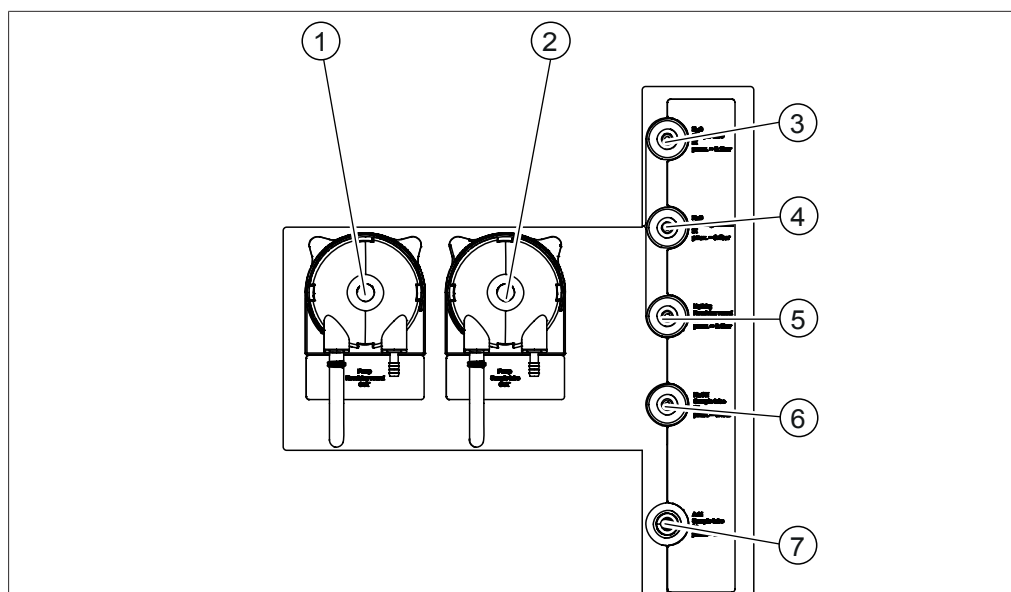


Fig. 14: Connections on the rear side

- | | | | |
|---|---|---|--|
| 1 | Waste pump receiving vessel | 2 | Waste pump for sample tube |
| 3 | H ₂ O supply for sample tube | 4 | H ₂ O supply for steam generation |
| 5 | Boric acid supply | 6 | NaOH supply for sample tube |
| 7 | Acid supply for sample tube | | |

3.3 Scope of delivery



NOTE

The scope of delivery depends of the configuration of the purchase order.

Accessories are delivered as per the purchase order, order confirmation, and delivery note.

3.4 Type plate

The type plate identifies the instrument. The type plate is located at the rear of the instrument.

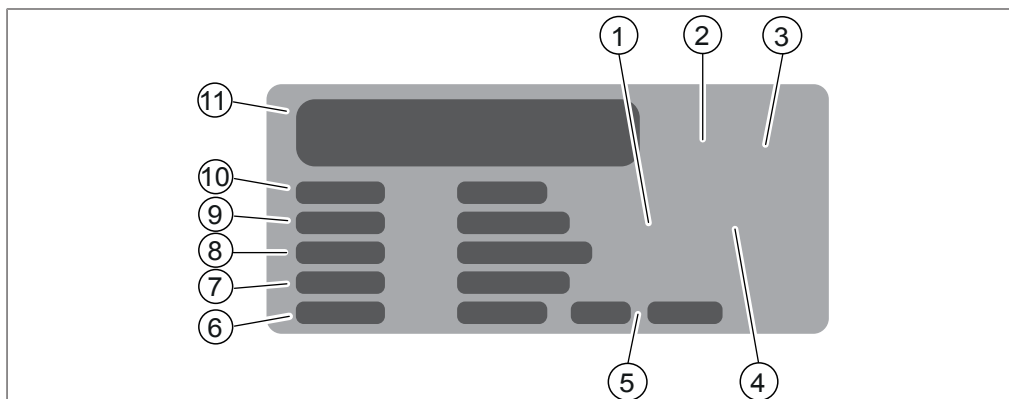


Fig. 15: Type plate

- | | | | |
|----|------------------------------------|----|---|
| 1 | Symbol for "electronics recycling" | 2 | Initial product code |
| 3 | Approvals | 4 | Symbol for "Do not dispose of as household waste" |
| 5 | Year of manufacture | 6 | Power consumption maximum |
| 7 | Frequency | 8 | Input voltage range |
| 9 | Serial number | 10 | Instrument name |
| 11 | Company name and address | | |

3.5 Technical data

3.5.1 K-365 Dist Line

	EasyDist	BasicDist	MultiDist
Dimensions (W x D x H)	320 x 400 x 730 mm	320 x 400 x 730 mm	320 x 400 x 730 mm
Weight	23 kg	23 kg	23 kg
Connection voltage	220 - 240 ± 10 % VAC	220 - 240 ± 10 % VAC	220 - 240 ± 10 % VAC
Power consumption	2100 W	2100 W	2100 W
Frequency	50 / 60 Hz	50 / 60 Hz	50 / 60 Hz
IP Code	IP20	IP20	IP20
Pollution degree	2	2	2
Overvoltage category	II	II	II
Tap water pressure	1 - 10 bar	1 - 10 bar	1 - 10 bar
Tap water condition	1- 25°C	1 - 25°C	1 - 25°C
(If the temperature is more than 25°C use a recirculating chiller)			
Tap water consumption	~ 1.2 L/min	~ 1.2 L/min	~ 1.2 L/min
Approvals	CE, CSA	CE, CSA	CE, CSA

Analysis specifications

	EasyDist	BasicDist	MultiDist
Measuring range	analyte specific	analyte specific	0.02 – 220 mg N
Analysis time for 30 mg N (with Metrohm Eco Titrator without titrator preparation)	-	-	3.5 min
Analysis time for 200 mg N (with Metrohm Eco Titrator without titrator preparation)	-	-	5 min
Recovery	analyte specific	analyte specific	N: >98% with digestion process N: >99.5% (direct distillation)
Measurement reproducibility (Direct distillation of 1 mg N absolute and 300 sec. distillation)	analyte specific	analyte specific	N: <0.8%

Instrument specifications

	EasyDist	BasicDist	MultiDist
Distillation capacity (100%)	~ 40 mL/min	~ 40 mL/min	~ 40 mL/min
Distillation capacity (10%)	~ 12.5 mL/min	~ 12.5 mL/min	~ 12.5 mL/min
Adjustable distillation power	10 - 100%	10 - 100%	10 - 100%
MaxAccuracy Mode (Compensation of power grid fluctuations)	Yes	Yes	Yes
Distillate reproducibility (RSD) (At 300 seconds distillation time)	<1%	<1%	<1%
AutoDist Mode (Automatic detection of the condensation start)	No	Yes	Yes
NaOH pump	No	No	Yes
H ₂ O pump	No	No	Yes
Acid pump	No	Yes	Yes
Boric acid pump	No	No	Yes
Sample disposal	No	Optional	Yes
Receiver disposal	No	No	Optional
Safety sensors	Yes	Yes	Yes
Titration vessel	No	No	Optional
Titrator connection	No	No	Optional
IQ/OQ	No	Yes	Yes

Interface specifications

	Interface EasyDist, Ba- sicDist	Interface pro BasicDist (option), MultiDist
Display	LCD color, 4.3 in	LCD color, 7.0 in
Control elements	Buttons / control knob	Touch-screen / control knob
Max. number of methods	8	96
Max. number of determinations	36	512
Max. number of determinations per series	16	32
Max. number of templates	4	16
Max. number of users	4	32
Max. number of determination results	40	256
Language	en, de, fr, it, es, pt, ja, zh, ru, pl, ko, id	en, de, fr, it, es, pt, ja, zh, ru, pl, ko, id

Connections

	EasyDist	BasicDist	MultiDist
LAN port	-	-	1
BUCHI standard communication ports (COM)	2	2	3
Stirrer port	-	-	1
Level detection sensor port	1	1	1
H ₂ O tank level sensor port	1	1	1
H ₃ BO ₃ tank level sensor port	-	-	1
Acid tank level sensor port	-	-	-
NaOH tank level sensor port	1	1	1
Waste tank level sensor port	-	option	2
Dispenser port (RS-232)	-	-	1
Titration port (RS-232)	-	-	1
Titration port (RS-232)	-	-	1

3.5.2 Ambient conditions

For indoor use only.

Max. altitude above sea level	2,000 m
Ambient and storage temperature	5–40 °C
Maximum relative humidity	80% for temperatures up to 31 °C decreasing linearly to 50 % relative humidity at 40 °C, non-condensing

3.5.3 Materials

Component	Materials of construction
Housing	Polyurethane
Housing electric	Stainless steel
Steam generator housing	Stainless steel
Housing stirrer	Glass fiber reinforced polypropylene
Rear cover	Polyvinyl chloride
Glass parts	Borosilicate 3.3
Plastic splash protector	Glass fiber reinforced polypropylene
Protective cover	PMMA
Sealing	CSM
NaOH hose	EPDM
Three-way valve	PP / PE

3.5.4 Installation site

- The installation site meets the safety requirements. See Chapter 2 “Safety”, page 9
- The installation site has a firm, level and nonslip surface.
- The installation site has no obstacles (e.g. water taps, drains, etc.).
- The installation site has enough space that the canisters can be placed safely.
- The installation site has an own mains outlet socket for the instrument.
- The installation site allows that the power supply can be disconnected at any time in an emergency.
- The installation site is not exposed to external thermal loads, such as direct solar radiation.
- The installation site has enough space that cables / tubes can be routed safely.
- The installation site meets the requirements for the connected devices. See related documentation
- The installation site meets the specifications according to the technical data (e.g. weight, dimension, etc.). See Chapter 3.5 “Technical data”, page 24

4 Transport and storage

4.1 Transport



NOTICE

Risk of breakage due to incorrect transportation

Make sure that the instrument is fully dismantled.

Pack every instrument components properly to prevent breakage. Use the original packaging whenever possible.

Avoid sharp movements during transit.

- ▶ After transporting, check the instrument and all glass components for damage.
- ▶ Damage that has occurred in transit should be reported to the carrier.
- ▶ Keep packaging for future transportation.

4.2 Storage

- ▶ Make sure that the ambient conditions are complied with (see Chapter 3.5 “Technical data”, page 24).
- ▶ Wherever possible, store the device in its original packaging.
- ▶ After storage, check the device, all glass components, seals and tubing for damage and replace if necessary.

4.3 Lifting the instrument




NOTICE

Dragging the instrument can damage the feet of the instrument.

- ▶ Lift the instrument when positioning or re-locating.

- ▶ Lift the instrument at the points indicated.

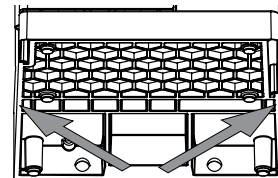

5 Installation

5.1 Before installation



NOTICE

Instrument damage due to switching it on too early.

Switching on the instrument too early after transportation can cause damage.

- ▶ Climatize the instrument after transportation.

5.2 Establishing electrical connections



NOTICE


Risk of instrument damage because of not suitable power supply cables.

Not suitable power supply cables can cause bad performance or an instrument damage

- ▶ Use only BUCHI power supply cables.

Precondition:


- The electrical installation is as specified on the type plate.
 - The electrical installation is equipped with a proper grounding system.
 - The electrical installation is equipped with suitable fuses and electrical safety features.
 - The installation site is as specified in the technical data. See Chapter 3.5 "Technical data", page 24
- ▶ Connect the power supply cable to the connection on the instrument. See Chapter 3.2 "Configuration", page 14
 - ▶ Connect the mains plug to an own mains outlet socket.

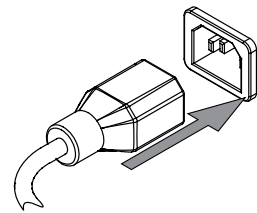

5.3 Securing against earthquakes

The instrument has an earthquake fixing point to protect the device against falling.

- ▶ Tie the lashing mount to a fixed point using strong cord or a wire.


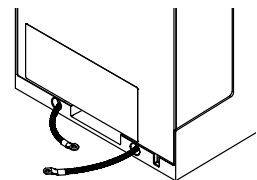


5.4 Installing the cooling water supply

5.4.1 Installing a tap water supply (option)

Precondition:

- ☑ The tap water supply complies with the specified parameters. See Chapter 3.5 “Technical data”, page 24
- ☑ Make sure that the instrument is not connected to the power supply.
- ▶ Install the inlet hose to the connection marked *Cooling Water / Chiller IN.*
- ▶ Install the drain hose to the connection marked *OUT Cooling Water / Chiller.*

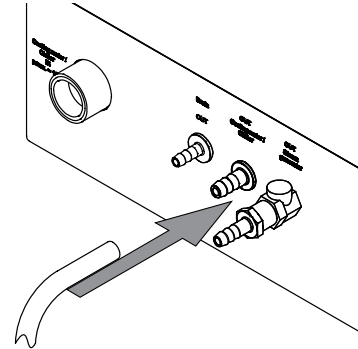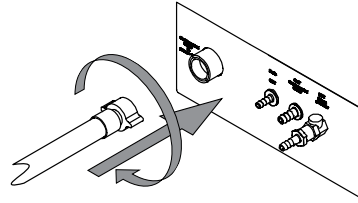

- ▶ Put the other end of the hose in a sink.

5.4.2 Installing the recirculating chiller on the condenser (option)





NOTE


Place the hose in the rear cable duct.

Precondition:

- ☑ The water supply complies with the tap water specifications. See Chapter 3.5 “Technical data”, page 24
- ▶ Make sure that the instrument is not connected to the power supply.
- ▶ Remove the connection at the condenser inlet.
- ▶ Attach the cooling water hose from the chiller to the condenser.

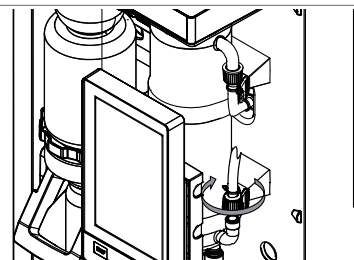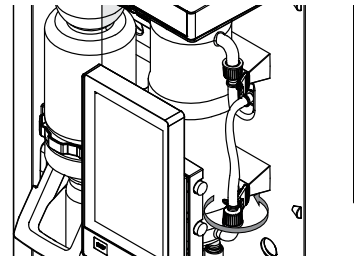


- ▶ Install the drain hose to the connection marked *OUT Cooling Water / Chiller* on the rear of the instrument.

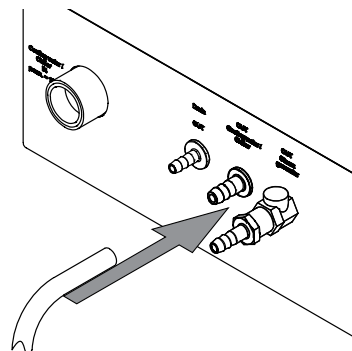

- ▶ Make sure that the hoses are not bend.


5.5 Installing the drain tubing (option)

- ▶ Install the drain hose to the connection marked *Drain*.
- ▶ Place the other end of the hose in an aqueous waste container located lower than the instrument.

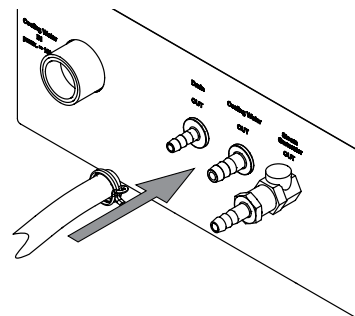

5.6 Installing the waste pump connection (MultiDist option only)


- ▶ Attach the waste hose onto the connection marked *Pump Receiving Vessel OUT*.
- ▶ Secure the hose in place with a hose clip.

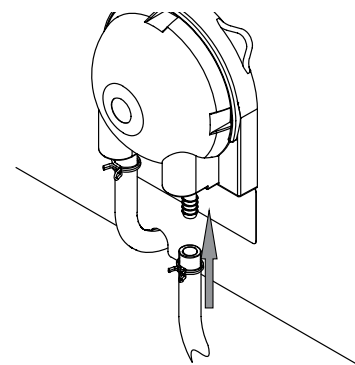

- ▶ Put the other end of the hose in a suitable collection device.

5.7 Installing the sample waste pump connection (BasicDist option, MultiDist only)

- ▶ Attach the waste hose to the connection marked *Pump Sample Tube OUT*.
- ▶ Secure the hose in place with a hose clip.

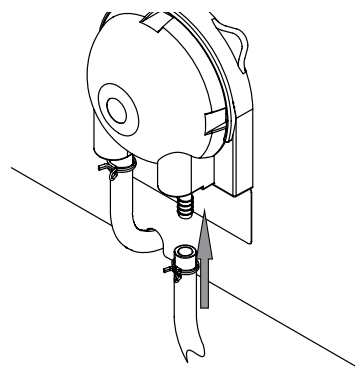

- ▶ Put the other end of the hose in a suitable collection device.

5.8 Installing the H₂O supply for steam generation




NOTE

MultiDist only


The water supply comes from one canister.

Precondition:


- ☑ The water complies with the specified parameters. See Chapter 3.5 “Technical data”, page 24
- ☑ Make sure that the instrument is not connected to the power supply.
- ▶ Install the inlet hose to the connection marked *H₂O Steam Gen. IN*.
- ▶ Attach the inlet hose in place with a hose clip.

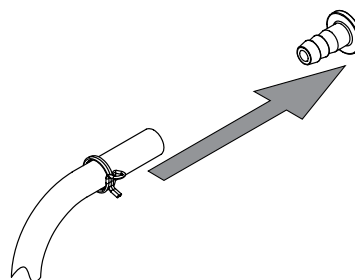


- ▶ Moist the aspiration hose.
- ▶ Push the hoses together.

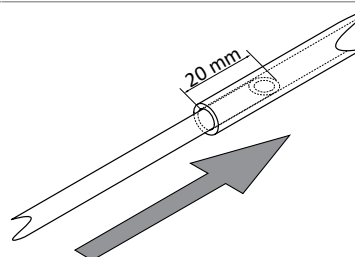

- ▶ Attach the hose grommet to the canister lid.

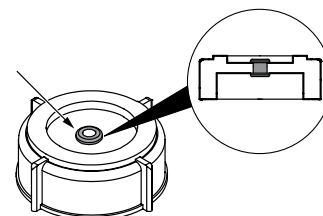


- ▶ Moist the hose.
- ▶ Insert the inlet hose through the opening of the cap nut and the screw cap of the water tank.
- ▶ Submerge the inlet hose into the water.

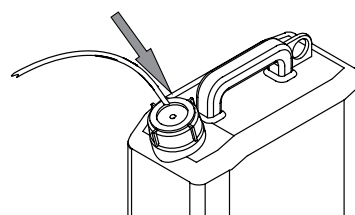


5.9 Installing NaOH supply (MultiDist only)


- ▶ Install the inlet hose onto the connection marked *NaOH Sample Tube IN.*
- ▶ Attach the inlet hose in place with a hose clip.

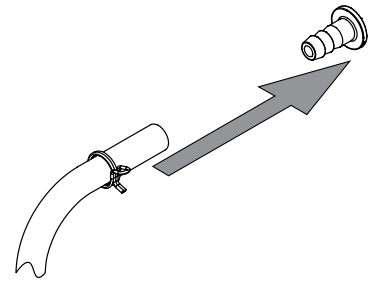


- ▶ Moist the aspiration hose.
- ▶ Push the hoses together.

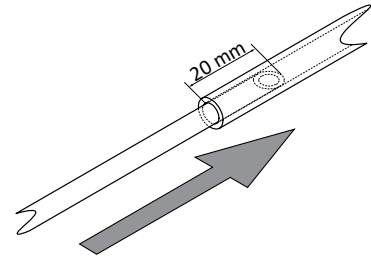

- ▶ Attach the hose grommet to the canister lid.

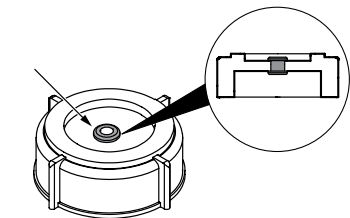


- ▶ Submerge the inlet hose into the NaOH.

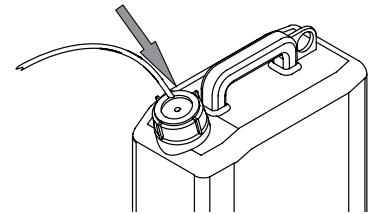


5.10 Installing H₃BO₃ supply (MultiDist only)


- ▶ Install the inlet hose to the connection marked *H₃BO₃ Receiving Vessel IN.*
- ▶ Attach the inlet hose in place with a hose clip.

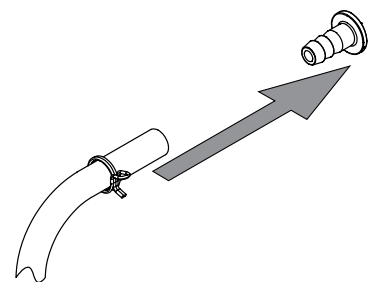


- ▶ Moist the aspiration hose.
- ▶ Push the hoses together.

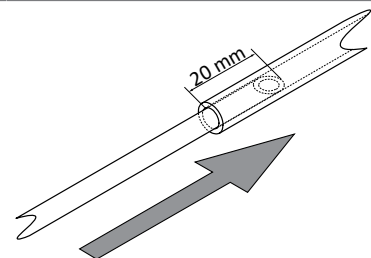

- ▶ Attach the hose grommet to the canister lid.

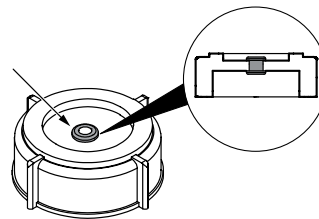

- ▶ Submerge the inlet hose into the H₃BO₃.

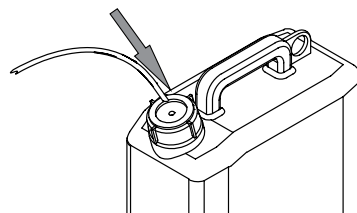

5.11 Installing the H₂O supply for the sample tube (MultiDist only)




NOTE


MultiDist only

The water supply comes from one canister.


- ▶ Install the inlet hose to the connection marked *H₂O Sample Tube IN.*
- ▶ Attach the inlet hose in place with a hose clip.

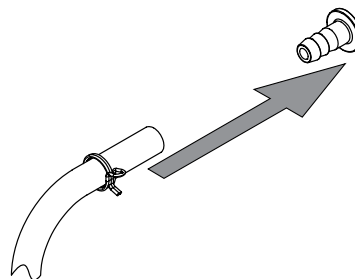


- ▶ Moist the aspiration hose.
- ▶ Push the hoses together.

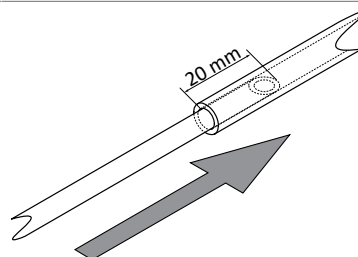

- ▶ Attach the hose grommet to the canister lid.

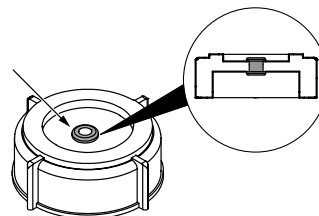

- ▶ Submerge the inlet hose into the H₂O.

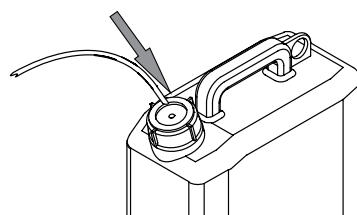


5.12 Installing the level sensor on the canister (option)




NOTE

Do not expose the tank level sensor to an electromagnetic field in the frequency range of 2 to 10 MHz.


- ▶ Attach the hook-and-loop fastener sticker to the sensor.

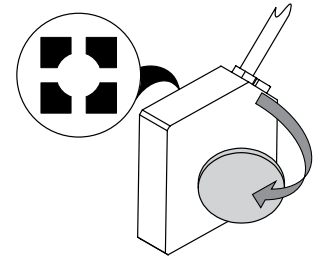


- ▶ Attach the strap to the canister.

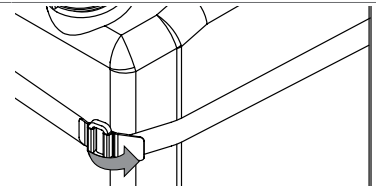

- ▶ Fix the sensor on the canister.

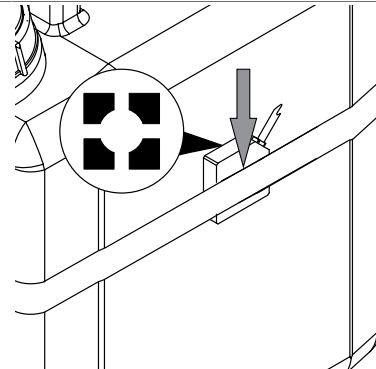

- ▶ Connect the sensor cable to the related connection on the instrument. See Chapter 3.2 "Configuration", page 14

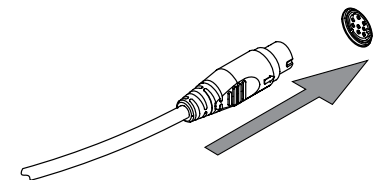

5.13 Installing a titrator with LAN (MultiDist option with Eco Titrator only)


Settings




NOTE

Make sure, that the IP on both instruments is 192.168.10.3


- ▶ Connect the LAN cable to the related connection on the instrument. See Chapter 3.2 "Configuration", page 14

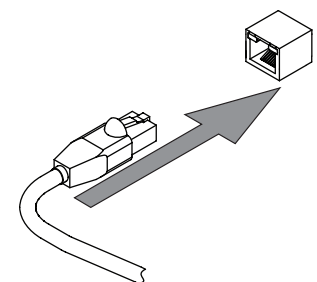


Specifications on the display:	
Precondition:	
<ul style="list-style-type: none"> ☑ The titrator is prepared. See related manual ▶ Navigate to the specified submenu. 	<div style="text-align: center;">↓</div> <div style="border: 1px solid black; padding: 2px; display: inline-block;">Peripherals ></div>
	<div style="text-align: center;">↓</div> <div style="border: 1px solid black; padding: 2px; display: inline-block;">Titrator Model ></div>
▶ Select the specified action.	Titrator Model ...
▶ Tap the specified function on the function bar.	EDIT
▶ Select the titrator model you wish to use.	<div style="display: flex; align-items: center;"> <div style="border: 1px solid black; border-radius: 10px; padding: 2px 10px; margin-right: 5px;">...</div> <div style="border: 1px solid black; padding: 2px;"> <div style="border: 1px solid black; border-radius: 5px; padding: 2px; margin-bottom: 2px;">...</div> <div style="border: 1px solid black; border-radius: 5px; padding: 2px; margin-bottom: 2px;">...</div> <div style="border: 1px solid black; border-radius: 5px; padding: 2px;">...</div> </div> </div>
▶ Tap the specified function on the function bar.	SAVE
⇒ The titrator model is saved.	

5.14 Installing a titrator with RS232 (MultiDist option only)

- Precondition:
- ☑ The dosing unit is prepared. See related documentation
 - ▶ Attach the plug onto the connection marked *Titrator (RS-232)*.
 - ▶ Secure the plug in place.

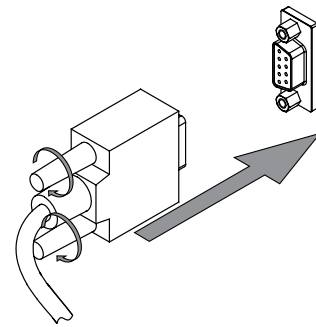


Specifications on the display:	
Precondition:	
<ul style="list-style-type: none"> ☑ The titrator is prepared. See related manual ☑ The instrument is connected to the LAN network. ▶ Navigate to the specified submenu. 	<div style="text-align: center;">↓</div> <div style="border: 1px solid black; padding: 2px; display: inline-block;">Peripherals ></div>
	<div style="text-align: center;">↓</div> <div style="border: 1px solid black; padding: 2px; display: inline-block;">Titrator Model ></div>
▶ Select the specified action.	Titrator Model ...
▶ Tap the specified function on the function bar.	EDIT
▶ Select the titrator model you wish to use.	<div style="display: flex; align-items: center;"> <div style="border: 1px solid black; border-radius: 10px; padding: 2px 10px; margin-right: 5px;">...</div> <div style="border: 1px solid black; padding: 2px;"> <div style="border: 1px solid black; border-radius: 5px; padding: 2px; margin-bottom: 2px;">...</div> <div style="border: 1px solid black; border-radius: 5px; padding: 2px; margin-bottom: 2px;">...</div> <div style="border: 1px solid black; border-radius: 5px; padding: 2px;">...</div> </div> </div>
▶ Tap the specified function on the function bar.	SAVE
⇒ The titrator model is saved.	

5.15 Installing a dispenser (MultiDist only)


Precondition:

- ☑ The dosing unit is prepared. See related Documentation
- ▶ Attach the plug onto the connection marked *Dosing Unit*.
- ▶ Secure the plug in place.

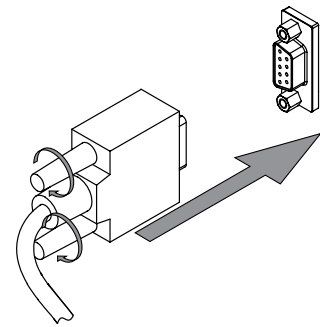


5.16 Installing the reaction detection sensor (MultiDist option only)

- ▶ Attach the reaction detection sensor to the instrument.

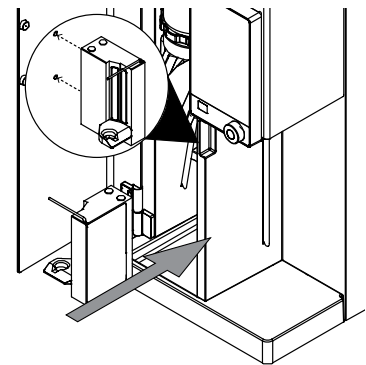


- ▶ Connect the sensor cable to the related connection on the instrument. See Chapter 3.2 "Configuration", page 14

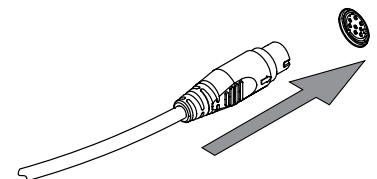

5.17 Installing a SD card (Interface pro only)

- ▶ Insert the SD card.
- ▶ Restart the instrument.
- ⇒ The status bar shows the SD card symbol.

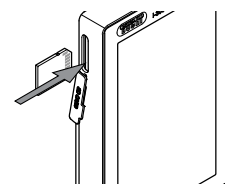


5.18 Installing the OnLevel sensor (option)




NOTE

Using the OnLevel sensor. See Chapter 8.11 "Preparing the OnLevel sensor (option)", page 71


- ▶ Attach the metal plate to the magnetic area of the instrument.

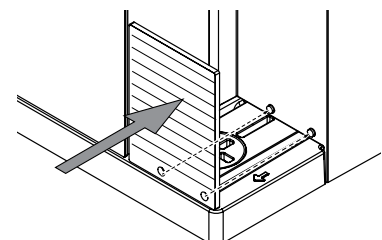

- ▶ Push the cable through the rear cable duct.

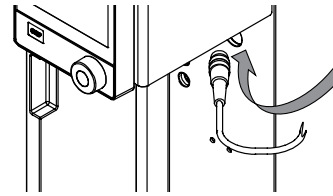


- ▶ Connect the sensor cable to the related connection on the instrument. See Chapter 3.2 “Configuration”, page 14

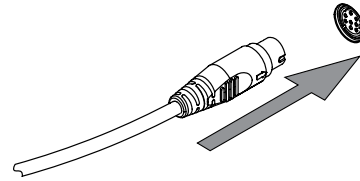


5.19 Installing the stirrer (MultiDist option only)


- ▶ Attach the stirrer to the receiving vessel.

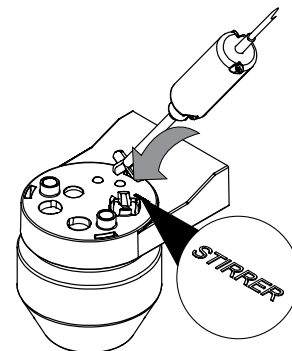

- ▶ Push the cable through the rear cable duct.

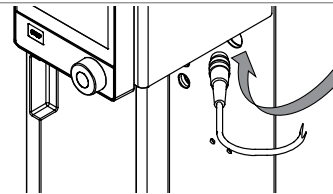

- ▶ Connect the sensor cable to the related connection on the instrument. See Chapter 3.2 “Configuration”, page 14

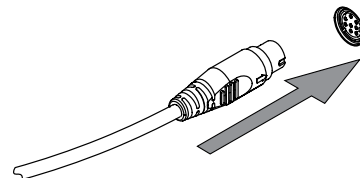

6 Description of the interface (EasyDist and BasicDist)

6.1 Layout of the interface


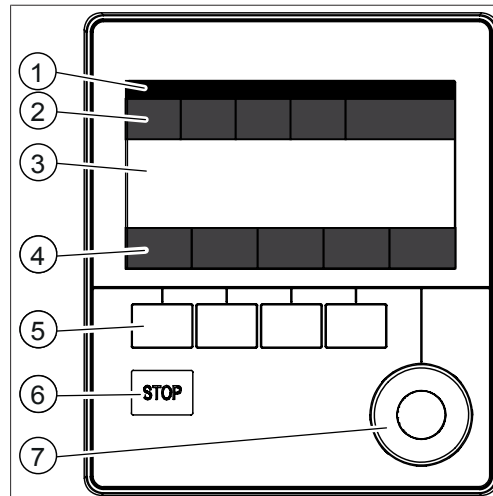


Fig. 16: Interface


No.	Description	Function
1	Status bar	Shows the determination name and status of the instrument. See Chapter 6.4 "Status bar interface", page 43
2	Menu bar	Shows symbols representing the menus. See Chapter 6.3 "Menu bar interface", page 40
3	Content area	Shows current settings, submenus or actions depending on the current operation.
4	Function bar	Shows functions that can be performed according to the current operation. See Chapter 6.2 "Function bar interface", page 40
5	Function buttons	Pressing a function button performs the assigned function on the function bar.
6	Stop button	
7	Navigation control	Used for navigating the user interface.

Using the navigation control

- Select an entry.

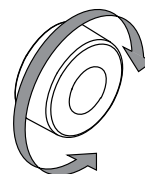




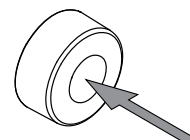





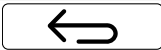

- Confirm the selected entry.








6.2 Function bar interface

The function bar shows available functions according to the current operation. The functions on the function bar are executed by tapping the relevant function buttons.

Symbol	Description	Meaning
	[Back]	The display reverts to the previous view.
	[Confirm]	Confirms an entry.
	[Start]	Starts a steam distillation process.
	[Copy]	Copies the selected method.
	[Edit]	Allows the selected item to be edited.
	[Ready]	Steam generator is ready to distill.
	[Preparation]	Carries out the method selected in the [Priming] submenu. See Chapter 8.2 "Editing the priming function", page 50
	[Template]	Saves the selected series as a template. Creates a series from the selected template.
	[Delete]	Deletes the selected entry.
	[Delete all]	Deletes all entries.
	[Standby]	The steam generator changes to standby mode.

6.3 Menu bar interface

Menu symbol	Name	Explanation
	Home menu	<ul style="list-style-type: none"> • Process control parameters
	Manual control menu	<ul style="list-style-type: none"> • Start various processes manually
	Process menu	<ul style="list-style-type: none"> • Editing and saving: <ul style="list-style-type: none"> ◦ Methods ◦ Series ◦ Templates ◦ Determinations
	Configuration menu	<ul style="list-style-type: none"> • Change settings • Service menu • System information
	Determination data menu	<ul style="list-style-type: none"> • View processed determinations

6.3.1 Home menu

6.3.2 Manual control menu

The manual control menu contains the following submenus:

Submenu	Explanation
Dose manually	See Chapter 8.13 "Dosing manually", page 72
Aspirate manually	See Chapter 8.14 "Aspirating manually", page 72
<i>[Preheating]</i>	See Chapter 8.1 "Editing the preheating function", page 50
<i>[Priming]</i>	See Chapter 8.2 "Editing the priming function", page 50
<i>[Cleaning]</i>	See Chapter 10.3 "Cleaning the glass components", page 79
<i>[Stirring]</i>	See Chapter 8.15 "Changing the stirring speed manually", page 73
<i>[Pump Calibration]</i>	See Chapter 10.14 "Calibrating the pumps", page 81

6.3.3 Method menu

The method menu contains the following submenus:

Submenu	Explanation
<i>[Single Determination]</i>	See Chapter 8.3 "Editing the single determination", page 50
<i>[Series]</i>	See Chapter 8.4 "Editing a series", page 53
<i>[Templates]</i>	See Chapter 8.5 "Editing a template", page 57
<i>[Methods]</i>	See Chapter 8.6 "Editing a method", page 58

6.3.4 Configuration menu

Settings submenu

Action	Option	Explanation
<i>[Language]</i>	Choice of display language on the interface	The following languages are available: English/German/French/Italian/Spanish/Japanese/Chinese/Russian/Polish
<i>[Current Date]</i>	Date input	Enter in sequence: Day, month, year. Apply the settings by pressing <i>[Save]</i> .
<i>[Current Time]</i>	Time input	Enter in sequence: Minutes, hours. Apply the settings by pressing <i>[Save]</i> .
Time zone	Choice of time zone	Select your time zone
<i>[Keyboard Layout]</i>	Choice of display keyboard on the interface	Select the keyboard layout
<i>[Key Tones]</i>	Off/On	Setting for audible signal in response to input controls.
<i>[Beep on Finish]</i>	Off/On	Setting for audible signal by the end of a determination.

Action	Option	Explanation
<i>[Beep on Error]</i>	Off/On	Setting for audible signal by the end of a determination.
<i>[Display Brightness]</i>	Enter setting	Display illumination level in %: 0 - 100
<i>[Zero Amount Warning]</i>	On/Off	The instrument gives a warning when the entry for the weight is zero.
<i>[Default Amount Unit]</i>	Select value	The following values are selectable: g, mL
<i>[MaxAccuracy Mode]</i>	On / Off	Compensation of the steam power caused by voltage fluctuations.
<i>[AutoDist Mode]</i>	On/Off	Conditioning and distillation are performed automatically.
<i>[Level Detect. Max. Time]</i>	Enter value	The time after which the distillation stops without level detection sensor triggering.
<i>[Pump Maintenance Interval]</i>	Select value	Selecting a frequency to carry out a pump rinsing. See Chapter 10.21 "Rinsing a pump", page 86
<i>[H₂O Dosage per Keypress]</i>	Enter value	Dosage volume which is applied by pressing the related button.
<i>[Acid Dosage per Keypress]</i>	Enter value	Dosage volume which is applied by pressing the related button.
<i>[NaOH Dosage per Keypress]</i>	Enter value	Dosage volume which is applied by pressing the related button.
<i>[Aspiration Time Receiver]</i>	Select value	Aspiration time for the related button.
<i>[Aspiration Time Sample Tube]</i>	Select value	Aspiration time for the related button.
<i>[H₃BO₃ Dosage per Keypress]</i>	Enter value	Dosage volume which is applied by pressing the related button.
<i>[Steam Generator Standby]</i>	Enter value	Enter the time until the instrument turns automatically into standby.
<i>[Demo Mode]</i>	On/Off	Simulate a distillation.
<i>[Service Data Logger]</i>	On/Off	Saves all data on the instrument for 14 months.

Peripherals submenu

Shows information about the connected peripherals.

Users submenu

Creating and setting users. See Chapter 8.7 "Editing user setting", page 68

Network submenu

Action	Option	Explanation
[Network]	Enter value	The following parameters can be edited: Device name/MAC address/DHCP/ System IP address/Subnet mask/ Gateway/DNS server/BUCHI Cloud/ Server IP address

Service submenu



NOTE

Only super user can carry out changes in the [Service] submenu.

System Information submenu

Technical information about the connected devices (e.g. serial number, firmware version).

6.3.5 Determination data menu

Shows the history of the determination data.

6.4 Status bar interface

The status bar shows actual information about the instrument.

View	Status
	The instrument is in standby mode.
	Process is running.
	The instrument is heating up.
	The BLE dongle is connected to the instrument.
	A level detection sensor is connected to the instrument.
	Reaction detection sensor attached.
	A tank level sensor is connected to the instrument.
	Dosing unit attached.
	Titration unit attached.
	White: A SD card is in the instrument. Red: The SD card in the instrument is in read only mode.
	A USB input device is connected.
	A USB memory stick is connected to the instrument.
	A recirculating chiller is connected.

7 Description of the interface pro (BasicDist and MultiDist)



⚠ CAUTION

Risk of injury from glass splinters

Sharp objects can damage the display.

- ▶ Keep sharp objects away from the display.

7.1 Layout of the interface pro


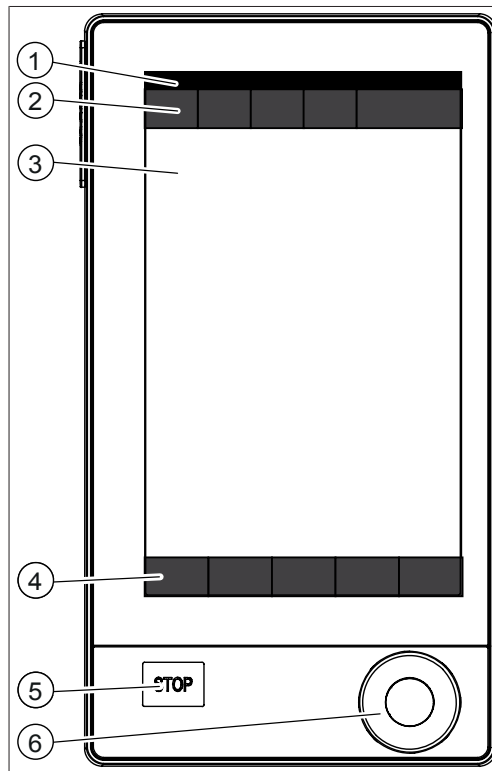


Fig. 17: Interface pro


No.	Description	Function
1	Status bar	Shows the determination name and status of the instrument. See Chapter 7.4 "Status bar interface pro", page 49
2	Menu bar	Shows symbols representing the menus.
3	Content area	Shows current settings, submenus or actions depending on the current operation.
4	Function bar	Shows functions that can be performed according to the current operation.
5	Stop button	
6	Navigation control	Used for navigating the user interface.

Using the navigation control

- ▶ Select an entry.

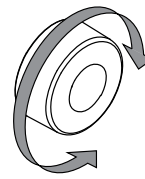



- ▶ Confirm the selected entry.

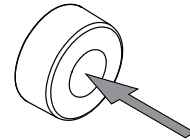



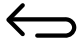


7.2 Function bar interface pro

The function bar shows available functions according to the current operation. The functions on the function bar are executed by tapping the relevant function buttons.

Symbol	Description	Meaning
	<i>[Back]</i>	The display reverts to the previous view.
OK	<i>[Confirm]</i>	Confirms an entry.
START	<i>[Start]</i>	Starts a steam distillation process.
COPY	<i>[Copy]</i>	Copies the selected method.
EDIT	<i>[Edit]</i>	Allows the selected item to be edited.
READY	<i>[Ready]</i>	Steam generator is ready to distill.
PREP	<i>[Preparation]</i>	Carries out the method selected in the <i>[Priming]</i> submenu. See Chapter 8.2 "Editing the priming function", page 50
	<i>[Template]</i>	Saves the selected series as a template. Creates a series from the selected template.
DELETE	<i>[Delete]</i>	Deletes the selected entry.
DEL ALL	<i>[Delete all]</i>	Deletes all entries.
STANDBY	<i>[Standby]</i>	The steam generator changes to standby mode.

7.3 Menu bar interface pro

Menu sym- bol	Name	Explanation
	Home menu	<ul style="list-style-type: none"> • Process control parameters

Menu sym- bol	Name	Explanation
	<i>Manual control menu</i>	<ul style="list-style-type: none"> • Start various processes manually
	<i>Process menu</i>	<ul style="list-style-type: none"> • Editing and saving: <ul style="list-style-type: none"> ◦ Methods ◦ Series ◦ Templates ◦ Determinations
	<i>Configuration menu</i>	<ul style="list-style-type: none"> • Change settings • Service menu • System information
	<i>Determination data menu</i>	<ul style="list-style-type: none"> • View processed determinations

7.3.1 Home menu



Fig. 18: Home menu during a determination

- | | | | |
|---|--|---|---|
| 1 | Select and shows operation mode | 2 | Preview of next determination |
| 3 | Short-cut to determination data of the running determination | 4 | Status of the determination (incl. corresponding icons) |
| 5 | Remaining time | | |

7.3.2 Manual control menu

The manual control menu contains the following submenus:

Submenu	Explanation
Dose manually	See Chapter 8.13 "Dosing manually", page 72
Aspirate manually	See Chapter 8.14 "Aspirating manually", page 72
[Preheating]	See Chapter 8.1 "Editing the preheating function", page 50
[Priming]	See Chapter 8.2 "Editing the priming function", page 50
[Cleaning]	See Chapter 10.3 "Cleaning the glass components", page 79
[Stirring]	See Chapter 8.15 "Changing the stirring speed manually", page 73
[Pump Calibration]	See Chapter 10.14 "Calibrating the pumps", page 81

7.3.3 Method menu

The method menu contains the following submenus:

Submenu	Explanation
<i>[Single Determination]</i>	See Chapter 8.3 “Editing the single determination”, page 50
<i>[Series]</i>	See Chapter 8.4 “Editing a series”, page 53
<i>[Templates]</i>	See Chapter 8.5 “Editing a template”, page 57
<i>[Methods]</i>	See Chapter 8.6 “Editing a method”, page 58

7.3.4 Configuration menu

Settings submenu

Action	Option	Explanation
<i>[Language]</i>	Choice of display language on the interface	The following languages are available: English/German/French/Italian/ Spanish/Japanese/Chinese/Russian/ Polish
<i>[Current Date]</i>	Date input	Enter in sequence: Day, month, year. Apply the settings by pressing <i>[Save]</i> .
<i>[Current Time]</i>	Time input	Enter in sequence: Minutes, hours. Apply the settings by pressing <i>[Save]</i> .
Time zone	Choice of time zone	Select your time zone
<i>[Keyboard Layout]</i>	Choice of display keyboard on the interface	Select the keyboard layout
<i>[Key Tones]</i>	Off/On	Setting for audible signal in response to input controls.
<i>[Beep on Finish]</i>	Off/On	Setting for audible signal by the end of a determination.
<i>[Beep on Error]</i>	Off/On	Setting for audible signal by the end of a determination.
<i>[Display Brightness]</i>	Enter setting	Display illumination level in %: 0 - 100
<i>[Zero Amount Warning]</i>	On/Off	The instrument gives a warning when the entry for the weight is zero.
<i>[Default Amount Unit]</i>	Select value	The following values are selectable: g, mL
<i>[MaxAccuracy Mode]</i>	On / Off	Compensation of the steam power caused by voltage fluctuations.
<i>[AutoDist Mode]</i>	On/Off	Conditioning and distillation are performed automatically.
<i>[Level Detect. Max. Time]</i>	Enter value	The time after which the distillation stops without level detection sensor triggering.

Action	Option	Explanation
[Pump Maintenance Interval]	Select value	Selecting a frequency to carry out a pump rinsing. See Chapter 10.21 “Rinsing a pump”, page 86
[H ₂ O Dosage per Keypress]	Enter value	Dosage volume which is applied by pressing the related button.
[NaOH Dosage per Keypress]	Enter value	Dosage volume which is applied by pressing the related button.
[Acid Dosage per Keypress]	Enter value	Dosage volume which is applied by pressing the related button.
[H ₃ BO ₃ Dosage per Keypress]	Enter value	Dosage volume which is applied by pressing the related button.
[Aspiration Time Receiver]	Select value	Aspiration time for the related button.
[Aspiration Time Sample Tube]	Select value	Aspiration time for the related button.
[Steam Generator Standby]	Enter value	Enter the time until the instrument turns automatically into standby.
[Demo Mode]	On/Off	Simulate a distillation.
[Service Data Logger]	On/Off	Saves all data on the instrument for 14 months.
[Export Device]	Select value	Select between the following data export devices: USB stick/SD card

Peripherals submenu

Shows information about the connected peripherals.

User administration submenu

Creating and setting users. See Chapter 8.7 “Editing user setting”, page 68

Network submenu

Action	Option	Explanation
[Network]	Enter value	The following parameters can be edited: Device name/DHCP/MAC address/ System IP address/Subnet mask/ Gateway

Service submenu



NOTE

Only super user can carry out changes in the [Service] submenu.

System Information submenu














Technical information about the connected devices (e.g. serial number, firmware version).

7.3.5 Determination data menu

Shows the history of the determination data.

7.4 Status bar interface pro

The status bar shows actual information about the instrument.

View	Status
	The instrument is in standby mode.
	Process is running.
	The instrument is heating up.
	The BLE dongle is connected to the instrument.
	A level detection sensor is connected to the instrument.
	Reaction detection sensor attached.
	A tank level sensor is connected to the instrument.
	Dosing unit attached.
	Titration unit attached.
	White: A SD card is in the instrument. Red: The SD card in the instrument is in read only mode.
	A USB input device is connected.
	A USB memory stick is connected to the instrument.
	A recirculating chiller is connected.

8 Preparations for a determination







8.1 Editing the preheating function



NOTE

If the AutoDist Mode is active, the preheating is automated.

The preheating function defines how long the instrument parts are conditioned with steam.

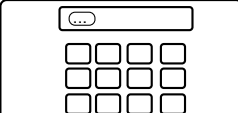	Specifications on the display:
▶ Navigate to the specified submenu.	 ↓
	
▶ Select the preheating function for which you want to change the preheating time.	
▶ Tap the specified function on the function bar.	
▶ Enter the preheating time in seconds.	
▶ Tap the specified function on the function bar. ⇒ The time is saved.	




8.2 Editing the priming function



NOTE

The method selected in this submenu is used for preparation.

The priming function puts the instrument in the ready to use status for a determination.

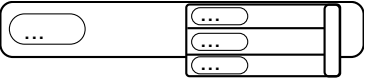	Specifications on the display:
▶ Navigate to the specified submenu.	 ↓
	
▶ Select the method you wish to use. ⇒ The priming method is selected.	






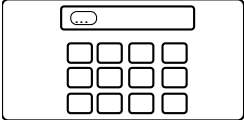

8.3 Editing the single determination

8.3.1 Changing the name of the single determination



NOTE


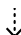
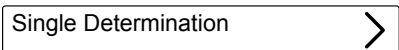


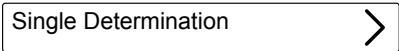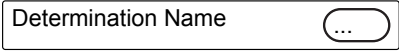

The name of the single determination is displayed in the *[Determination data]* menu.

Specifications on the display:	
▶ Navigate to the specified submenu.	 
	
▶ Select the specified action.	
▶ Tap the specified function on the function bar.	
▶ Enter the name for the single determination.	
▶ Tap the specified function on the function bar.	
⇒ The name for the determination is saved.	






8.3.2 Selecting the determination type for a single determination

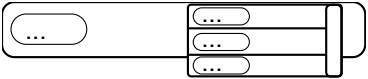
The following determination types are selectable:

Determination type	Explanation
<i>[Blank]</i>	For determinations without sample.
<i>[Reference Substance]</i>	For determinations with a defined reference substance.
<i>[Sample]</i>	For determinations with unknown sample.


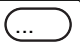
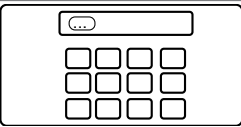
Specifications on the display:	
▶ Navigate to the specified submenu.	 
	
▶ Select the specified action.	
▶ Tap the specified function on the function bar.	
▶ Select the determination type you wish to use.	
⇒ The determination type is selected.	
▶ Tap the specified function on the function bar.	

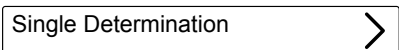8.3.3 Selecting the method for a single determination

Specifications on the display:	
▶ Navigate to the specified submenu.	 
	
▶ Select the specified action.	
▶ Tap the specified function on the function bar.	

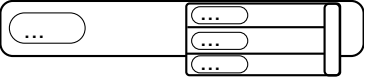Specifications on the display:	
▶ Select the method you wish to use. ⇒ The method is selected.	
▶ Tap the specified function on the function bar.	SAVE


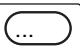
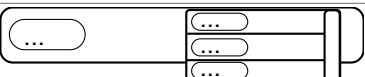
8.3.4 Changing the sample quantity for single determination

Specifications on the display:	
Precondition: ☑ The determination type is set to <i>[Sample]</i> . See Chapter 8.3.2 “Selecting the determination type for a single determination”, page 51	 ↓
▶ Navigate to the specified submenu.	Single Determination >
▶ Select the specified action.	
▶ Tap the specified function on the function bar.	EDIT
▶ Enter a value for the sample weight.	
▶ Tap the specified function on the function bar. ⇒ The sample weight is saved.	SAVE

8.3.5 Changing the unit for a single determination

The following units are available:

Unit	Explanation
<i>[g]</i>	Enter the weight for the determination in gram.
<i>[ml]</i>	Enter the weight for the determination in mL.

Specifications on the display:	
Precondition: ☑ The determination type is set to <i>[Sample]</i> or <i>[Reference Substance]</i> . See Chapter 8.3.2 “Selecting the determination type for a single determination”, page 51	 ↓
▶ Navigate to the specified submenu.	Single Determination >
▶ Select the specified action.	Unit 
▶ Tap the specified function on the function bar.	EDIT
▶ Select the unit you wish to use. ⇒ The unit is saved.	
▶ Tap the specified function on the function bar.	SAVE


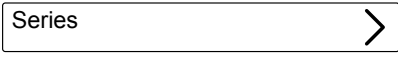

8.4 Editing a series

8.4.1 Creating a new series





There are two possibilities to create new series:

- Chapter “Creating a new series”, page 53
- Chapter “Creating a new series from a template”, page 53






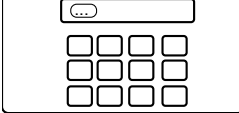

Creating a new series

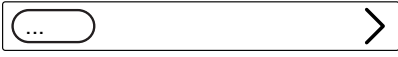	Specifications on the display:
▶ Navigate to the specified submenu.	 ↓ 
▶ Tap the specified function on the function bar. ⇒ The new series is created.	

Creating a new series from a template





	Specifications on the display:
▶ Navigate to the specified submenu.	 ↓ 
▶ Select the template from which you want to create a series.	
▶ Tap the specified function on the function bar. ⇒ The series is created.	

8.4.2 Changing the name of a series




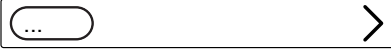

	Specifications on the display:
▶ Navigate to the specified submenu.	 ↓ 
▶ Select the name of the series that you wish to edit.	
▶ Select the specified action.	
▶ Tap the specified function on the function bar.	
▶ Enter a name for the series.	
▶ Tap the specified function on the function bar. ⇒ The name is saved.	

8.4.3 Adding a determination to a series






Add the first determination to a series

	Specifications on the display:
▶ Navigate to the specified submenu.	
	↓
	
▶ Select the series you wish to edit.	
▶ Tap the specified function on the function bar. ⇒ The determination is created.	


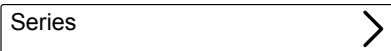
Add more determinations to a series by copy existing




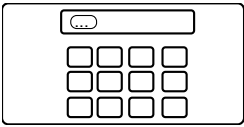
	Specifications on the display:
▶ Navigate to the specified submenu.	
	↓
	
▶ Select the series you wish to edit.	
▶ Select the determination you wish to copy.	
▶ Tap the specified function on the function bar. ⇒ The determination is created.	

8.4.4 Deleting a determination from a series

	Specifications on the display:
▶ Navigate to the specified submenu.	
	↓
	
▶ Select the series you wish to edit.	
▶ Select the determination you wish to delete.	
▶ Tap the specified function on the function bar. ⇒ The determination is deleted.	

8.4.5 Changing the determination name for a series







	Specifications on the display:
▶ Navigate to the specified submenu.	
	↓
	

Specifications on the display:	
▶ Select the series you wish to edit.	
▶ Select the determination you wish to edit.	
▶ Select the specified action.	Determination Name 
▶ Tap the specified function on the function bar.	EDIT
▶ Enter the name for the single determination.	
▶ Tap the specified function on the function bar. ⇒ The name for the determination is saved.	SAVE


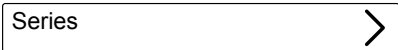
8.4.6 Selecting the determination type within a series







The following determination types are selectable:

Determination type	Explanation
<i>[Blank]</i>	For determinations without sample.
<i>[Reference Substance]</i>	For determinations with a defined reference substance.
<i>[Sample]</i>	For determinations with unknown sample.



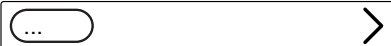





Specifications on the display:	
▶ Navigate to the specified submenu.	 ↓ 
▶ Select the series you wish to edit.	
▶ Select the determination you wish to edit.	
▶ Select the specified action.	Type 
▶ Tap the specified function on the function bar.	EDIT
▶ Select the determination type you wish to use. ⇒ The determination type is selected.	
▶ Tap the specified function on the function bar. ⇒ The name for the determination is saved.	SAVE

8.4.7 Selecting the method for a determination within a series

Specifications on the display:	
▶ Navigate to the specified submenu.	 ↓ 

	Specifications on the display:
▶ Select the series you wish to edit.	
▶ Select the determination you wish to edit.	
▶ Select the specified action.	Method Name 
▶ Tap the specified function on the function bar.	
▶ Select the determination type you wish to use. ⇒ The determination type is selected.	
▶ Tap the specified function on the function bar.	



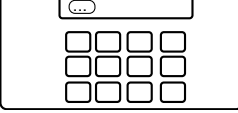





8.4.8 Changing the sample quantity for determination within a series

	Specifications on the display:
Precondition: <input checked="" type="checkbox"/> The determination type is set to <i>[Sample]</i> or <i>[Reference Substance]</i> . See Chapter 8.4.6 “Selecting the determination type within a series”, page 55	 ↓ Series 
▶ Navigate to the specified submenu.	
▶ Select the series you wish to edit.	
▶ Select the determination you wish to edit.	
▶ Select the specified action.	
▶ Tap the specified function on the function bar.	
▶ Enter a value for the sample weight.	
▶ Tap the specified function on the function bar. ⇒ The value for the sample weight is saved.	

8.4.9 Changing the amount unit for a determination within a series

The following units are available:

Unit	Explanation
<i>[g]</i>	Enter the weight for the determination in gram. Chapter 8.3.4 “Changing the sample quantity for single determination”, page 52
<i>[ml]</i>	Enter the weight for the determination in mL. See Chapter 8.3.4 “Changing the sample quantity for single determination”, page 52

	Specifications on the display:
Precondition: <input checked="" type="checkbox"/> The determination type is set to <i>[Sample]</i> or <i>[Reference Substance]</i> . See Chapter 8.4.6 "Selecting the determination type within a series", page 55	 ↓ 
▶ Navigate to the specified submenu.	
▶ Select the series you wish to edit.	
▶ Select the determination you wish to edit.	
▶ Select the specified action.	
▶ Select the determination type you wish to use. ⇒ The determination type is selected.	
▶ Tap the specified function on the function bar.	





8.5 Editing a template

8.5.1 Creating a new template




There are two possibilities to create a template:

- Chapter "Creating a new template", page 57
- Chapter "Creating a new template from an existing series", page 57




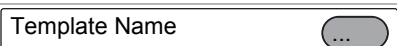

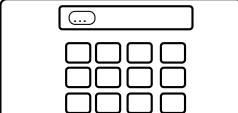

Creating a new template from an existing series

	Specifications on the display:
▶ Navigate to the specified submenu.	 ↓ 
▶ Select the series from which you want to create a template.	
▶ Tap the specified function on the function bar. ⇒ The template is created.	

Creating a new template


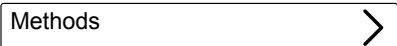


	Specifications on the display:
▶ Navigate to the specified submenu.	 ↓ 
▶ Tap the specified function on the function bar. ⇒ The template is created.	

8.5.2 Changing the name of a template

	Specifications on the display:
▶ Navigate to the specified submenu.	 ↓
	
▶ Select the name of the template that you wish to edit.	
▶ Select the specified action.	
▶ Tap the specified function on the function bar.	
▶ Enter a name for the template.	
▶ Tap the specified function on the function bar. ⇒ The template name is saved.	

8.6 Editing a method

8.6.1 Creating a new method





	Specifications on the display:
▶ Navigate to the specified submenu.	 ↓
	
▶ Select a method you wish to copy.	
▶ Tap the <i>[Copy]</i> function on the function bar. ⇒ A new method is created.	

8.6.2 Deleting a method






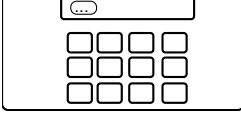


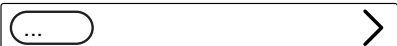
NOTE

Default methods *[Priming Method]* and *[Standard Method]* cannot be deleted.

	Specifications on the display:
▶ Navigate to the specified submenu.	 ↓
	
▶ Select a method you wish to delete.	
▶ Tap the specified function on the function bar.	
▶ Confirm the secure question. ⇒ The method is deleted.	




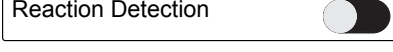

8.6.3 Changing the name of a method

	Specifications on the display:
▶ Navigate to the specified submenu.	 ↓
	
▶ Select the method that you wish to edit.	
▶ Select the specified action.	
▶ Tap the specified function on the function bar.	
▶ Enter a name for the method.	
▶ Tap the specified function on the function bar. ⇒ The name is saved.	

8.6.4 Changing the reaction detection for a method (option for MultiDist only)

The reaction detection sensor monitors the alkalization and doses the NaOH accordingly. The instrument switches automatically off when the alkalization is completed.

Status	Explanation
On	Automatic alkalization during the determination.
Off	Enter the parameters for the alkalization manually.

	Specifications on the display:
▶ Navigate to the specified submenu.	 ↓
	
▶ Select the method that you wish to edit.	
▶ Select the specified action.	
▶ Tap the specified function on the function bar. ⇒ The status is changed.	

8.6.5 Changing the H₂O volume for a method (MultiDist only)






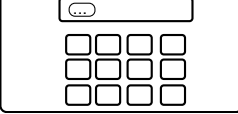



NOTE

The *[Reaction Detection]* action is set to Off. See Chapter 8.6.4 “Changing the reaction detection for a method (option for MultiDist only)”, page 59

Manually entering the quantity of H₂O to dilute the sample.

The input values can be found in the corresponding *Application note*.

Specifications on the display:	
▶ Navigate to the specified submenu.	
	↓
	
▶ Select the method that you wish to edit.	
▶ Select the specified action.	
▶ Tap the specified function on the function bar.	
▶ Enter the volume quantity.	
▶ Tap the specified function on the function bar.	
⇒ The quantity for the dosing is saved.	




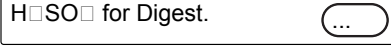

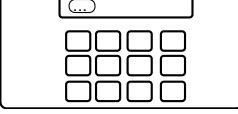

8.6.6 Changing the H₂SO₄ volume for a method (MultiDist only)



NOTE

The *[Reaction Detection]* action is set to On. See Chapter 8.6.4 “Changing the reaction detection for a method (option for MultiDist only)”, page 59

Enter manually the quantity of sulfuric acid.
The value is known from the previous digestion process.

Specifications on the display:	
▶ Navigate to the specified submenu.	
	↓
	
▶ Select the method that you wish to edit.	
▶ Select the specified action.	
▶ Tap the specified function on the function bar.	
▶ Enter the quantity of the sulfuric acid.	
▶ Tap the specified function on the function bar.	
⇒ The quantity for the sulfuric acid is saved.	






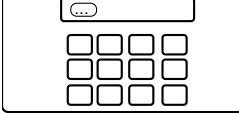

8.6.7 Changing the NaOH volume for a method (MultiDist only)



NOTE




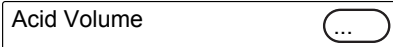

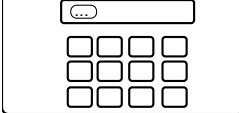

The *[Reaction Detection]* action is set to Off. See Chapter 8.6.4 “Changing the reaction detection for a method (option for MultiDist only)”, page 59

Enter manually the quantity of NaOH for alkalizing the sample.
For the correct quantity check related *Application note* or use the KjelOptimizer App.

	Specifications on the display:
▶ Navigate to the specified submenu.	 ↓
	
▶ Select the method that you wish to edit.	
▶ Select the specified action.	
▶ Tap the specified function on the function bar.	
▶ Enter the quantity for the dosing.	
▶ Tap the specified function on the function bar. ⇒ The quantity for the dosing is saved.	

8.6.8 Changing the acid volume for a method (BasicDist, MultiDist only)

Enter manually the quantity of reagent to be dosed.

	Specifications on the display:
▶ Navigate to the specified submenu.	 ↓
	
▶ Select the method that you wish to edit.	
▶ Select the specified action.	
▶ Tap the specified function on the function bar.	
▶ Enter the quantity for the dosing.	
▶ Tap the specified function on the function bar. ⇒ The quantity for the dosing is saved.	






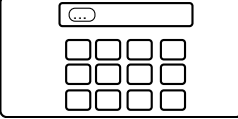

8.6.9 Changing the reaction time for a method (BasicDist, MultiDist only)



NOTE

The *[Reaction Detection]* action is set to Off. See Chapter 8.6.4 “Changing the reaction detection for a method (option for MultiDist only)”, page 59








The reaction time is the time between the addition of the NaOH and the start of the distillation.

Specifications on the display:	
▶ Navigate to the specified submenu.	 ↓
	
▶ Select the method that you wish to edit.	
▶ Select the specified action.	
▶ Tap the specified function on the function bar.	
▶ Enter a value for the reaction time.	
▶ Tap the specified function on the function bar.	
⇒ The reaction time is saved.	

8.6.10 Changing the number of steam steps for a method

Select steps to get the steam power gradually.

Selection	Explanation
[None]	Steam power in % selectable. See Chapter 8.6.12 “Changing the steam power for a method”, page 63
[2]	Steam power in % selectable. See Chapter 8.6.12 “Changing the steam power for a method”, page 63 Duration time of the steam step selectable. See Chapter 8.6.11 “Changing the time a steam step is carried out for a method”, page 63
[3]	Steam power in % selectable. See Chapter 8.6.12 “Changing the steam power for a method”, page 63 and Duration time of the steam step selectable. See Chapter 8.6.11 “Changing the time a steam step is carried out for a method”, page 63

Specifications on the display:	
▶ Navigate to the specified submenu.	 ↓
	
▶ Select the method that you wish to edit.	
▶ Select the specified action.	
▶ Tap the specified function on the function bar.	
▶ Select the number of steps you wish to use.	
▶ Tap the specified function on the function bar.	

8.6.11 Changing the time a steam step is carried out for a method



NOTE

Calculating the starting time for a steam step

- ☑ Make sure that the distillation time is set longer than the calculated time for the steam steps. See Chapter 8.6.14 “Changing the distillation time for a method”, page 64
- ▶ Starting time for steam step two:
 - ⇒ The duration time of the first steam step.
- ▶ Starting time for steam step three:
 - ⇒ The duration time of the first steam step.
 - ⇒ +
 - ⇒ The duration time of the second steam step.

	Specifications on the display:
Precondition:	
☑ More than one steam step is selected. See Chapter 8.6.10 “Changing the number of steam steps for a method”, page 62	↓
▶ Navigate to the specified submenu.	Methods >
▶ Select the method that you wish to edit.	... >
▶ Select the specified action.	Calibration NaOH ...
▶ Tap the specified function on the function bar.	EDIT
▶ Enter the time for the steam step.	
▶ Tap the specified function on the function bar.	SAVE
⇒ The time for the steam step is saved.	

8.6.12 Changing the steam power for a method

The steam power in % during a steam step.

	Specifications on the display:
▶ Navigate to the specified submenu.	
	↓
	Methods >
▶ Select the method that you wish to edit.	... >
▶ Select the specified action.	Steam Power ...
▶ Tap the specified function on the function bar.	EDIT
▶ Enter the level for the steam power.	

Specifications on the display:

- ▶ Tap the specified function on the function bar. SAVE
- ⇒ The level for the steam power is saved.

8.6.13 Changing the level detection sensor settings (option) for a method

The level detection stops the distillation when a defined distillate volume is reached.



NOTE

The maximum level detection time can be changed in the *Settings* submenu.

Selection	Explanation
On	The OnLevel sensor is active.
Off	The OnLevel sensor is off.

Specifications on the display:

- ▶ Navigate to the specified submenu. ✎
↓
- Methods >
- ▶ Select the method that you wish to edit. ... >
- ▶ Select the specified action. Level Detection
- ▶ Tap the specified function on the function bar. EDIT
- ⇒ The status changed.

8.6.14 Changing the distillation time for a method






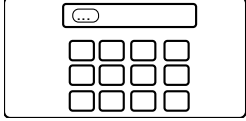

Enter manually the time for the determination.

Specifications on the display:

- Precondition:
- The *[Level Detection]* action is set to Off. See Chapter 8.6.13 “Changing the level detection sensor settings (option) for a method”, page 64
- ▶ Navigate to the specified submenu. ✎
↓
 - Methods >
 - ▶ Select the method that you wish to edit. ... >
 - ▶ Select the specified action. Distillation Time
 - ▶ Tap the specified function on the function bar. EDIT
 - ▶ Enter a value for the distillation time.





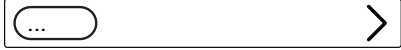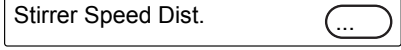


0	1	2	3
4	5	6	7
8	9	.	=
 - ▶ Tap the specified function on the function bar. SAVE
 - ⇒ The distillation time is saved.

8.6.15 Changing the distillation stirring speed for a method (MultiDist only)

	Specifications on the display:
▶ Navigate to the specified submenu.	 ↓
	
▶ Select the method that you wish to edit.	
▶ Select the specified action.	
▶ Tap the specified function on the function bar.	
▶ Enter the stirring speed.	
▶ Tap the specified function on the function bar.	
⇒ The stirring speed is saved.	

8.6.16 Changing the titration type for a method (MultiDist option only)

Titration type	Explanation
[None]	No titration
[Boric Acid Titration]	Titration
[Back Titration]	Back titration






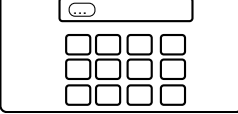

	Specifications on the display:
▶ Navigate to the specified submenu.	 ↓
	
▶ Select the method that you wish to edit.	
▶ Select the specified action.	
▶ Tap the specified function on the function bar.	
▶ Select the titration type you wish to use.	
▶ Tap the specified function on the function bar.	

8.6.17 Changing the H₃BO₃ volume for a method (MultiDist only)



NOTE

The availability depends on the selected titration type.


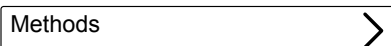

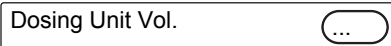

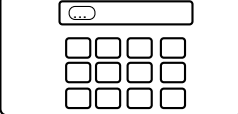

Specifications on the display:	
▶ Navigate to the specified submenu.	 ↓
	
▶ Select the method that you wish to edit.	
▶ Select the specified action.	
▶ Tap the specified function on the function bar.	
▶ Enter the quantity for the volume.	
▶ Tap the specified function on the function bar. ⇒ The volume is saved.	

8.6.18 Changing the dosing unit volume for a method (MultiDist option only)



NOTE

The availability depends on the selected titration type.





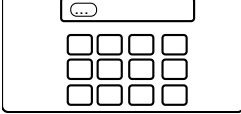

Specifications on the display:	
▶ Navigate to the specified submenu.	 ↓
	
▶ Select the method that you wish to edit.	
▶ Select the specified action.	
▶ Tap the specified function on the function bar.	
▶ Enter the quantity for the volume.	
▶ Tap the specified function on the function bar. ⇒ The volume is saved.	

8.6.19 Changing the titration stirring speed for a method (MultiDist option only)



NOTE

The availability depends on the selected titration type.

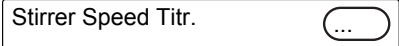
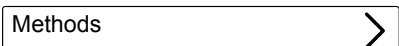
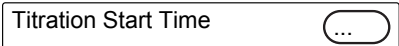

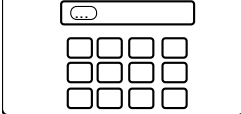

	Specifications on the display:
Precondition: <input checked="" type="checkbox"/> A titration type is activated. See ▶ Navigate to the specified submenu.	 ↓
▶ Select the method that you wish to edit.	
▶ Select the specified action.	
▶ Tap the specified function on the function bar.	
▶ Enter the stirring speed.	
▶ Tap the specified function on the function bar. ⇒ The stirring speed is saved.	

8.6.20 Changing the titration start time for a method (MultiDist option only)





NOTE

The availability depends on the selected titration type.

	Specifications on the display:
▶ Navigate to the specified submenu.	 ↓
▶ Select the method that you wish to edit.	
▶ Select the specified action.	
▶ Tap the specified function on the function bar.	
▶ Enter a value for the titration start time.	
▶ Tap the specified function on the function bar. ⇒ The start time is saved.	

8.6.21 Changing the aspiration time for the sample tube for a method (BasicDist, MultiDist only)

	Specifications on the display:
▶ Navigate to the specified submenu.	 ↓
	

Specifications on the display:	
▶ Select the method that you wish to edit.	
▶ Select the specified action.	Sample Tube Aspiration
▶ Tap the specified function on the function bar.	EDIT
▶ Enter a value for the aspiration time.	
▶ Tap the specified function on the function bar. ⇒ The aspiration time is saved.	SAVE

8.6.22 Changing the aspiration time for the receiving vessel for a method (MultiDist only)

Specifications on the display:	
▶ Navigate to the specified submenu.	 ↓ Methods
▶ Select the method that you wish to edit.	
▶ Select the specified action.	Receiver Aspiration
▶ Tap the specified function on the function bar.	EDIT
▶ Enter a value for the aspiration time.	
▶ Tap the specified function on the function bar. ⇒ The aspiration time is saved.	SAVE

8.7 Editing user setting



NOTE

The user *[Service User]* is a default user and password protected.

8.7.1 Creating a new user

Specifications on the display:	
▶ Navigate to the specified submenu.	 ↓ User Administration
▶ Tap the specified function on the function bar.	NEW
▶ Enter a name for the user.	

Specifications on the display:


▶ Tap the specified function on the function bar.



8.7.2 Deleting a user

Specifications on the display:

▶ Navigate to the specified submenu.

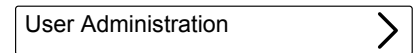

▶ Select the user you wish to delete.



▶ Tap the specified function on the function bar.





▶ Confirm the secure question.

⇒ The user is deleted.

8.7.3 Changing the name of a user

Specifications on the display:

▶ Navigate to the specified submenu.

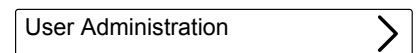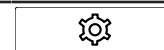


▶ Select the user that you wish to edit.



▶ Tap the specified function on the function bar.



▶ Enter a new name for the user.

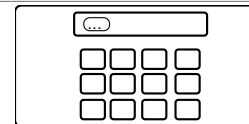

▶ Tap the specified function on the function bar.




⇒ The new name is saved.

8.7.4 Selecting a user

Specifications on the display:

▶ Navigate to the specified submenu.

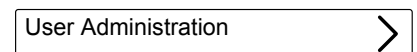

▶ Select the user you wish to use.



▶ Tap the specified function on the function bar.



⇒ The user is active.

⇒ The user is marked up green.

8.8 Installing the sample tube




NOTE


Removing is done in reverse sequence.

Precondition:


- ☑ Make sure, that the end stop is in the correct position for the sample tube in use. See Chapter 8.10 “Preparing the end stop for the sample tube size”, page 71
- ▶ Hold the sample tube with the tongs.

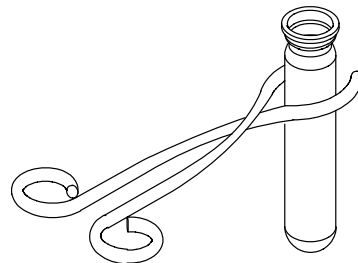


- ▶ Insert the sample tube.

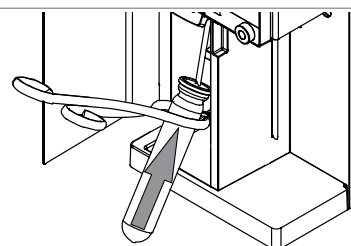

- ▶ Apply a light force to push the sample tube against the end stop and the sample tube sealing.

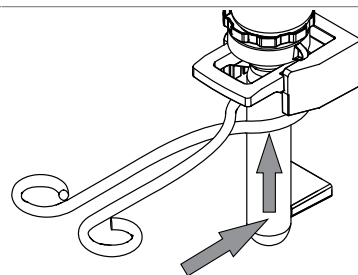


- ▶ Secure the sample tube in place with the handle.

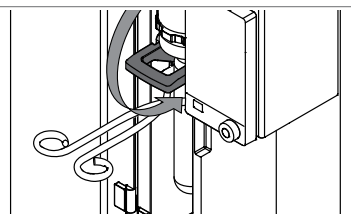


8.9 Installing a receiving vessel

- ▶ Insert the sample receiving vessel to the instrument.

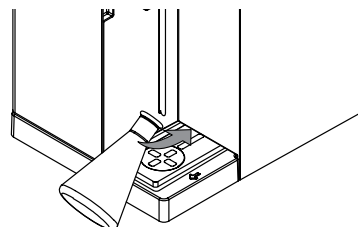


- ▶ Place the sample receiving vessel on the receiving vessel area.

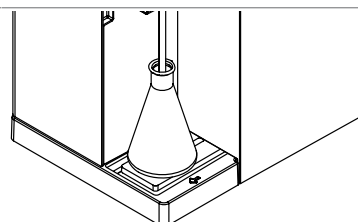

8.10 Preparing the end stop for the sample tube size


- ▶ Pull the end stop.
- ▶ Twist the end stop.

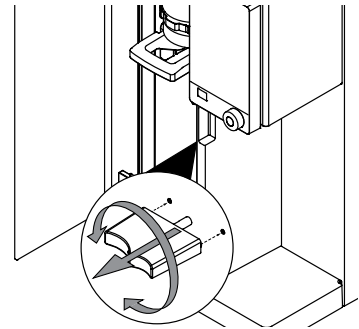


8.11 Preparing the OnLevel sensor (option)

Precondition:

- The installations for the OnLevel sensor are completed. See Chapter 5.18 “Installing the OnLevel sensor (option)”, page 37
- ▶ Attach the magnetic sensor to the metal plate.

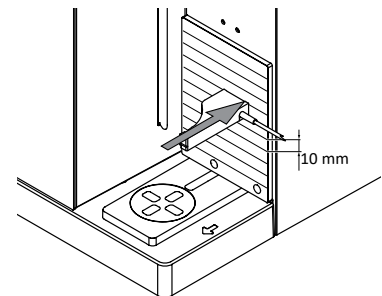

- ▶ Adjust the sensor near the sample receiving vessel.

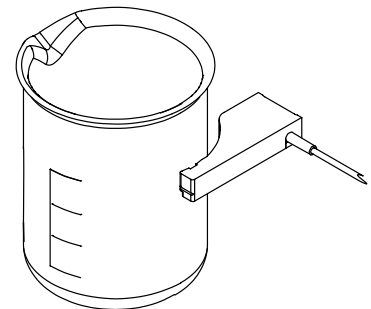

8.12 Selecting the operation mode

Select one of the following operation modes:

Operation mode	Explanation
[Single Determination]	For carrying out one determination. Editing settings see Chapter 8.3 “Editing the single determination”, page 50
[Series]	For planning and carrying out a series of determinations. Editing settings see Chapter 8.4 “Editing a series”, page 53

Specifications on the display:

▶ Navigate to the specified menu.	
▶ Select the specified action.	Operation Mode
▶ Tap the specified function on the function bar.	
▶ Select the operation mode according to your requirements.	
▶ Tap the specified function on the function bar.	

8.13 Dosing manually

8.13.1 Dosing H₂O manually



NOTE

The quantity of liquid per keystroke can be changed in the *[Settings]* submenu.

Specifications on the display:

Precondition:

- A sample tube is installed. See Chapter 8.8
"Installing the sample tube", page 70

▶ Navigate to the specified menu.

▶ Tap the specified button.

⇒ The pump applies a defined quantity of H₂O.



8.13.2 Dosing NaOH manually



NOTE

The quantity of liquid per keystroke can be changed in the *[Settings]* submenu.

Specifications on the display:

Precondition:

- A sample tube is installed. See Chapter 8.8
"Installing the sample tube", page 70

▶ Navigate to the specified menu.

▶ Tap the specified button.

⇒ The pump applies a defined quantity of NaOH.



8.13.3 Dosing H₃BO₃ manually



NOTE

The quantity of liquid per keystroke can be changed in the *[Settings]* submenu.

Specifications on the display:

Precondition:

- A receiving vessel is installed. See Chapter 8.9
"Installing a receiving vessel", page 70

▶ Navigate to the specified menu.

▶ Tap the specified button.

⇒ The pump applies a defined quantity of H₃BO₃.



8.14 Aspirating manually

8.14.1 Aspirating the sample tube manually



NOTE

The aspiration time per keystroke can be changed in the *Settings* submenu.

Specifications on the display:

▶ Navigate to the specified menu.



Specifications on the display:

- ▶ Tap the specified button. [Sample Tube Aspiration]
- ⇒ The pump aspirates liquid from the sample tube.

8.14.2 Aspirating the receiving vessel manually



NOTE

The aspiration time per keystroke can be changed in the *Settings* submenu.

Specifications on the display:

- ▶ Navigate to the specified menu.
- ▶ Tap the specified button. [Receiver Aspiration]
- ⇒ The pump aspirates liquid from the receiving vessel.

8.15 Changing the stirring speed manually

Specifications on the display:

- ▶ Navigate to the specified submenu.
↓
Stirring >
- ▶ Select the specified action. Stirrer Speed ...
- ▶ Tap the specified function on the function bar. EDIT
- ▶ Enter the stirring speed.
- ▶ Tap the specified function on the function bar. SAVE
- ⇒ The stirring speed is saved.

8.16 Switching valve adjustment (MultiDist only)



NOTE


After switching the valve, the instrument automatically rinses the pumps.

The following valve positions are available:

Valve position	Explanation
<i>Acid</i>	The system dose acid to the sample tube.
<i>NaOH</i>	The system dose NaOH to the sample tube.

Precondition:

- The On /Off master switch is set to on.
- ▶ Rinse the pump you want to use. See Chapter 10.21 “Rinsing a pump”, page 86
- ▶ Turn the valve in the correct position.
- ▶ Follow the instructions on the interface.

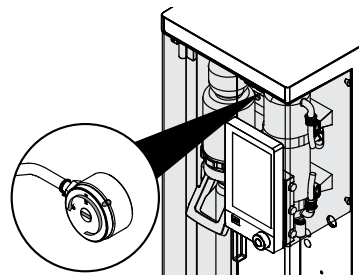

9 Carrying out a determination

9.1 Preparing the instrument

Navigation path



Precondition:

- All commissioning operations have been completed. See Chapter 5 “Installation”, page 29
- ▶ Set the On/Off master switch to On.
 - ⇒ The instrument is starting up.
- ▶ Make sure that no defective sealings or glass parts are used.
- ▶ Navigate to the *Home* menu according the navigation path.
- ▶ Tap the **[READY]** button on the function bar.
 - ⇒ The instrument is heating up.
- ▶ Install a receiving vessel (EasyDist and BasicDist only). See Chapter 8.9 “Installing a receiving vessel”, page 70
- ▶ Install the sample tube. See Chapter 8.8 “Installing the sample tube”, page 70
- ▶ Close the protection shield.
- ▶ Flush the tubing. See Chapter 8.13 “Dosing manually”, page 72
- ▶ Aspirate the receiving and the sample vessels. See Chapter 8.14 “Aspirating manually”, page 72

9.2 Starting a determination

Navigation path



Precondition:

- The instrument is prepared. See Chapter 9.1 “Preparing the instrument”, page 75
- The sample is prepared.
- The connected instruments are prepared (MultiDist only). See related documentation.
- ▶ Navigate to the *Home* menu according the navigation path.
- ▶ Press the **[PREP]** button on the function bar.
 - ⇒ Wait until the status bar shows the status **READY**.
- ▶ Select the operation mode. See Chapter 8.12 “Selecting the operation mode”, page 71
- ▶ Tap the **[START]** button on the function bar.

9.3 Ending a determination

Navigation path



Precondition:


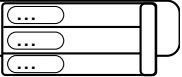
- The display shows the *Finished* dialog.
- ▶ Confirm the message on the dialog.

9.4 Shutting down the instrument


Precondition:

- ☑ The distillation process has ended.
- ▶ Shut down the connected instruments (MultiDist only). See related documentation.
- ▶ Clean the instrument. See Chapter 10.3 “Cleaning the glass components”, page 79
- ▶ Install an empty sample tube. See Chapter 8.8 “Installing the sample tube”, page 70
- ▶ Install an empty receiving vessel (EasyDist and BasicDist only). See Chapter 8.9 “Installing a receiving vessel”, page 70
- ▶ Set the On/Off master switch to off.

9.5 Filtering determination data


	Specifications on the display:
▶ Navigate to the specified menu.	
▶ Select the specified action.	Filter by ...
▶ Select the filter you wish to use.	... 

9.6 Showing determination data


▶ Navigate to the specified submenu.	
▶ Select the determination you wish to view. ⇒ The display shows the data of the determination.	... >

9.7 Deleting determination data

9.7.1 Deleting one determination




	Specifications on the display:
▶ Navigate to the specified menu.	
▶ Select the determination you wish to delete.	... >
▶ Tap the specified function on the function bar.	DELETE
▶ Confirm the secure question. ⇒ The determination data is deleted.	

9.7.2 Deleting all determination data



	Specifications on the display:
▶ Navigate to the specified menu.	
▶ Tap the specified function on the function bar.	DEL ALL
▶ Confirm the secure question. ⇒ All determination data are deleted.	

9.8 Exporting determination data (Interface pro only)

9.8.1 Exporting one determination (Interface pro only)

Specifications on the display:	
Precondition:	
<input checked="" type="checkbox"/> A storage medium is available. See Chapter "Settings submenu", page 47	
▶ Navigate to the specified menu.	
▶ Select the determination you wish to export.	
▶ Tap the specified function on the function bar.	

9.8.2 Exporting all determination data (Interface pro only)

Specifications on the display:	
▶ Navigate to the specified menu.	
▶ Tap the specified function on the function bar.	

10 Cleaning and servicing



NOTE

Users may only carry out the servicing and cleaning operations described in this section.

Any servicing and repair work which involves opening up the casing may only be carried out by BUCHI service technicians.

- ▶ Use only genuine BUCHI consumables and spare parts in order to ensure correct operation of the device and preserve the warranty.

10.1 Regular maintenance work

Action	Daily	Weekly	Monthly	Yearly	Additional information
10.2 Cleaning and servicing the sample tube	1				Carry out this action before every instrument use.
10.3 Cleaning the glass components	1				Carry out this action after every instrument use.
10.4 Cleaning and servicing the hoses and hose connections		2			
10.5 Cleaning and servicing the sample tube seal		2			Replace the sample tube seal twice per year
10.6 Cleaning and servicing the splash protector		2			Replace the glass splash protector after ~ 3500 distillations (twice per year). Replace the plastic splash protector once in 2 years.
10.7 Cleaning and servicing the steam inlet and condenser outlet tube		2			Replace the tube once per year.
10.8 Cleaning the housing		2			
10.9 Cleaning and servicing the warning and directive symbols		2			
10.10 Cleaning and servicing the bridge splash protector to condenser		2			Replace the bridge once per year.
10.11 Cleaning and servicing the dosing pumps		2			Replace the NaOH pump once per year
10.15 Cleaning and servicing the waste pumps			1		
10.12 Cleaning and servicing the condenser			2		
10.13 Cleaning and servicing the steam generator			2		
10.14.1 Calibrating the H ₂ O pump			2		

Action	Daily	Weekly	Monthly	Yearly	Additional information
10.14.2 Calibrating the acid pump			2		
10.14.3 Calibrating the NaOH pump			2		
10.14.4 Calibrating the H ₃ BO ₃ Pump			2		
10.16 Decalcifying the instrument				2	

1 - User; 2 - Operator

10.2 Cleaning and servicing the sample tube

- ▶ Before using, check the sample tube for defects (cracks/splintering).
- ▶ Prevent temperature shocks on the sample tubes.
 - ⇒ Temperature shocks may lead to breakage.

Cleaning

If cleaning with a dishwashing machine:

- ▶ Make sure, that there is no contact with other glassware.





Storage (e.g. drawer)

- ▶ Prevent rolling and small hits.
 - ⇒ This can cause hair cracks which may lead to breakages.

10.3 Cleaning the glass components

The following values for the cleaning can be adjusted (the list differs according to the instrument configuration):

- [H₂O Volume]
- [Steam Power]
- [Distillation Time]
- [Sample Tube Aspiration]
- [Receiver Aspiration]

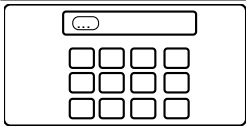	Specifications on the display:
Precondition:	
<input checked="" type="checkbox"/> A sample tube is installed. See Chapter 8.8 "Installing the sample tube", page 70	↓
<input checked="" type="checkbox"/> A sample receiving vessel is installed. See Chapter 8.9 "Installing a receiving vessel", page 70	
▶ Navigate to the specified submenu.	
▶ Adjust the values according to your needs.	
▶ Confirm the entries with the specified function in the function bar.	

Specifications on the display:

- ▶ Tap the specified function on the function bar.

START**10.4 Cleaning and servicing the hoses and hose connections**

- ▶ Check the hoses and hose connections for defects (cracks, brittle areas).
- ▶ If defective, replace the hoses.

10.5 Cleaning and servicing the sample tube seal

- ▶ Check the seal for the following damage that may cause leakage:
 - Deterioration (surface roughness)
 - Deformation (rubber shape)
 - Residues (dirt, white crystals)
- ▶ Rinse the seal with water.
- ▶ If necessary, replace the seal. See Chapter 10.22 “Replacing the sample tube seal”, page 87

10.6 Cleaning and servicing the splash protector

- ▶ Check for leaks on the connectors.
- ▶ Check for residues inside the splash protector.
- ▶ Check the glass wall of the main body and the connectors.
 - ⇒ If the part is dirty, clean it with detergents.
 - ⇒ If the glass part is showing deterioration (transparency loss/reams in the glass wall) or leakages (white residues), replace it.

10.7 Cleaning and servicing the steam inlet and condenser outlet tube

- ▶ Check the tubes concerning dirt or residues.
- ▶ Clean and replace periodically depending on the use.

10.8 Cleaning the housing

- ▶ Wipe down the housing with a damp cloth.
- ▶ If heavily soiled, use ethanol or a mild detergent.
- ▶ Wipe down the display with a damp cloth.

10.9 Cleaning and servicing the warning and directive symbols

- ▶ Check that the warning symbols on the instrument are legible.
- ▶ If they are dirty, clean them.

10.10 Cleaning and servicing the bridge splash protector to condenser

- ▶ Check if it is tight.
- ▶ Check for residues inside the tube.
- ▶ If necessary, clean it with a brush and water.

10.11 Cleaning and servicing the dosing pumps

- ▶ Check the pump connectors in the dosing pump window for:
 - Leakages
 - Crystallization
- ▶ If necessary, contact BUCHI customer service.

10.12 Cleaning and servicing the condenser

Decalcification solution:

160 g citric acid or 80 g amidosulfuric acid dissolved in 0.8 L water.






- ▶ Check if the cooling media of the condenser is transparent.
- ▶ Check if the inside of the coil is showing condensation droplets.
 - ⇒ If droplets appear:
 - Flush the condenser coil through the condensate expansion vessel with an 1% hydrochloric acid.
 - Flush the condenser coil through the condensate expansion vessel with water and water/ethanol mixture.
- ▶ Clean the condenser on the outside of the coil (water connectors) with decalcification solution to remove chalk, algae, and rust.

10.13 Cleaning and servicing the steam generator


- ▶ Check the distillation volume.
 - ⇒ If the condensate amount is less than 160 mL water/per 5 minutes distillation (after preheating), decalcify the instrument. See Chapter 10.16 “Decalcifying the instrument”, page 83

10.14 Calibrating the pumps

10.14.1 Calibrating the H₂O pump

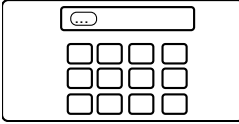	Specifications on the display:
Precondition:	
<input checked="" type="checkbox"/> A sample tube is installed. See Chapter 8.8 “Installing the sample tube”, page 70	⋮
▶ Navigate to the specified submenu.	Pump Calibration 
	⋮
	H ₂ O 
▶ Select the specified action.	Nominal Volume 
▶ Tap the specified function on the function bar.	EDIT
▶ Enter the calibration volume you wish to use.	
▶ Tap the specified function on the function bar.	START


Specifications on the display:

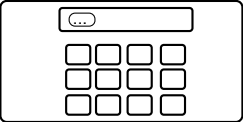- ▶ Wait until the display shows a numeric dialog. 

- ▶ Remove the sample tube.


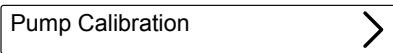
- ▶ Measure the volume in the sample tube.


- ▶ Enter the value in the numeric dialog.


- ▶ Tap the specified function on the function bar. 


10.14.2 Calibrating the acid pump

Specifications on the display:

- Precondition:
- A sample tube is installed. See Chapter 8.9 “Installing a receiving vessel”, page 70
- ▶ Navigate to the specified submenu.  ↓ 

 - ▶ Select the specified action. 


 - ▶ Tap the specified function on the function bar. 

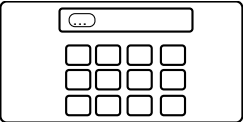 - ▶ Wait until the display shows a numeric dialog. 

 - ▶ Remove the sample tube.


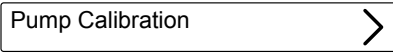
 - ▶ Measure the volume in the sample tube.


 - ▶ Enter the value in the numeric dialog.


 - ▶ Tap the specified function on the function bar. 


10.14.3 Calibrating the NaOH pump


Specifications on the display:


- Precondition:
- A sample tube is installed. See Chapter 8.9 “Installing a receiving vessel”, page 70
- ▶ Navigate to the specified submenu.  ↓ 

 - ▶ Select the specified action. 

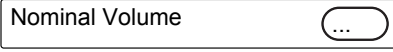 - ▶ Tap the specified function on the function bar. 


 - ▶ Tap the specified function on the function bar. 

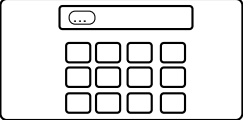 - ▶ Enter the calibration volume you wish to use. 

 - ▶ Tap the specified function on the function bar. 

Specifications on the display:

▶ Wait until the display shows a numeric dialog.

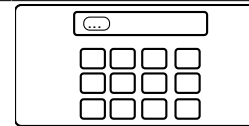

▶ Remove the sample tube.

▶ Measure the volume in the sample tube.

▶ Enter the value in the numeric dialog.

▶ Tap the specified function on the function bar.




10.14.4 Calibrating the H₃BO₃ Pump

Specifications on the display:

Precondition:

A receiving vessel is installed. See Chapter 8.9 “Installing a receiving vessel”, page 70

▶ Navigate to the specified submenu.

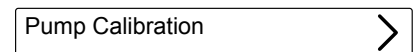


▶ Select the specified action.



▶ Tap the specified function on the function bar.




▶ Enter the calibration volume you wish to use.

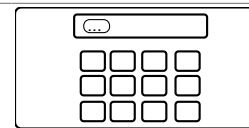

▶ Tap the specified function on the function bar.



▶ Wait until the display shows a numeric dialog.

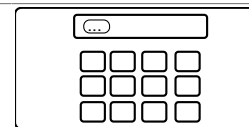

▶ Remove the sample receiving vessel.

▶ Measure the volume in the sample receiving vessel.

▶ Enter the measured volume.

▶ Tap the specified function on the function bar.



10.15 Cleaning and servicing the waste pumps

- ▶ Check for leaks on the connectors.
- ▶ Check the tubing to peristaltic pump.
- ▶ If necessary, replace the tubing to peristaltic pump.
- ▶ If necessary, contact BUCHI customer service for exchanging the pump.

10.16 Decalcifying the instrument

Time required: approximately 2.5 hours

Decalcification solution:

160 g citric acid or 80 g amidosulfuric acid dissolved in 0.8 L water.

Precondition:


- ☑ The instrument temperature is the same as the ambient temperature.
- ▶ Remove the water from the instrument. See Chapter 12.4 “Removing water from the steam generator”, page 93
- ▶ Prepare the decalcification solution.
- ▶ Install a suitable hose to the H₂O supply for steam generation.
- ▶ Put the other end of the hose in the decalcification solution.
- ▶ Set the On/Off master switch to On.
- ▶ Press the *[READY]* button on the function bar.
- ▶ Wait until the pumps stop working.
- ▶ Set the On/Off master switch to Off.
- ▶ Wait 30 minutes.
- ▶ Remove the decalcification solution from the instrument. See Chapter 12.4 “Removing water from the steam generator”, page 93
- ▶ Do steps (5) thru (10) again.
- ▶ Install the H₂O supply for steam generation.
- ▶ Set the On/Off master switch to On.
- ▶ Press the *[READY]* button on the function bar.
- ▶ Wait until the pumps stop working.
- ▶ Set the On/Off master switch to Off.
- ▶ Remove the water from the instrument. See Chapter 12.4 “Removing water from the steam generator”, page 93
- ▶ Do steps (11) thru (16) five times.

10.17 Replacing the splash protector


**NOTE**

Installing is done in reverse sequence.


- ▶ Remove the protection shield. See Chapter 10.19 “Attaching and removing the protection shield”, page 86
- ▶ Loosen the cap nut at the splash protector.

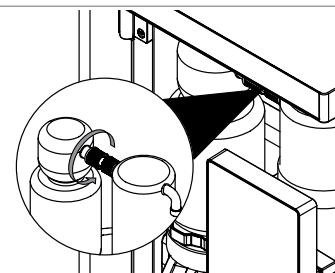


- ▶ Open the splash protector nut.

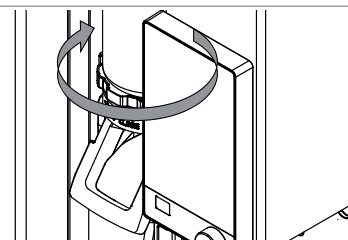

- ▶ Remove the splash protector.

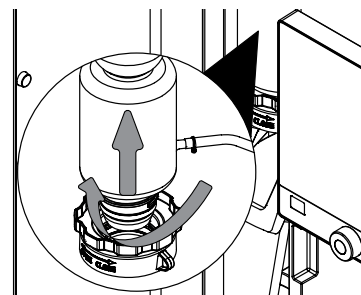

- ▶ Remove the steam connection.
- ▶ Remove the NaOH/acid connection (according to the instrument configuration).

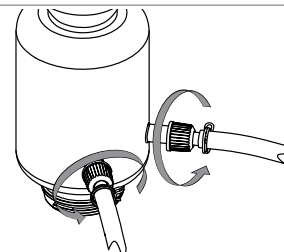

10.18 Replacing the condenser




NOTE


Installing is done in reverse sequence.

Precondition:


- A one liter beaker flask is available.
- ▶ Remove the protection shield. See Chapter 10.19 “Attaching and removing the protection shield”, page 86
- ▶ Loosen the cap nut at the condenser.

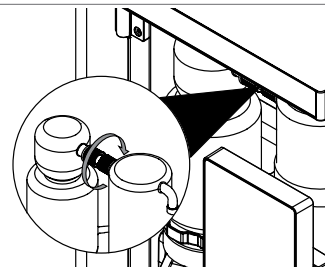


- ▶ Remove the lower cooling water hose from the instrument.

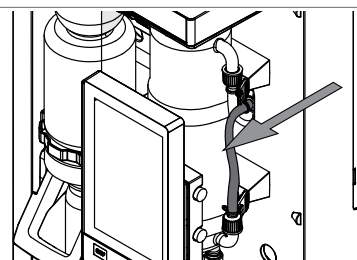

- ▶ Put the cooling water hose in the beaker flask.
- ▶ Loosen the upper cooling hose from the condenser.

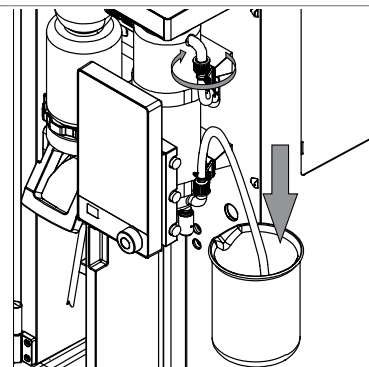


- ▶ Wait until the condenser is empty.
- ▶ Remove the rubber band.

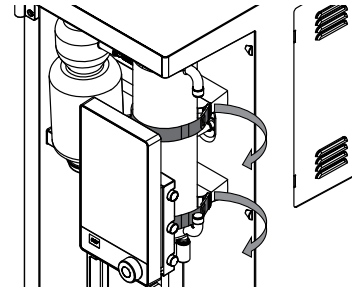

10.19 Attaching and removing the protection shield


10.19.1 Removing the protection shield

- ▶ Pull the protection shield from the instrument.

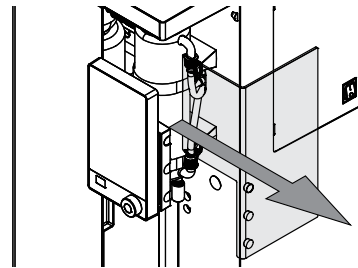


10.19.2 Attaching the protection shield

- ▶ Attach the protection shield onto the instrument.

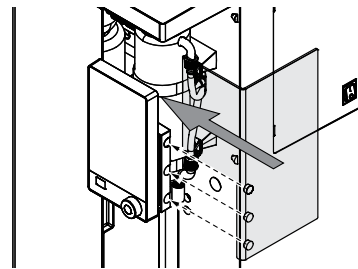

10.20 Cleaning the drip tray

- ▶ Pull the drip tray.
- ▶ Rinse the trip tray with water.

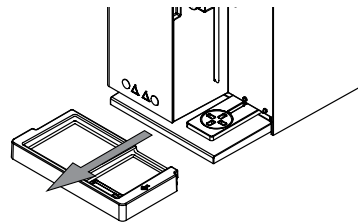

10.21 Rinsing a pump

Precondition:

- A sample tube is installed. See Chapter 8.8 “Installing the sample tube”, page 70
- ▶ Put the side of the hose in distilled water.
- ▶ Apply 100 mL with the manual functions. See Chapter 8.13 “Dosing manually”, page 72

10.22 Replacing the sample tube seal

**NOTE**


Installing is done in reverse sequence.

**NOTE**


Make the seal moist with water before installation.

- ▶ Remove the protection shield. See Chapter 10.19.1 “Removing the protection shield”, page 86


- ▶ Loosen the cap nut.

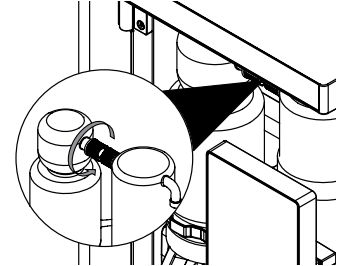



- ▶ Open the splash protector nut.

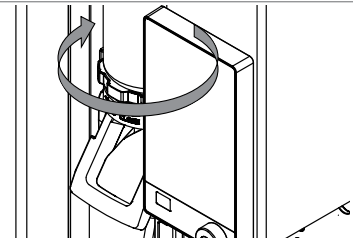

- ▶ Remove the splash protector.

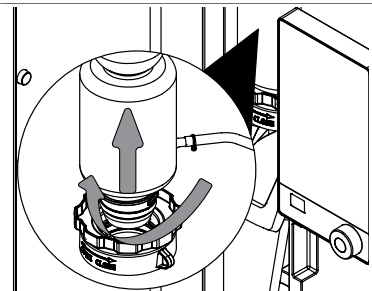

- ▶ Remove the seal.

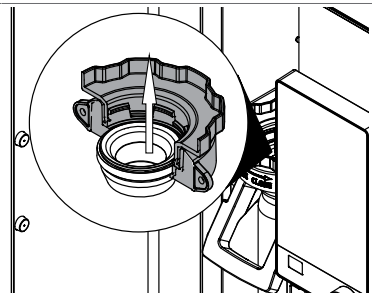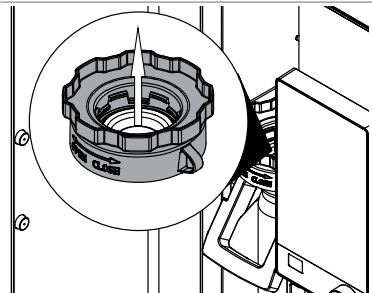

11 Help with faults

11.1 Troubleshooting digestion

Problem	Possible cause	Action
Crystallization after digestion	False ratio of H ₂ SO ₄ to catalyst.	▶ Use KjelOptimizer software to optimize the digestion (download available on BUCHI website).
	Digestion time too long.	▶ Increase digestion time. ▶ Compare to conditions to similar applications.
	Suction capacity of scrubber too strong.	▶ Lower the suction capacity on the scrubber. See <i>Scrubber</i> operation manual
Samples do not get clear	No or not enough catalyst used.	▶ Use KjelOptimizer software to optimize the digestion (download available on BUCHI website).
	Digestion temperature is too low.	▶ Increase digestion time. ▶ Compare to conditions to similar applications.
	Temperature is too high.	▶ Decrease digestion temperature.
	Sealing material was flushed into the sample.	▶ Check the sealing.
Fume leakage	The seals are defective.	▶ Replace the seals.
	Suction capacity of scrubber is too weak.	▶ Carry out the scrubber maintenance. See <i>Scrubber</i> operation manual
	Leakage in the system, e.g. hose connector not tight.	▶ Check connections between digester and scrubber.
	Blocked hoses.	▶ Clean the hoses.
	Reduced suction on the bypass valve.	
Boiling retardation/ bumping/foaming	Missing digestion rods or boiling stones.	▶ Add boiling rods during digestion.
	Missing antifoam tablet or other anti-foaming agent.	▶ Add antifoam tablet.
Samples do not get dark blue/brown after addition of NaOH	Air in NaOH hose.	▶ Flush the system. See Chapter 8.13 "Dosing manually", page 72
	No catalyst used for digestion (only H ₂ O ₂ or Kjeldahl tablets ECO).	

11.2 Troubleshooting distillation

Problem	Possible cause	Action
Splashing during distillation or addition of chemicals	The sample tube is too small.	▶ Select a bigger sample tube.
	Sample volume too high.	▶ Lower sample volume.
	Amount of water used for dilution is too low.	▶ Increase dilution volume.
Measured nitrogen content is too high	Air in titration system, burette or tubes.	▶ Tighten the connections. ▶ Flush the tubing with titrant. ▶ Refill burette.
	Carry over during distillation.	▶ Insert less sample. ▶ Avoid big concentration differences in measurement series.
	Titrant concentration too high.	▶ Lower the titrant concentration.
	Error in calculation.	▶ Check calculation. ▶ Check titration parameters. ▶ Check the titer for the titrant in use.
	pH electrode is defective.	▶ Calibrate the electrode. See related documentation ▶ Maintain electrode. See related documentation ▶ If necessary, replace it.
	The glassware is dirty.	▶ Clean the glassware. See Chapter 10.2 "Cleaning and servicing the sample tube", page 79

Problem	Possible cause	Action
Measured nitrogen content is too low	Incomplete digestion.	<ul style="list-style-type: none"> ▶ Increase digestion time. ▶ Use different Kjeldahl tablet.
	The H ₂ SO ₄ volume is too low.	▶ Use KjelOptimizer software to optimize the digestion (download available on BUCHI website).
	Kjeldahl Tablets and H ₂ SO ₄ in wrong ratio.	<ul style="list-style-type: none"> ▶ Correct ratio of Kjeldahl Tablets and H₂SO₄. ▶ Use KjelOptimizer software to optimize the digestion (download available on BUCHI website).
	Nitrogen content per sample tube is too high.	<ul style="list-style-type: none"> ▶ Not apply more than 200 mg Nitrogen per sample tube. ▶ Use KjelOptimizer software to optimize the digestion (download available on BUCHI website).
	Not enough NaOH or incorrect concentration of NaOH used (required is 32 %)	▶ Correct volume for complete alkalization of the digested sample.
	Leakage during distillation.	<ul style="list-style-type: none"> ▶ Check connection between condenser and splash protector. ▶ Tighten the connection. ▶ If necessary, replace it.
	Titrant solution	▶ Check titer of titrant.
	pH electrode is defective.	<ul style="list-style-type: none"> ▶ Calibrate the electrode. See related documentation ▶ Maintain electrode. See related documentation ▶ If necessary, replace it.
	The glassware is dirty.	▶ Clean the glassware. See Chapter 10.2 "Cleaning and servicing the sample tube", page 79
	Incorrect weighing.	<ul style="list-style-type: none"> ▶ Use weighing boats (easy sample transfer from balance to sample tube). ▶ Use anti-static equipment. ▶ Use larger sample sizes.

Problem	Possible cause	Action
Poor repeatability	Air bubbles in titration system, burette, tubes.	<ul style="list-style-type: none"> ▶ Tighten the connections. ▶ Flush the tubing with titrant. ▶ Refill burette.
	Aspiration not working properly.	<ul style="list-style-type: none"> ▶ Check for leaks. ▶ Tighten the connections.
	Sample is inhomogeneous.	<ul style="list-style-type: none"> ▶ Homogenize the sample.
	Sample weighing problems.	<ul style="list-style-type: none"> ▶ Use weighing boats (easy sample transfer from balance to sample tube). ▶ Use anti-static equipment. ▶ To decrease the degree of error, keep the weighed sample portion as high as possible.
	Incomplete digestion, digestion time too short.	<ul style="list-style-type: none"> ▶ Choose digestion time accordingly. ▶ Check color of samples during digestion. ⇒ Solution should be transparent by the end of the digestion.
	Stirrer is defective.	<ul style="list-style-type: none"> ▶ Clean the stirrer. ▶ If necessary, replace the stirrer.
	Loose contact of the sensor cables.	<ul style="list-style-type: none"> ▶ Check the sensor cables.


11.3 Troubleshooting instrument

Problem	Possible cause	Action
The instrument does not work	The installation site has no power supply.	<ul style="list-style-type: none"> ▶ Check the power supply of the installation site.
	The instrument is not connected to the power supply.	<ul style="list-style-type: none"> ▶ Connect the instrument to the power supply. See Chapter 5.2 “Establishing electrical connections”, page 29
	The power supply cable is defect.	<ul style="list-style-type: none"> ▶ Replace the power supply cable.
	The fuse was triggered.	<ul style="list-style-type: none"> ▶ Reset the fuse
	The switch is defect.	<ul style="list-style-type: none"> ▶ Contact BUCHI Customer Service.

Problem	Possible cause	Action
Steam generator is not reaching the status <i>Ready</i>	Not enough water in the steam generator.	<ul style="list-style-type: none"> ▶ Make sure, that the installation H₂O supply for steam generation is correct. ▶ Make sure, that there is enough water in the canister. ▶ Make sure, that the hose is submerged in the water. ▶ Check for leaks between instrument and canister. ▶ Check water level status in the steam generator.
Display is black	The connection between instrument and display is interrupted.	<ul style="list-style-type: none"> ▶ Check the connection cable from the instrument to the display. ▶ Contact BUCHI Customer Service.
No cooling water flow	The cooling water supply is blocked.	<ul style="list-style-type: none"> ▶ Make sure, that the hoses are not bend. ▶ Check for leaks between instrument and cooling water source.
	Water flow sensor blocked.	<ul style="list-style-type: none"> ▶ Clean the hoses.
Dosing pump is not feeding	Not enough liquid for feeding.	<ul style="list-style-type: none"> ▶ Make sure, that the installation is correct. ▶ Make sure, that there is enough liquid in the canister. ▶ Make sure, that the hose is submerged in the liquid. ▶ Check for leaks between instrument and canister. ▶ Make sure that liquid is inside the pump.
Aspiration is not working	Leaks	<ul style="list-style-type: none"> ▶ Check the hoses connected with the pump for leaks and deterioration.
	The aspiration pump is defect.	<ul style="list-style-type: none"> ▶ Check if the wheel inside the pump is turning. ▶ Contact BUCHI Customer Service.

11.4 Tighten the sample tube sealing

- ▶ Turn the splash protector nut.

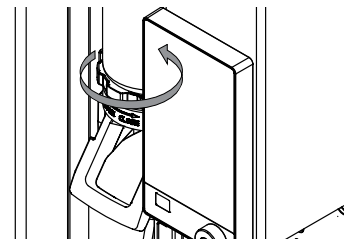

12 Taking out of service and disposal

12.1 Taking out of service

- ▶ Rinse all pumps. See Chapter 10.21 “Rinsing a pump”, page 86
- ▶ Remove the water from the steam generator. See Chapter 12.4 “Removing water from the steam generator”, page 93
- ▶ Remove all reagents and coolants.
- ▶ Clean the instrument.
- ▶ Set the On/Off master switch to Off.
- ▶ Disconnect the power supply.
- ▶ Remove all tubing and cables from the instrument.

12.2 Disposal

The operator is responsible for proper disposal of the instrument.

- ▶ When disposing of equipment observe the local regulations and statutory requirements regarding waste disposal.
- ▶ When disposing, observe the disposal regulations of the materials used. Materials used see Chapter 3.5 “Technical data”, page 24

12.3 Returning the instrument


Before returning the instrument, contact the BÜCHI Labortechnik AG Service Department.

<https://www.buchi.com/contact>


12.4 Removing water from the steam generator

Precondition:

- The On / Off master switch is set to Off.
- ▶ Install a drain hose to the connection marked *Steam Generator OUT*.
- ▶ Put the other end of the hose in a sink.

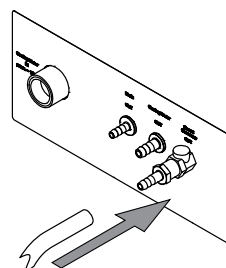







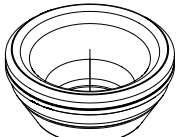
- ▶ Open the valve.

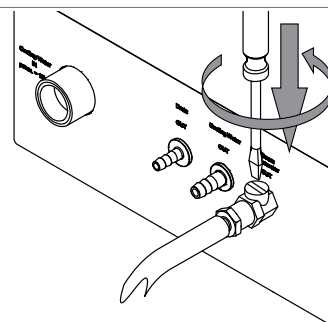


13 Appendix


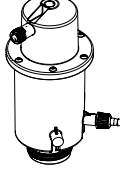

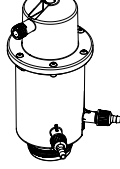

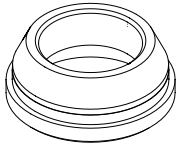
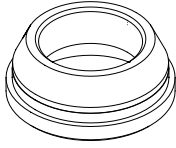
13.1 Spare parts and accessories

13.1.1 Accessories

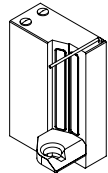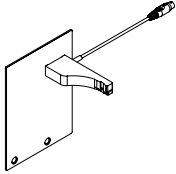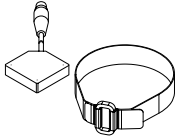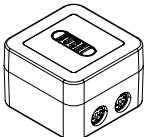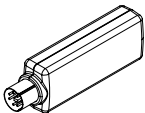
	Order no.	Image
<p>Reaction detection sensor</p> <p>Detects the optimal alkalization point, saves up to 30 % of NaOH in every determination, saves resources and reduces the ecological footprint</p>	11072666	
<p>OnLevel sensor</p> <p>Sensor to set the desired distilled volume in the receiving vessel, for most reproducible results in every determination</p>	11070270	
<p>Tank level sensor kit</p> <p>Detects the current reagent / water / waste level in the tank, for higher convenience and smooth operation during runs</p>	11072294	
<p>Eco Titrator</p> <p>Including pre-defined methods for MultiKjel / MultiDist and connection to enable automatic titration</p>	11072748	
<p>StatusLight cpl., incl. communication cable</p> <p>Indicates the status of the instrument (instrument is ready to use, has an error or is in operation).</p>	11068959	
<p>BUCHI Bluetooth® Dongle, connects instrument to smartphone via Bluetooth®</p>	11067770	
<p>Adapter for 3rd party sample tubes</p> <p>Adapter setup for other sample tubes (250 mL), sealing included</p>	11072398	
<p>Sealing for sample tube 3rd party adapter</p>	11072180	

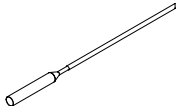	Order no.	Image
Condenser outlet for alcohol application	043096	







13.1.2 Splash protector spare parts

	Order no.	Image
Glass splash protector (EasyDist)	11071015	
Plastic splash protector (EasyDist)	11072055	
Glass splash protector For visible distillation processes.	11071013	
Plastic splash protector The most durable splash protector.	11070670	
Devarda splash protector Prevents from foaming into the condenser.	11071014	
Sealing with ring (CSM)	11073674	
Sealing with ring (CSM) 4 pcs.	11085448	






	Order no.	Image
Sealing with ring (EDPM)	11073675	A line drawing of a thick, circular O-ring or sealing ring with a slightly raised outer edge.
Steam inlet tubing to sample tube	043424	A line drawing of a thin, straight tube with a slight bend at one end.
Steam inlet tubing to sample tube (750 mL)	043119	A line drawing of a thin, straight tube with a slight bend at one end, similar to the previous one but possibly longer.
Connecting piece (CSM) Rubber piece between splash protector and steam outlet tube, alkali-resistant.	019002	A line drawing of a tapered, conical rubber fitting with a flange on the wider end.
Connecting piece (CSM) 4 pcs.	11085449	A line drawing of a tapered, conical rubber fitting, similar to the previous one.
Connecting piece EPDM Rubber piece between splash protector and steam outlet tube, acid-resistant.	11062478	A line drawing of a tapered, conical rubber fitting, similar to the previous ones.
Bridge splash protector to condenser	11070620	A line drawing of a cylindrical component with a flange and a textured, ribbed surface.
Cap nut GL14	033577	A line drawing of a cap nut with a flange and a textured, ribbed surface.
Hose barbs, set. 4 pcs, straight, GL 14, FPM seal Content: Hose barbs, cap nuts, seals.	040296	A line drawing showing a collection of small components: four hose barbs, two cap nuts, and several seals.

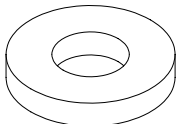	Order no.	Image
Seals, set. 10 pcs, for hose barbs GL 14, FPM, green	040040	



13.1.3 Tank spare parts

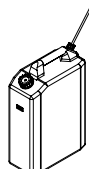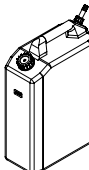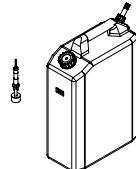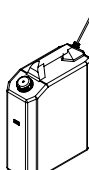
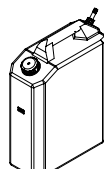
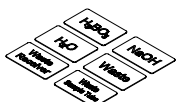

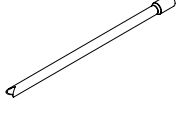
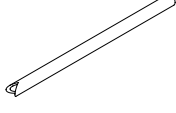
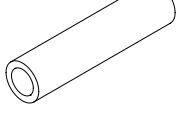
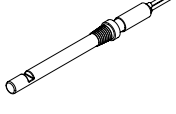
	Order no.	Image
Cap set, 10L tank	11073639	
Storage tank (10 L) for chemicals	043468	
Storage tank (10 L) for waste	043470	
Storage tank (10 L) with level sensor for waste	043474	
Cap Set, 20 L tank	11072174	
Storage tank (20 L) for chemicals	043469	
Caps included		
Storage tank (20 L) for waste	043471	
Caps included		
Tank labels	043434	


13.1.4 Condenser spare parts

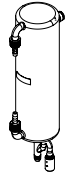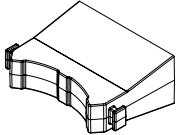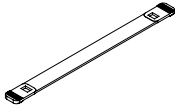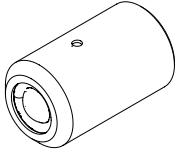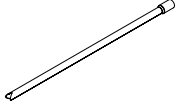	Order no.	Image
Glass condenser with check valve	11072183	
Clamp for condenser	11066868	
Rubber band Holding the glass condenser	11070669	
Check valve For overpressure safety	11071740	
Dist. Outlet PTFE L=300 for receiver side	11071940	



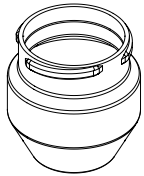
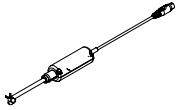



13.1.5 Titration related parts

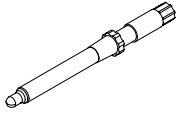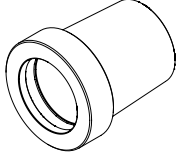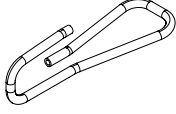
	Order no.	Image
Receiving vessel 420 mL	043390	
Stirrer For the receiving vessel.	11070246	
pH Electrode glass, refillable For potentiometric titrations.	11065834	

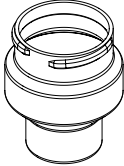	Order no.	Image
pH Electrode gel, maintenance-free For potentiometric titrations.	11056842	
Clamp cone pH-electrode	11069793	
Hose Tygon Ø 8.0 x 4.8 Hose to receiving module (waste aspiration, boric acid dosing)	043364	
Receiving support For attaching the receiving vessel to the distillation unit	11071003	
Condensate outlet to receiving module, PTFE	11071941	
Aspiration tubing receiving vessel	11072589	
H ₃ BO ₃ tubing receiving vessel For dosing H ₃ BO ₃	11072637	
Colorimetric sensor (Optrode) For colorimetric titrations.	11066601	
Kit colorimetric titration (Optrode) Includes all Optrode accessories (w/o Optrode sensor and receiving vessel).	11068260	

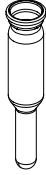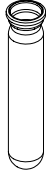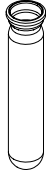
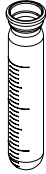



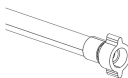
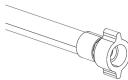
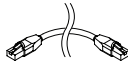
	Order no.	Image
Receiving vessel (Optrode) For colorimetric titrations with Optrode sensor.	11068263	

13.1.6 Sample tubes




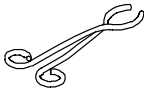
	Order no.	Image
Sample tubes micro (100 mL) 4 pcs.	11057442	
Sample tubes 300 mL 20 pcs.	11059690	
Sample tubes 300 mL For sample volumes up to 200 mL or 5 g in weight 4 pcs.	037377	
Sample tubes graduated 300 mL 4 pcs.	043049	
Sample tube 500 mL	026128	
Sample tubes 500 mL 4 pcs.	043982	
Set sample tube 750 mL with long suction tubing	11058999	

	Order no.	Image
Pair of glass tongs	002004	

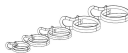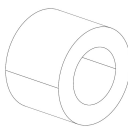13.1.7 Cable and tubing

Specification	Order no.	Image
Hose peristaltic pump out (2.5 m)	11071630	
Hose clips peristaltic pump hose	043586	
Hose acid pump (2 m)	043588	
Hose clip D7.6	043571	
Hose to NaOH/ H ₂ O/H ₃ BO ₃ pumps or drain (2 m)	11072687	
Hose clip to NaOH/H ₂ O/ H ₃ BO ₃ pumps or drain	043841	
Suction to tube tank (0.58 m)	043407	
Tap water hose cpl.	037780	
Set of sealings tap water hose	040043	
Connection cable to SI-Analytics TitroLine/ Titronic for K-365	043618	
Cable to Metrohm & Mettler (T-series) titrator To connect the Metrohm Titrimo plus, a Metrohm remote box is needed, see also quick guide.	043617	
Tubing to peristaltic pump	11070015	
Communication cable RJ45, 2 m Connection between vacuum controller/interface and recirculating chiller, vacuum controller/interface and vacuum pump or steam distillation unit and Eco titrator.	044989	

13.1.8 Other spare parts

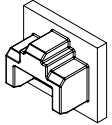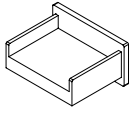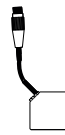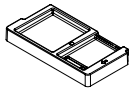	Order no.	Image
Cap RJ-45 cap	11055949	
USB cover	11069375	
Level sensor (capacitive) Part of tank level sensor kit	11065245	
Tank level sensor (Velcro point)	11070517	
Tank level sensor (Velcro strap)	11070516	
Collecting tray	11066465	

13.1.9 Consumables

	Order no.
2% boric acid pH 4.65 +/- 0.15 with Sher indicator, 5 L	11064972
4% boric acid pH 4.65 +/- 0.15 with Sher indicator, 5 L	11064973
4% boric acid pH 4.65 +/- 0.15 with bromocresol green / methly red mixed indicator, 5 L	11064976
Buffer solution pH 4.00 ± 0.01 twin-neck bottle, 250 mL	11064974
Buffer solution pH 7 ± 0.01 twin-neck bottle, 250 mL	11064975
KCl electrolyte, saturated Electrode storage solution, 250 mL	11064978
Sher indicator 100 mL Color indicator for colorimetric titrations.	003512
Ammonium phosphate monobasic, 25 g Reference substance for the verification and OQ of the Kjeldahl distillation units.	045527
Weighing boats Nitrogen-free weighing paper 58 x 10 x 10 mm 100 pcs	11060522
Kjeldahl Tablets Titanium For efficient digestions, 1000 tablets (3.5 g K ₂ SO ₄ / 0.105 g CuSO ₄ * 5 H ₂ O / 0.105 g TiO ₂)	11072627
Kjeldahl Tablets Titanium Micro For small sample sizes, 1000 tablets (1.5 g K ₂ SO ₄ / 0.045 g CuSO ₄ * 5 H ₂ O / 0.045 g TiO ₂)	11072628
Kjeldahl Tablets Missouri For efficient digestions without Titanium, 1000 tablets (4.98 g K ₂ SO ₄ / 0.02 g CuSO ₄ * 5 H ₂ O)	11072629
Kjeldahl Tablets ECO For eco-friendly digestions with minimum amount of copper, 1000 tablets (3.998 g K ₂ SO ₄ / 0.002 CuSO ₄ * 5 H ₂ O)	11072630
Kjeldahl Tablets Antifoam For foaming samples in digestions to avoid cross contamination due to foaming, 1000 tablets (0.97 g Na ₂ SO ₄ / 0.03 g Silicone)	11072631
Kjeldahl Tablets Copper Micro For small sample sizes without Titanium, 1000 tablets (1.5 g K ₂ SO ₄ / 0.15 g CuSO ₄ * 5 H ₂ O)	11072632

13.1.10 Maintenance kit

	Order no.
Customer Kit for Dist Line	11073025
Including suction tube with EPDM connecting piece, gaskets and EPDM seal with ring	

13.1.11 Upgrade kits

We offer the following upgrade kits:

- Upgrade Basic base unit with aspiration pump
- Upgrade Basic base unit and aspiration pump with I-300 Pro
- Upgrade Basic base unit with aspiration pump and I-300 Pro
- Upgrade Multi base unit with titration vessel

For ordering these upgrades please contact your local service support.

13.1.12 Documentation

	Order no.
IQ/OQ set MultiDist en	11073602
Including Eco Titrator if available	
Repeating OQ MultiDist en	11073603
Including Eco Titrator if available	
IQ/OQ set BasicDist en	11073631
Repeating OQ BasicDist en	11073632
Kjeldahl Knowledge Base (EN)	11595478
Comprehensive guide covering all aspects of the theoretical and practical know-how.	
Kjeldahl Practice Guide (EN)	11592548
Provides theoretical background information, useful hints and calculation tables for daily routine work	
Kjeldahl Practice Guide (DE)	11592547
Kjeldahl Practice Guide (CN)	11592549



11594193 | F en

We are represented by more than 100 distribution partners worldwide.
Find your local representative at:

www.buchi.com

Quality in your hands
