

# Installation Guide

## TKN Set Throughput

### Scope of Delivery

Product

TKN Set Throughput

Article No.

11085543

### Installation

1



1. Remove the plastic tubing between condenser and neutralization vessel from the scrubber K-415. Close the condenser with the cap nut.

2



1. Place the 5 L bottle below the scrubber and cover it with the screw cap with 2 connections GL14.
2. Connect both EPDM tubings to the hose barbs and then attach these to the GL14 connections in the screw cap.
3. Connect one of the open endings to the S35 condensate drain connector.

**3**

1. Attach the S35 condensate drain connector to the outlet of the condenser. Secure it with the catchpot clip.

**4**

1. Connect the other end to the neutralization vessel of the scrubber.
2. Ensure that the connections are tight (e.g. by the help of cable straps or hose clamps) and that all other ports of the connection are closed with an appropriate cap or plug.